


Firmware Update Guide

This procedure shows the screen shots of Lab Advisor B.02.07.


Latest firmware, tools and documents can be downloaded from the Agilent web site

[http://www.chem.agilent.com/ layouts/agilent/downloadFirmware.aspx?whid=69761](http://www.chem.agilent.com/layouts/agilent/downloadFirmware.aspx?whid=69761)


1. Open Lab Advisor and create new system..

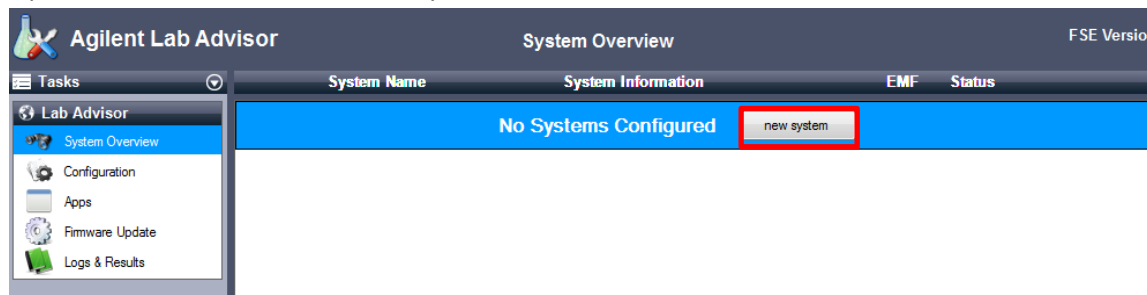

2. Enter **Instrument Name** and a valid IP address (default 192.168.254.11, fixed own IP or the MAC address of the LAN card/detector) and click on **OK**.

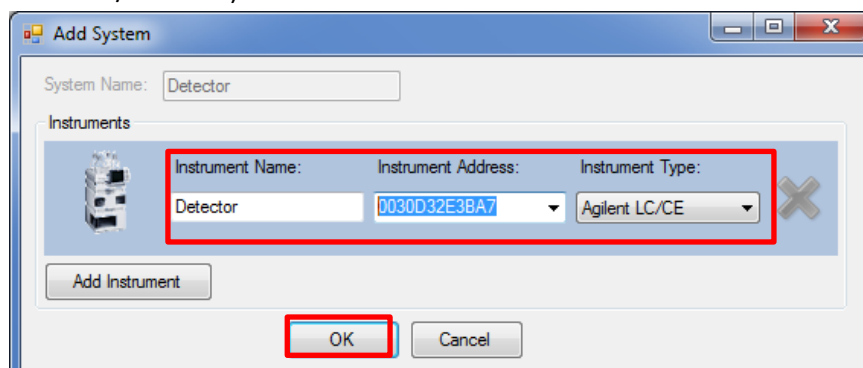

3. The System Overview shows the system with the modules. Click on **Firmware Update**.

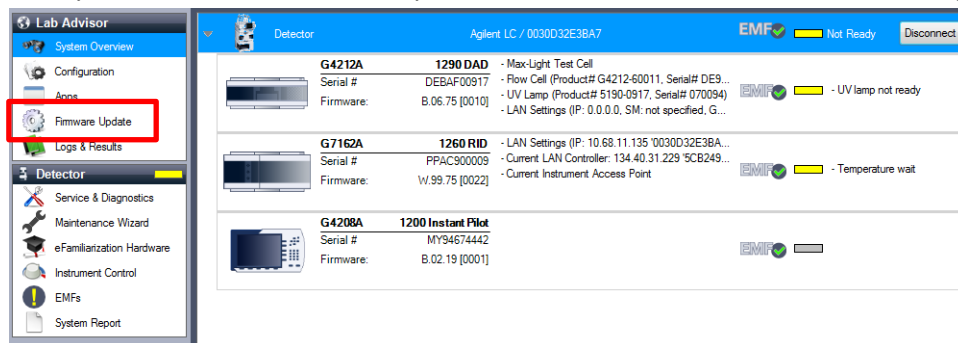

Firmware Update Guide

- Enter the **Firmware Update** and assure that the **firmware folder** is set correctly.

Caution: During the firmware update procedure, do not turn instrument off or disconnect AC power. Doing so may result in the instrument becoming unresponsive.

Detector [Update] Unlocked **Lock**

Device:	Resident Firmware:	Main Firmware:	Progress:
1290 DAD Serial # DEBAF00917		B.06.75 [0010]	
1260 RID Serial # PPAC900009		W.99.75 [0022]	
1200 Instant Pilot Serial # MY94674442		B.02.19 [0001]	

Firmware files are currently taken from: \\wadnas1\WAD-Firmware\Public\Firmware\Release-Archive\Release_A.06.50-B.06.50+Patches [Change folder](#)

- Lock** the system.
- Select the correct **Main Firmware** for all modules (if a newer resident version is available, update it as well) and click on **Update** to start the firmware update.

Detector [Update] Locked [Unlock]

Device:	Resident Firmware:	Main Firmware:	Progress:
G7162A 1260 RID Serial # PPAC900009 Keep current version	D.06.70 [0001]	W.99.75 [0022] Keep current version	
G4208A 1200 Instant Pilot Serial # MY94674442 Keep current version	B.02.19 [0001]	B.02.19 [0001]	
G4212A 1290 DAD Serial # DEBAF00917 B.06.70 [0001]	B.06.70 [0001]	B.06.75 [0010] B.06.75 [0010]	Update pending

Firmware files are currently taken from: \\wadnas1\WAD-Firmware\Public\Firmware\Release-Archive\Release_A.06.50-B.06.50+Patches [Change folder](#)

- When the firmware update has completed, the Lab Advisor shows **Finished**.

Caution: During the firmware update procedure, do not turn instrument off or disconnect AC power. Doing so may result in the instrument becoming unresponsive.

Detector [Update] Unlocked [Lock]

Device:	Resident Firmware:	Main Firmware:	Progress:
G7162A 1260 RID Serial # PPAC900009	D.06.70 [0001]	W.99.75 [0022]	
G4208A 1200 Instant Pilot Serial # MY94674442	B.02.19 [0001]	B.02.19 [0001]	
G4212A 1290 DAD Serial # DEBAF00917	B.06.70 [0001]	B.06.75 [0010]	Finished

Firmware files are currently taken from: \\wadnas1\WAD-Firmware\Public\Firmware\Release-Archive\Release_A.06.50-B.06.50+Patches [Change folder](#)

- Close the Lab Advisor.