

**Agilent UV-visible ChemStation and
Security Pack for Revision A.10.01****Site Preparation Specification****Purpose of Procedure**

To assure that the installation of Agilent instruments and systems can be completed successfully by careful preparation and evaluation of the installation site and by ensuring the availability of appropriate utilities, consumables and supplies.

Customer Responsibilities

Customers should ensure that all necessary operating supplies for the successful installation of instruments and systems are available. Installation sites should be prepared in accordance with the following specifications.

Important Information

If you have problems in providing any of the following, please contact your local Agilent office for assistance. Assistance with user specific applications may be provided but should be contracted separately. Users of the instrument should be present throughout the installation and familiarization otherwise important operational, maintenance and safety information may be missed.

Procedure Checklist**Minimum PC Hardware Requirements**

- Pentium III Processor 600 MHz
- Super VGA display (800 x 600 resolution)
- 4 GB hard disk
- CD-ROM drive

Minimum Memory Requirements

- 128 MB RAM for a single instrument configuration
- plus 32 MB for each additional instrument configured
- 256 MB RAM for a single-bath dissolution system (G1118AA) or kinetic measurements (G1117AA)
- plus 64 MB RAM for each additional bath in a multi-bath dissolution system (G1118AA)

Software Requirements Operating System

- Windows 2000 with service pack 3
- or**
- Windows XP with service pack 1a
- At least one NTFS partition
- User logon with administrative capabilities

ChemStation Software Requirements

- No other ChemStation software must be installed

**Additional Requirements for Workstation/Server
Installation**

- Make sure that the 'Prerequisites for the Server/Workstation Installation' of the G1813AA User's Guide are met
- or**
- Not applicable

Tick Boxes☐☐☐☐☐☐☐☐☐☐☐☐☐☐☐