
ChemStation E-Seminars

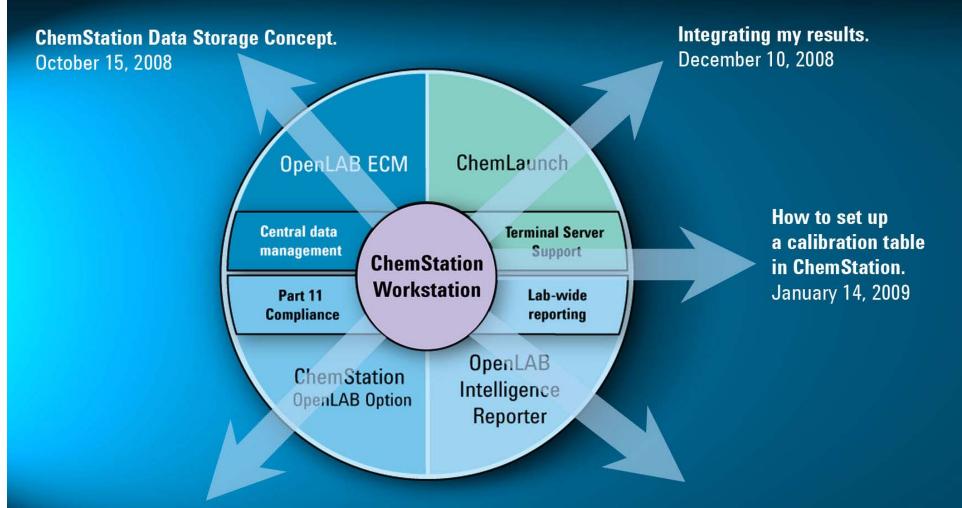

ChemStation Navigation Table — How to improve Review and Reprocess. November 12, 2008

Various Reporting in Chemstation. February 04, 2009



ChemStation E-Seminar Agenda

Schedule	Short description	Agilent Speaker
15 Oct 2008	ChemStation Data Storage Concept	Steven Brown Ortrud Emde
12 Nov 2008	ChemStation Navigation Table How to improve Review and Reprocess	Steven Brown Ortrud Emde
10 Dec 2008	Operational Functions for Data Generation, Data Analysis and Instrument Configuration	Steven Brown Ortrud Emde
14 Jan 2009	How to set up a calibration table in ChemStation	Steven Brown
04 Feb 2009	Various Reporting in Chemstation	Steven Brown Ortrud Emde

Seminar 1: ChemStation Data Storage Concept

Topics

ChemStation User Interface Enhancements Rev. B.02.01 or higher

The ChemStation "Preferences"

ChemStation Data Structure using Preference settings "Unique Folder Creation ON/OFF"


Advantages of the Sequence Container Concept "Unique Folder Creation ON"

Partial Acquisition Sequence in ChemStation B.04.01

Benefiz of OpenLAB with ChemStation


User Interface Enhancements ChemStation Rev. B.02.01 or higher


Method and Run Control - User Interface

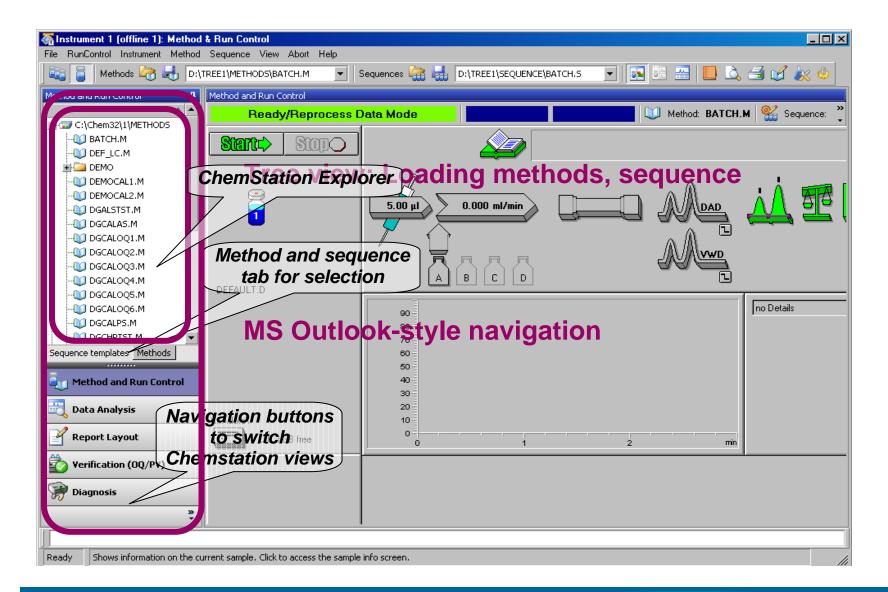


March. 2008

ChemStation Explorer - Views

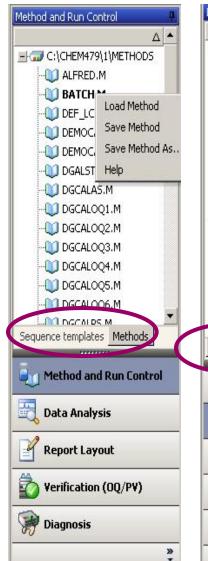

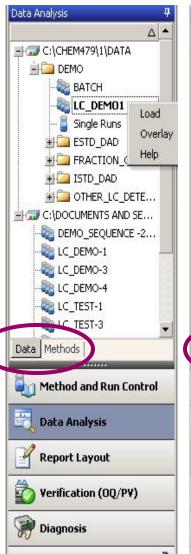


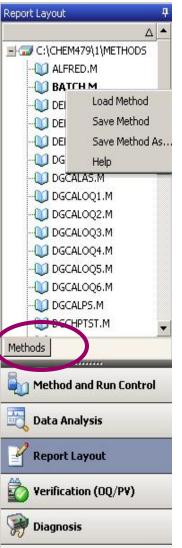







Navigation Pane – Customization

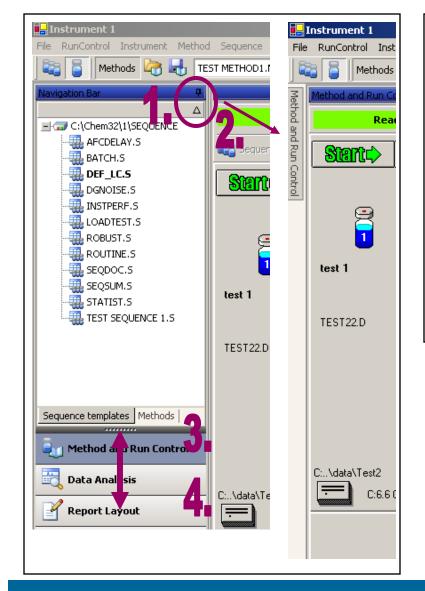


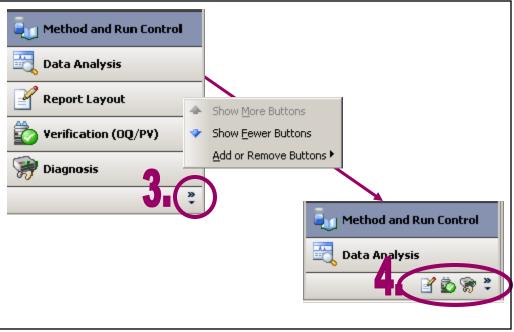

- ➤ Recommended Screen resolution: 1280x1024
- ➤ Customization features
- ➤ Sorting features

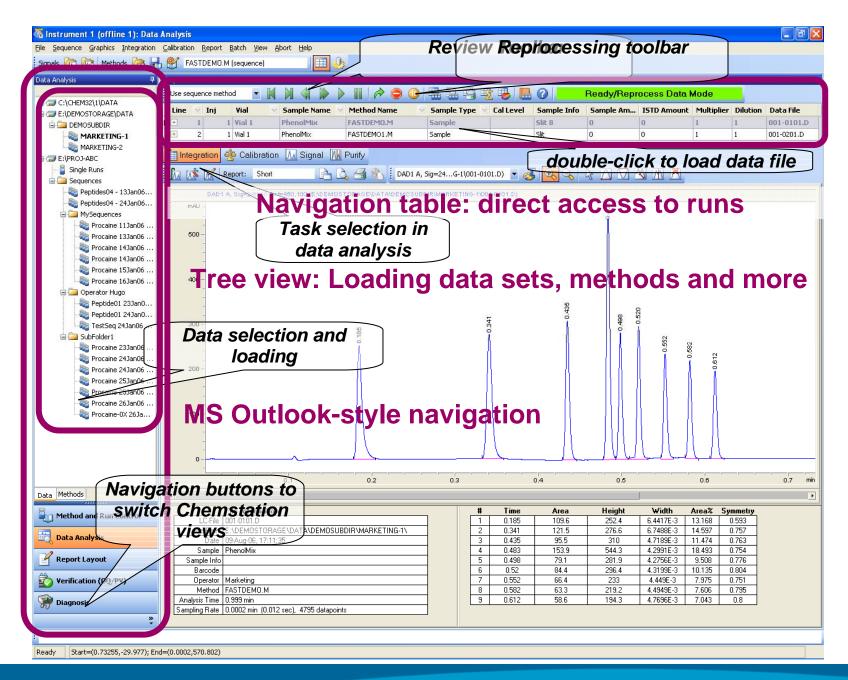


Loading Sequences and Single Runs from ChemStation Explorer

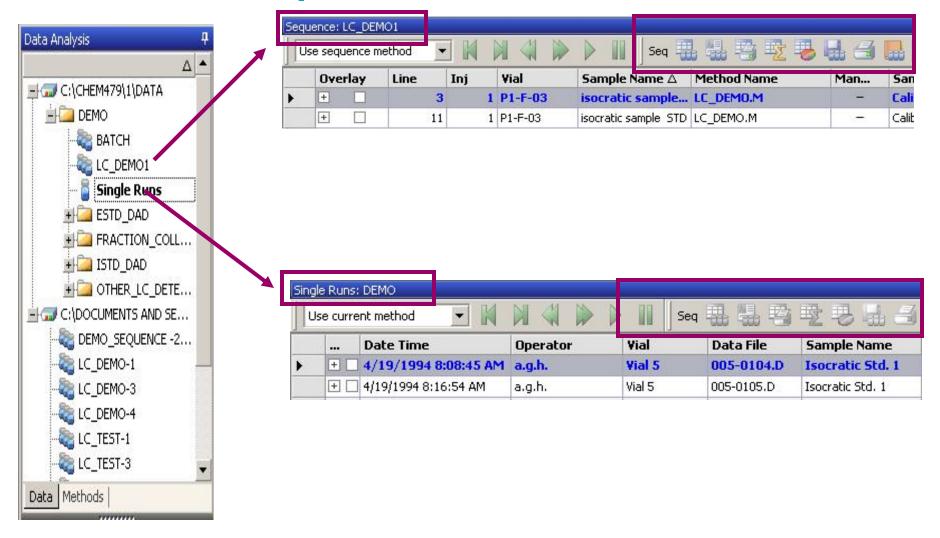

The ChemStation "Preferences"


Preferences Dialog – Flexible Path Settings

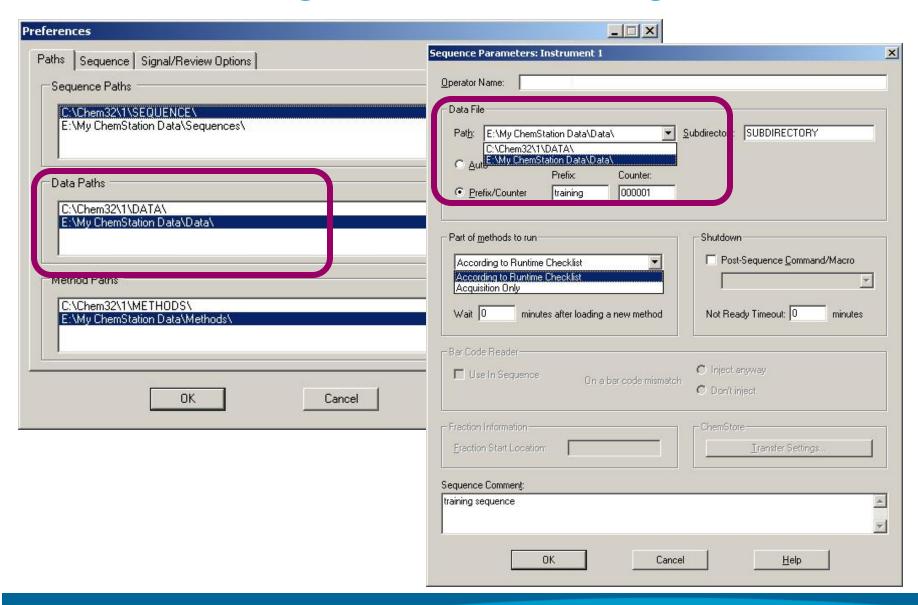

Choose your Data Storage Model

Unique Folder Creation OFF:


- ChemStation only
- Data are stored at a defined directory, the methods and sequence used to generate the data remain at their original location
- Possible to overwrite data
- Easy data review sequence possible
- Reprocessing of a complete sequence needs to be performed in Method and Run Control

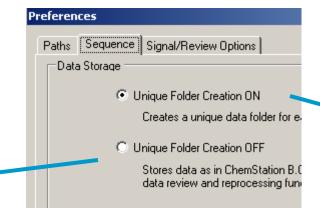

Unique Folder Creation ON:

- With ECM integration or ChemStore
 - Ensure data integrity:
 store related methods and
 sequence templates
 with sequence data
- Each sequence creates
 a uniquely named sequence container
 - Not possible to overwrite data
 - Easy data review or reprocessing of a complete sequence possible


Preferences Dialog: Sequence Tab

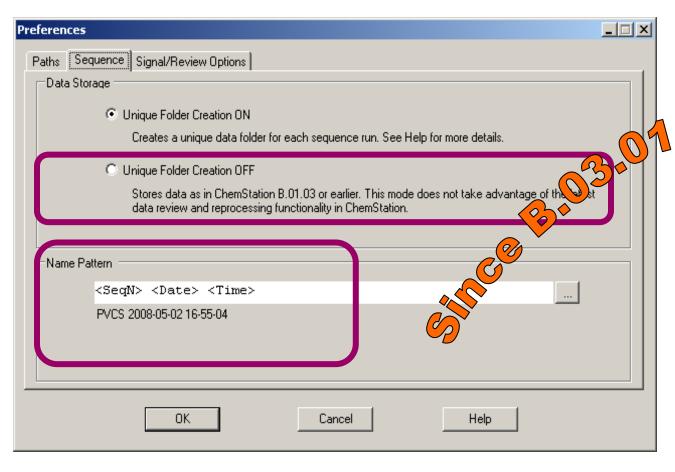

- Present in Online ChemStation only
- This setting influences only future data acquisition.

Sequence Name Pattern is grayed out when Unique Folder Creation is off.

Preferences Dialog: Sequence Tab

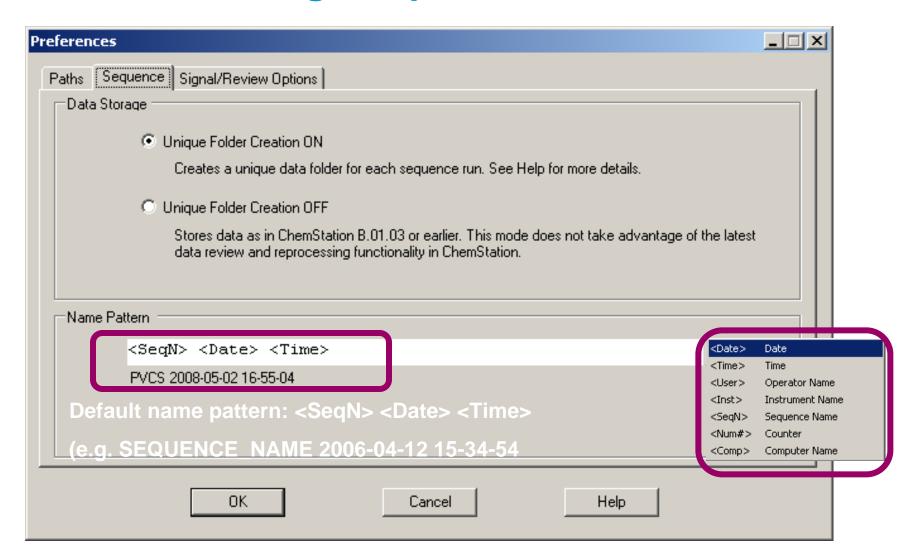

ChemStation Data Structure using Preferences Setting "Unique Folder Creation ON/OFF"

Comparison Start Prerequisites

ChemStation Rev. A – B.01.03 | ChemStation Rev. B.02.01

Methods: Master Methods:

C:\Chem32\1methods*.M C:\Chem32\1methods*.M

Sequence: Sequence Templates:

C:\Chem32\1\sequence*.S C:\Chem32\1\sequence*.S

Data Structure: Data Structure:

C:\Chem32\1\data\subdirectory\ C:\Chem32\1\data\subdirectory\

Press "Start Sequence" – What happens now?

Comparison - Single Run Acquisition

ChemStation Rev. A – B.01.03

ChemStation Rev. B.02.01

Run Method

Start the sequence – impact of the Data Storage chosen

Master Methods: C:\Chem32\1methods\Method1.M \Method2.M

Sequence Template: C:\Chem32\1\sequence\Sequence1.S

Data Structure: C:\ Chem32\1\data\subdirectory\


Chemstation Rev. A . B.01.03 Data1.D Data2.D

Data3.D

Data4.D


Data5.D

Sequence1.log Sequence Logbook Sequence1.b Sequence Batch File

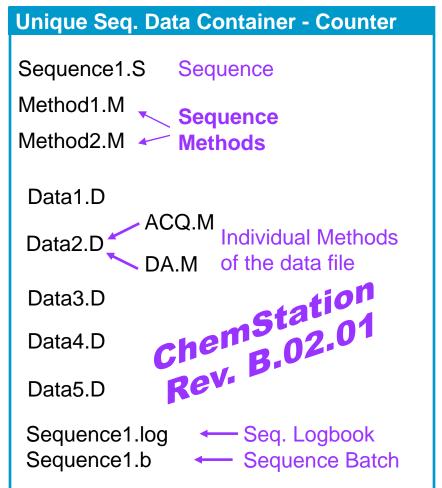


ChemStation - Data Integrity


- Unique Sequence Data Container
- No data is ever overwritten
- Data are filed along with relevant methods and sequence template
- Container includes sequence logbooks
- Each data file contains its own run logbook

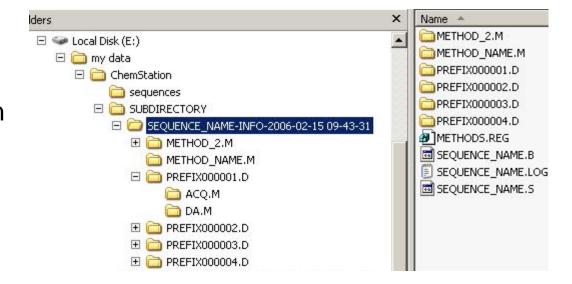


Saving the Method

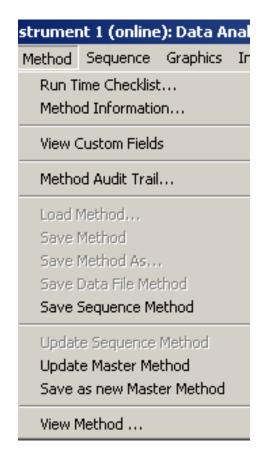

- Save Sequence Method:
 - Save the method with the same name, overwrite the previous parameters.
- Save as new Master Method:
 - Allows saving the method with a different name and to a different location; default target location is the master methods folder.

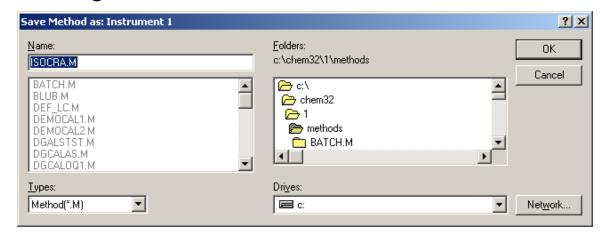

Saving the Sequence File

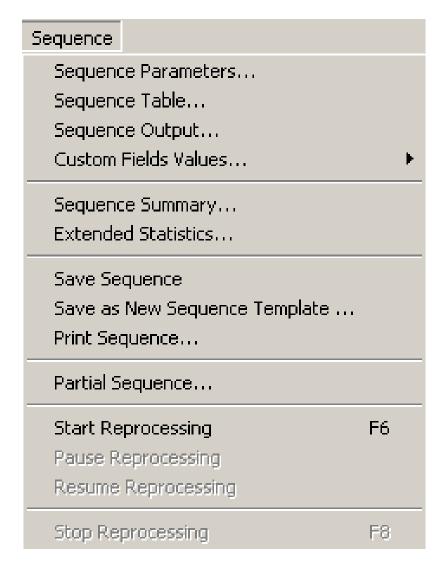

Save Current Sequence:


 Save the sequence file with the same name, overwrite the previous parameters.

Save as New Sequence Template:

• Allows saving the sequence file as a sequence template; converts the file to enable acquisition and editing of acquisition parameters.

Advantages of the Sequence Container Concept

Impact on Data Analysis Tasks – Review and Reprocess

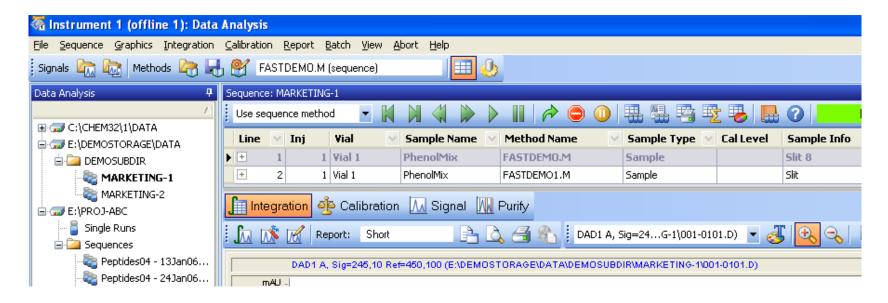

The data are stored with all necessary information needed for data analysis.

Your master method is now the sequence method of your newly acquired data container. All changes now are changes for the data set only, and remain only for this data set.

The sequence method used for the specific data file is automatically loaded.

Which Method is the Best to Use?


Working Methods:

For Data Acquisition – use the Master Method

For Data Analysis – use the Sequence Method

(which is a copy of your master method for the data set you are working on)

Master Methods:
C:\Chem32\1methods
\method A.M
\method B.M

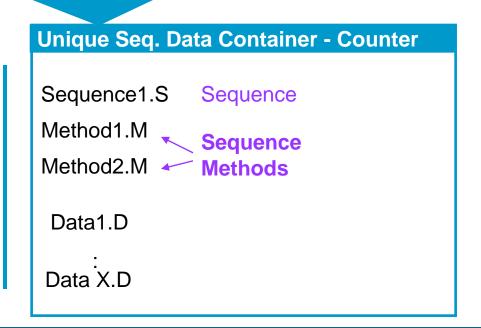

Unique Folder Creation switched OFF: Summary

- Methods are saved to the C:\Chem32\x\Methods directory
- > Sequences are saved to the C:\Chem\x\Sequence directory
- Data are saved to the C:\Chem32\x\Data directory
- Additional paths can be specified in the Preferences dialog
- ➤ Reprocessing is only possible from Method & Run Control Sequence Parameters\Part of the methods to run Reprocessing only
- Reprocessing is NOT possible from the Data Analysis view
- Data review IS possible from the Data Analysis view
- Possibility to perform partial acquisition into an existing folder
- Flexibility (for method development)
- Switching Unique Folder Creation off is not possible when ECM Integration or ChemStore are installed

March. 2008

New Data Organization - Key Advantages (Unique Folder Creation ON)

- All data is kept consistent, as related sets of data files, methods and sequences are kept together
- Data files created by the same sequence template are not accidentally overwritten, but automatically organized
- Results can be stored along with data files
- Easy data review and reprocessing of a complete sequence possible in Data Analysis
- Allows easier integration with data storage solutions such as Agilent ECM or ChemStore/Security Pack

March. 2008

History of Changes in the Sequence Container Concept

B.02.01	Introduction of sequence container concept	
B.02.01 SR1	Enhancements of UI in Data Analysis view	
B.02.01 SR2	Partial Sequence in Data Analysis view (Partial Reprocessing)	
B.03.01	Option to switch Unique Folder Creation off	
B.04.01	Partial acquisition into an existing sequence container	
	Storage of manual integration results in the data file	
	Introduction of a method menu in Data Analysis view	
	Possibility to store a sequence file inside a sequence container as a sequence template	


QUESTIONS?



Partial Acquistion Sequence in ChemStation B.04.01


Acquiring Data into an Existing Sequence Container (1)

Select Sequence / Partial Sequence

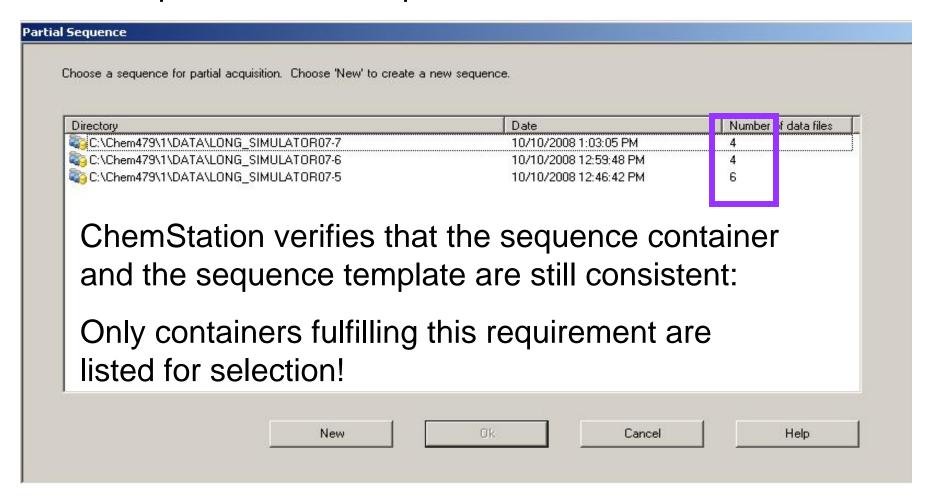

Acquiring Data into an Existing Sequence Container (2)

After either selecting an existing container or pressing New, the familiar Partial Sequence dialog comes up:

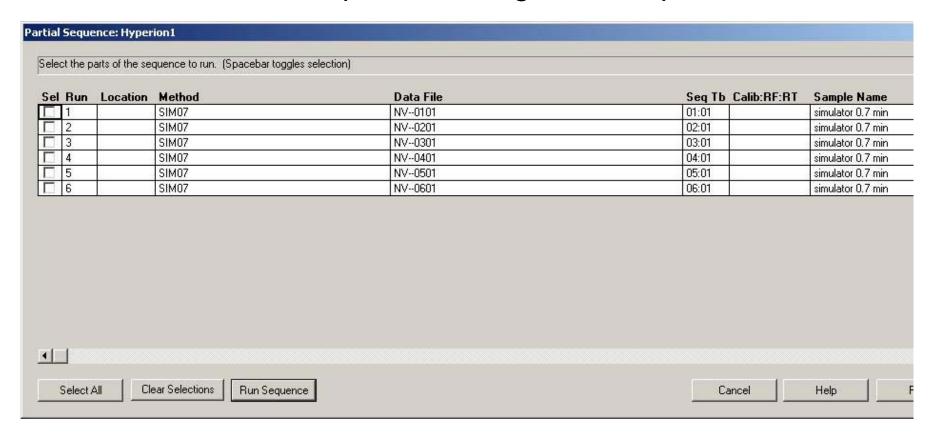


Information on ChemStation

Documentation can be found on the Agilent website http://www.chem.agilent.com/Scripts/PDS.asp?IPage=61233

- Product Datasheet
- Specification
- Application Notes
- Manuals

Manuals can be found on your ChemStation Installation DVD

- Getting Started with New ChemStation Workflow Part No. G2170-90042
- OpenLAB Option Part No. G2170-90233
- Configure and Maintain Your Agilent ChemStation Computer Part No. G2170-9013

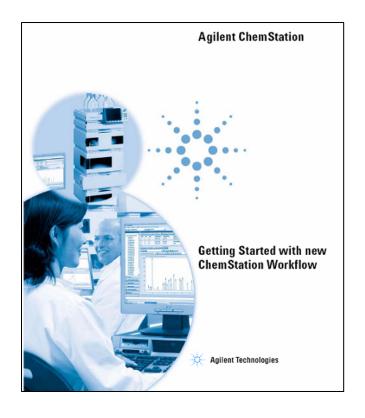

Software Status Bulletin

http://www.chem.agilent.com/en-US/Support/Downloads/Patches/Pages/default.aspx

Customer Trainings (NorthAmerica)

http://www.chem.agilent.com/cag/training/northam/index.asp

Learning Products – North America Course Catalog

Users needing to increase productivity in the lab by utilizing both standard and advanced features available in the Agilent GC/LC ChemStation may want to attend one of the following courses:

- ▶ H2606A ChemStation for GC Data Analysis and Reporting (2 days)
- ➤ H5928A Agilent HPLC (2D) Data Analysis and Reporting (2 days)
- ➤ H4039A Agilent HPLC (3D) Data Analysis and Reporting (3 days)

Course Features

- > Data acquisition and method creation
- > Data analysis including integration and calibration
- Sequencing
- Reporting

For more information concerning course content, dates and locations, please visit:

http://www.chem.agilent.com/en-US/education/en-us/classroomtraining/Pages/Courselisting.aspx

Learning Products – European Course Catalog

Users needing to increase productivity in the lab by utilizing both standard and advanced features available in the Agilent GC/LC ChemStation may want to attend one of the following courses:

- ➤ H4033A Agilent HPLC (3D) Method&Run Control, Data Analysis and Reporting (4 days)
- ➤ H8718A Agilent HPLC (3D) Data Analysis and Reporting (2 days)
- ➤ H5928A Advanced User Training, Quantification and Result Reporting (2 days)

Course Features

- Data acquisition and method creation
- > Data analysis including integration and calibration
- Sequencing
- Reporting

For more information concerning course content, dates and locations, please visit:

http://www.chem.agilent.com/en-US/education/pages/homepage.aspx

QUESTIONS?

