
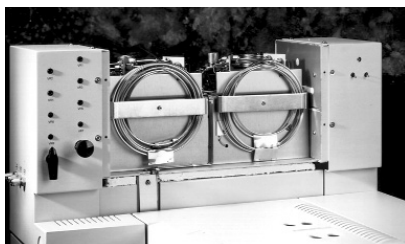


# Application 583D-00


## Agilent Fast Refinery Gas Analyzer

### Technical Overview

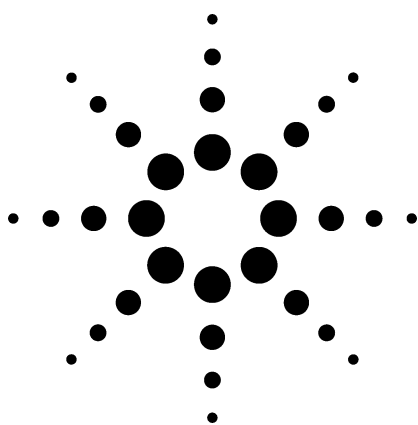



### Application Highlights

- Expect an analysis time of approximately 6 minutes.
- A Flame Ionization Detector (FID) to detect the C1 through C7 paraffins and olefins to a lower detection limit of 20 ppm, except for trace peaks eluting on the tail of a major component.
- A Thermal Conductivity Detector (TCD) is used to detect hydrogen in a nitrogen carrier to a lower detection limit of 100 ppm.
- A second TCD is used to detect carbon dioxide, ethane, ethylene, acetylene, hydrogen sulfide, O<sub>2</sub>/Ar, nitrogen, methane, and carbon monoxide to a lower detection limit of 200 ppm except for carbon monoxide (400 ppm), and hydrogen sulfide (500 ppm).

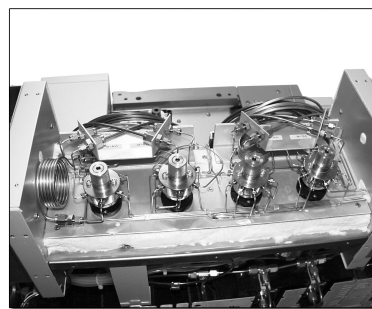

### Optional Configurations

- Refinery gas analysis with trace sulfurs by FPD or SCD
- Additional boiling point column for the analysis of heavy hydrocarbons (C1–C30)
- Standard analysis with the addition of trace CO by methanizer
- Custom analyzer for performing ASTM D2163, ASTM D1945, ASTM D2712, and ISO 7941
- High temperature injection for heavy fractions
- High temperature reactor effluent with percent level water
- TCD/TCD/MSD for the analysis of reactor effluent gases
- Valves for the injection of pressurized liquid samples.

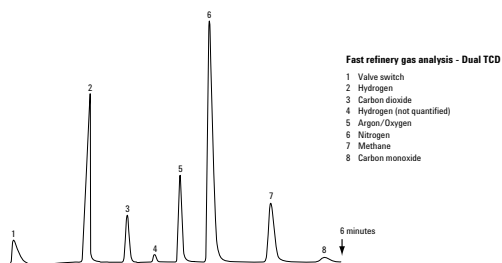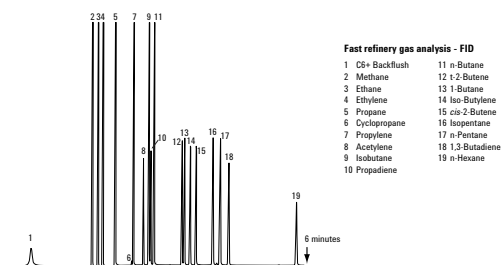

### For More Information


For more information on our products and services, visit our Web site at [www.agilent.com/chem](http://www.agilent.com/chem).



Agilent Technologies




WASSON·ECE  
INSTRUMENTATION

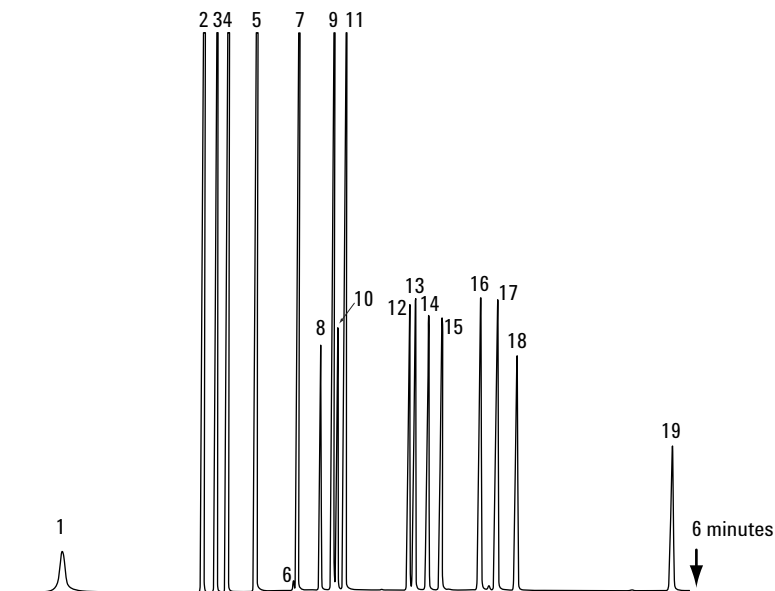

**Fast refinery gas analysis - FID**

- |                 |                         |
|-----------------|-------------------------|
| 1 C6+ Backflush | 11 n-Butane             |
| 2 Methane       | 12 t-2-Butene           |
| 3 Ethane        | 13 1-Butane             |
| 4 Ethylene      | 14 Iso-Butylene         |
| 5 Propane       | 15 <i>cis</i> -2-Butene |
| 6 Cyclopropane  | 16 Isopentane           |
| 7 Propylene     | 17 n-Pentane            |
| 8 Acetylene     | 18 1,3-Butadiene        |
| 9 Isobutane     | 19 n-Hexane             |
| 10 Propadiene   |                         |

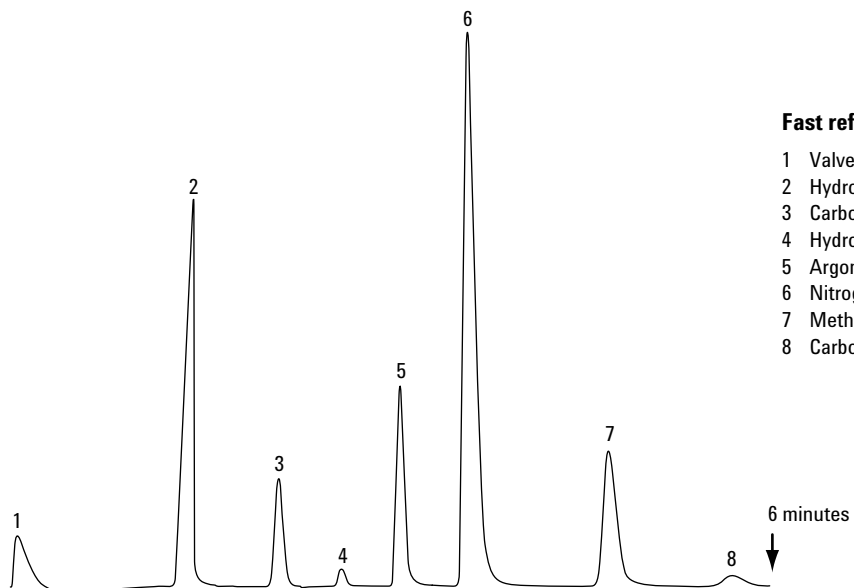

**Fast refinery gas analysis - Dual TCD**

- |                             |
|-----------------------------|
| 1 Valve switch              |
| 2 Hydrogen                  |
| 3 Carbon dioxide            |
| 4 Hydrogen (not quantified) |
| 5 Argon/Oxygen              |
| 6 Nitrogen                  |
| 7 Methane                   |
| 8 Carbon monoxide           |

**FID and TCD output from Agilent Fast Refinery Gas Analyzer.**

Agilent shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.

Information, descriptions, and specifications in this publication are subject to change without notice.

© Agilent Technologies, Inc. 2002

Printed in the USA  
December 6, 2002  
5988-6741EN