



Antiestrogen Drugs

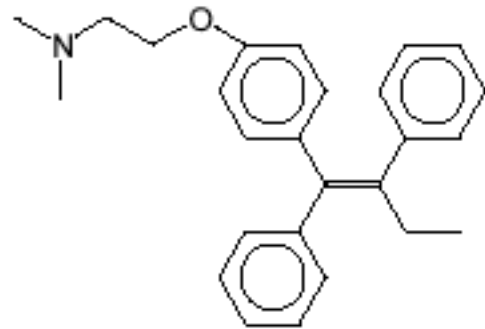

Tamoxifen

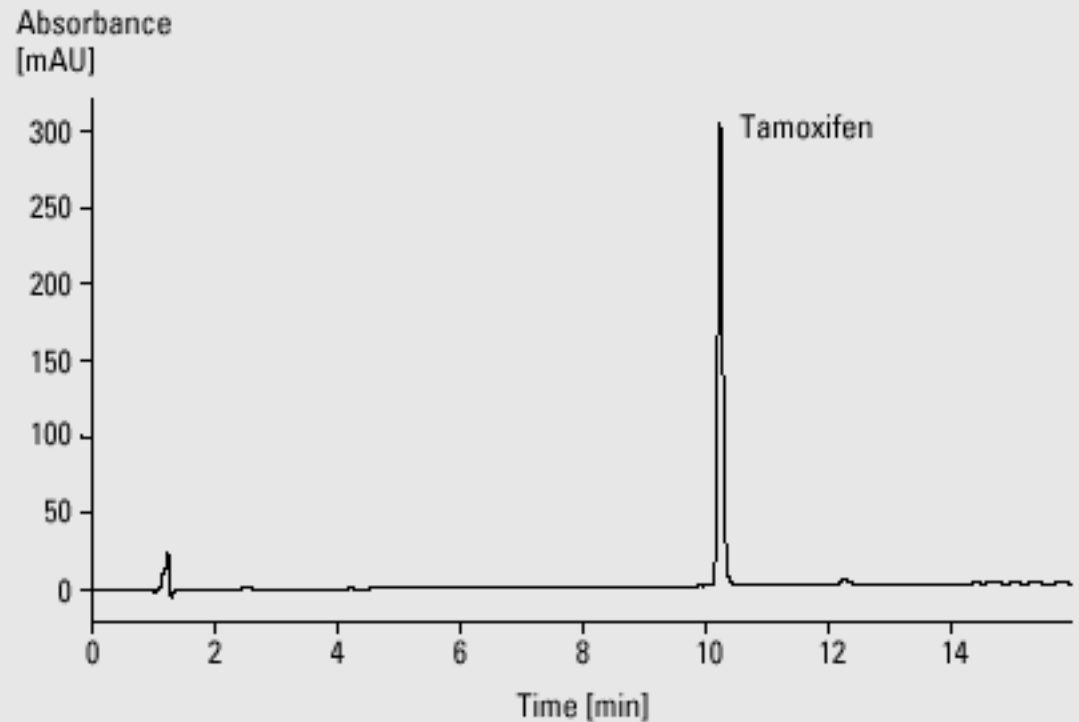

Analysis of antiestrogen drugs

Column	4 x 125 mm Purospher RP-18, 5 μ m
Mobile phase	A = 0.025 M KH_2PO_4 in water (pH = 3), B = acetonitrile
Flow rate	1.0 ml/min
Gradient	at 0 min 20 % B at 14 min 60 % B
Column wash	at 16 min 20 % B
UV detector	variable wavelength detector 204 nm, standard cell
Column compartment temperature	60 $^\circ\text{C}$
Stop time	16 min
Post time	5 min
Injection volume	5 μ l