

Fusel Oil Standard

DB-MTBE

124-0034

30 m x .45 mm I.D. x 2.55 μ m

DB-35

Custom column

30 m x .53 mm I.D. x 1.5 μ m

Inlet: 250°C, split

FID: 300°C

Carrier: H₂, 50 cm/sec

Oven: 40°C for 5 min.


10°C/min to 250°C

1. acetaldehyde

2. methanol

3. 3-methyl-butanol (isoamyl alcohol)

4. 2-methyl-butanol (active amyl alcohol)

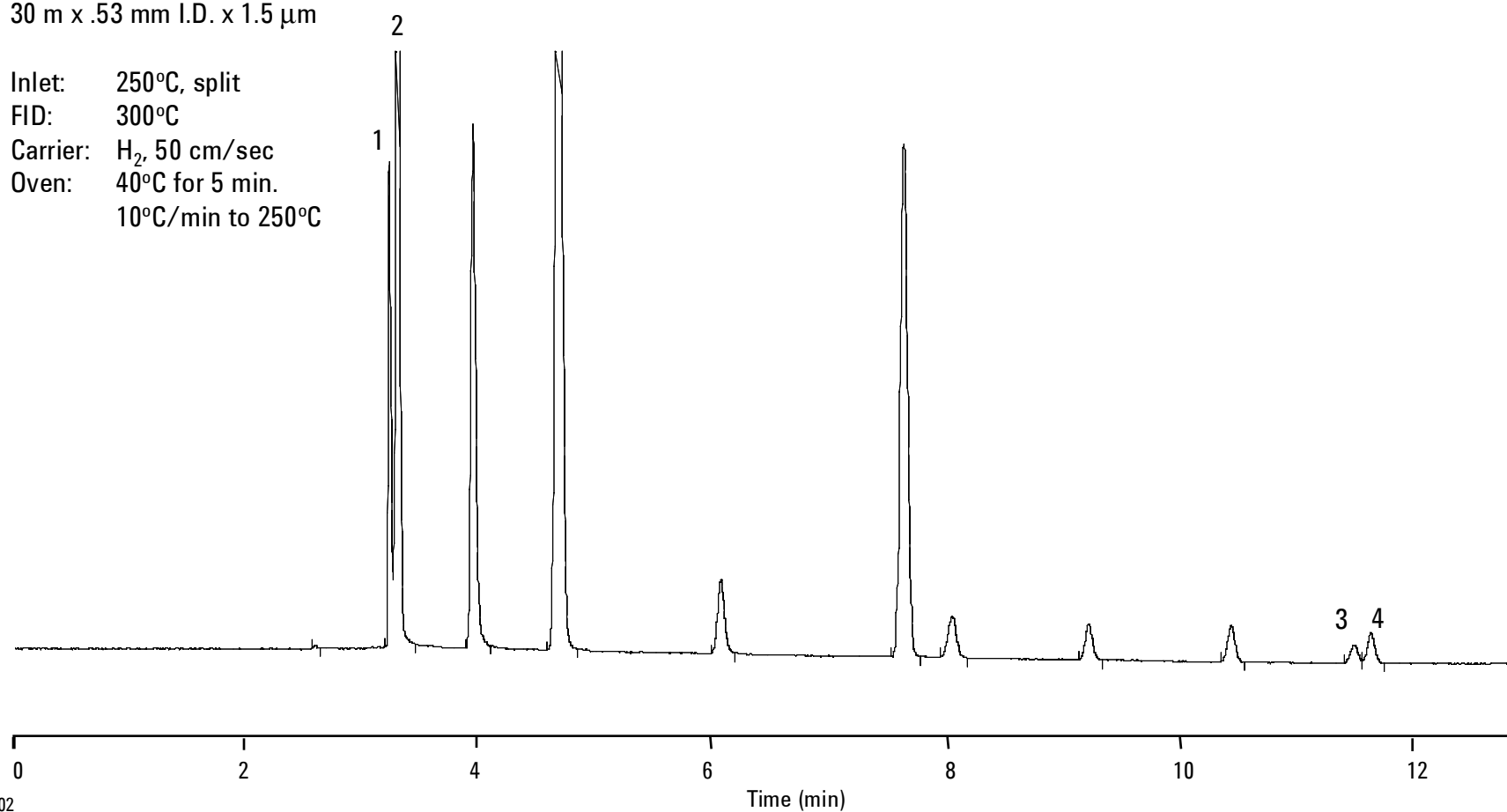

C1102

B-0457

Agilent Restricted



Agilent Technologies

Page 4