

CE Separation of Globin Chains With Urea

Column: μ SIL-FC
50 μ m I.D. x 75 cm (50 cm to detector)
J&W P/N: 291-8111 (MD-Phosphate Kit)
Buffer: 30mM phosphate, pH 3.0
+ 0.05% of FC-PN (J&W P/N: 590-4301) + 7M urea
Voltage: 25 kV
Injector: hydrostatic
Temperature: ambient
Detector: UV, 200 nm
Sample: diluted Hb standards (A, S, F, A₂) (Helena Lab)

