
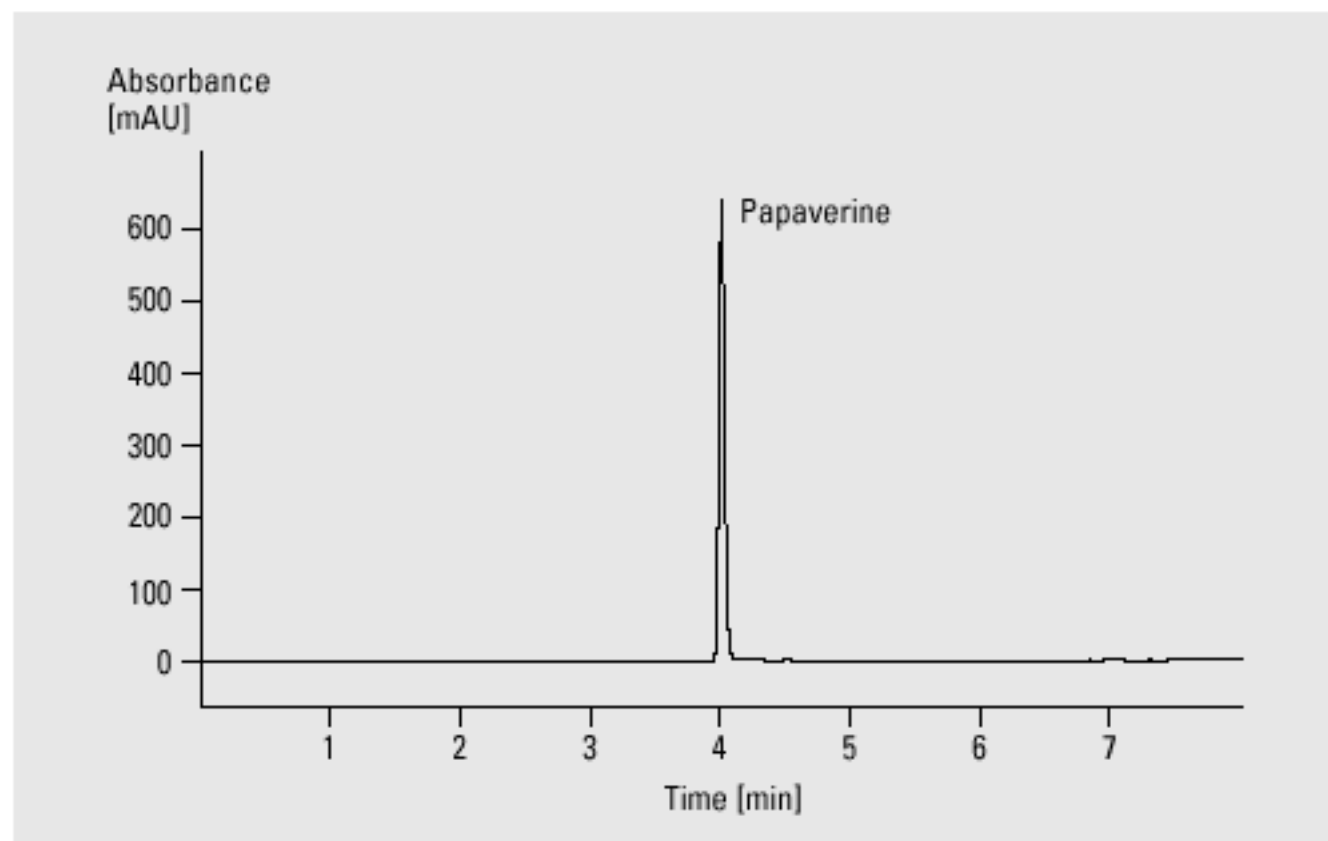


Muscle-relaxing Drugs

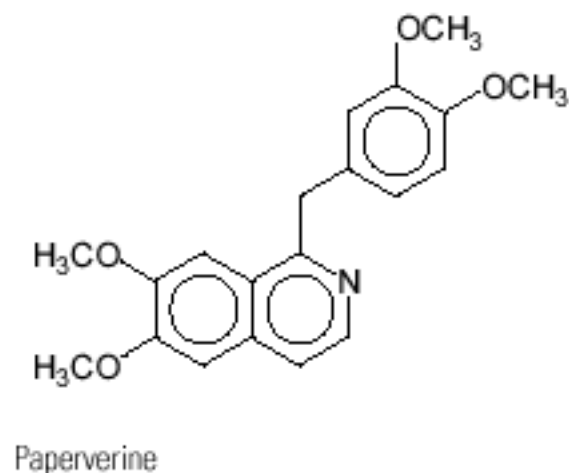

Analysis of muscle-relaxing drugs

Column	4.6 x 75 mm Zorbax SB-C18 , 3.5 μ m
Mobile phase	A = 0.025 M KH_2PO_4 in water (pH = 3), B = acetonitrile
Flow rate	1.0 ml/min
Gradient	at 0 min 10 % B at 8 min 80 % B
Column wash	at 10 min 10 % B
UV detector	variable wavelength detector 254 nm, standard cell
Column compartment temperature	60 $^\circ\text{C}$
Stop time	10 min
Post time	5 min
Injection volume	5 μ l