



**ZORBAX 300SB-C18**  
**Totally Porous Particle**

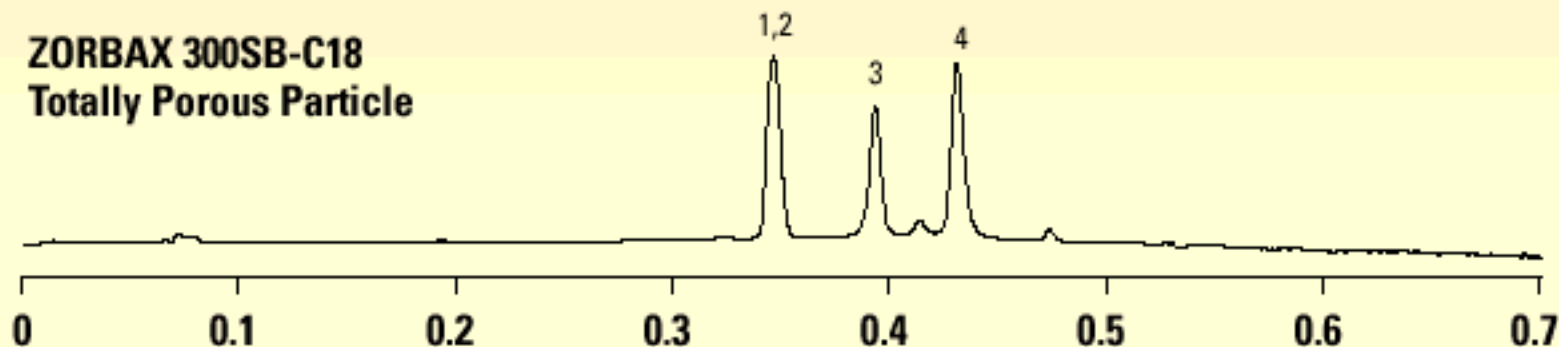

**Poroshell 300SB-C18**  
**Superficially Porous Particle**  
(Agilent Part No. 660750-902)

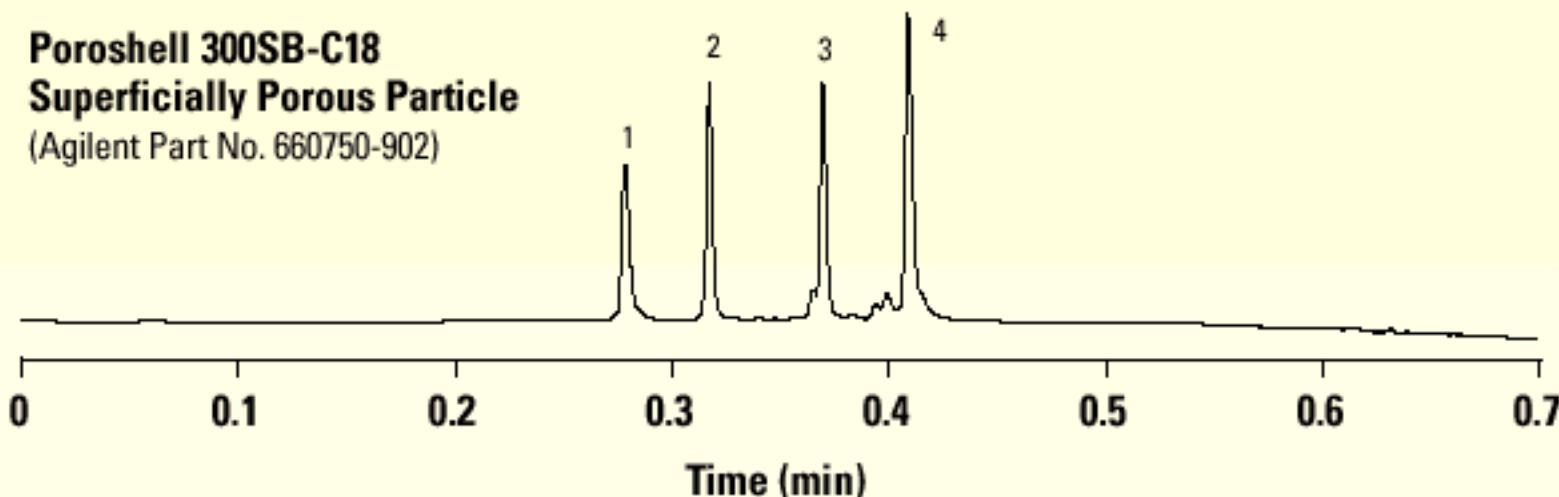

Column: 300SB-C18  
2.1 x 75 mm, 5  $\mu$ m

Mobile Phase:  
Gradient: 5 - 100% B in 0.67 min.  
A: 95% H<sub>2</sub>O, 5% ACN with 0.1% TFA  
B: 5% H<sub>2</sub>O, 95% ACN with 0.07% TFA

Flow: 3.0 mL/min  
Temperature: 70°C  
Detector: UV 215 nm  
Sample: 200  $\mu$ g each  
1. Neurotensin 3. Lysozyme  
2. RNaseA 4. Myoglobin