


Therapeutic Peptides

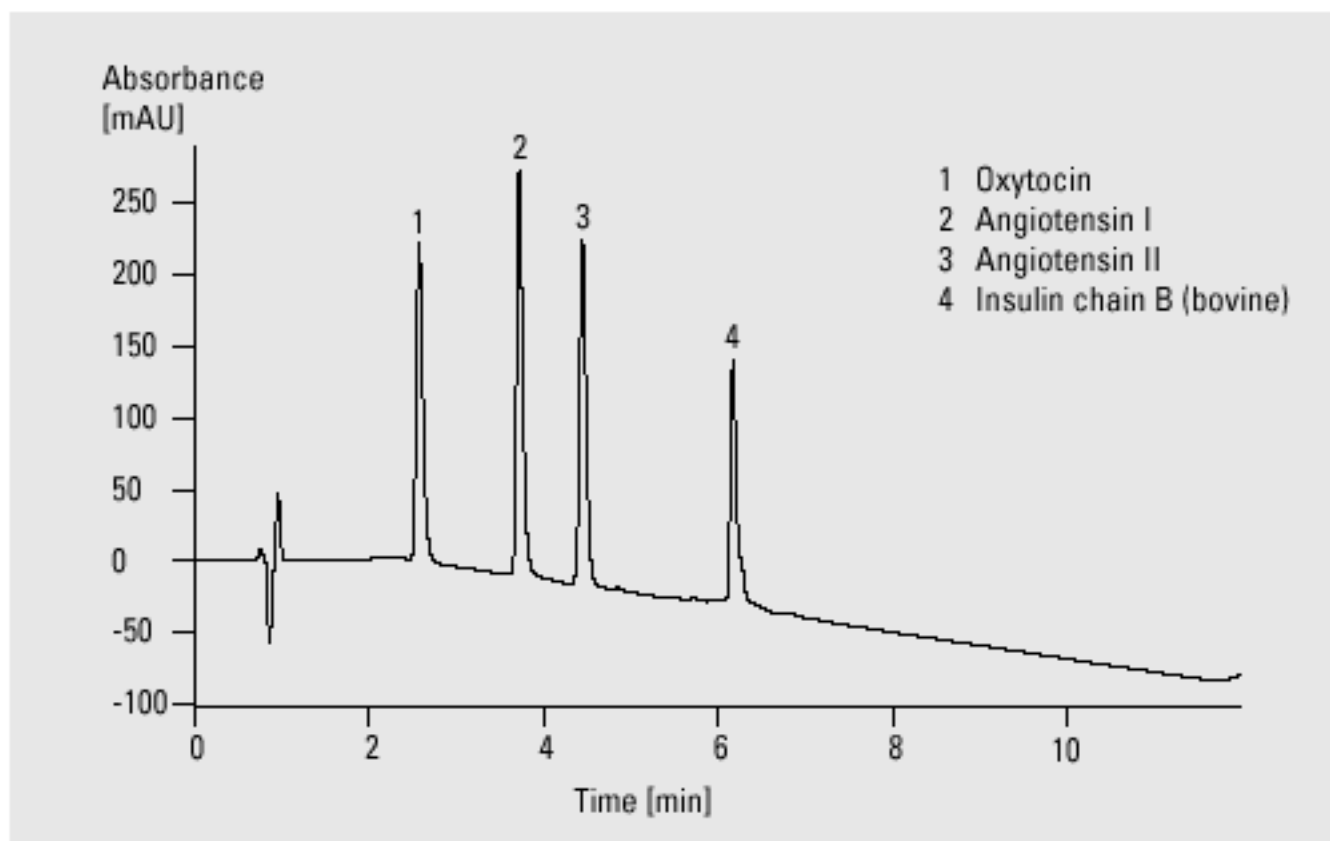

Analysis of therapeutic peptides

Column	4.6 x 75 mm Zorbax SB-C18 , 3.5 μ m
Mobile phase	A = 0.1 % TFA in water, B = acetonitrile
Flow rate	1.0 ml/min
Gradient	at 0 min 20 % B at 10 min 50 % B
Column wash	at 12 min 20 % B
UV detector	variable wavelength detector 210 nm, standard cell
Column compartment temperature	25 $^{\circ}$ C
Stop time	12 min
Post time	5 min
Injection volume	5 μ l