

Propylene Impurities Using GS-Alumina/KCl

Better resolution of isobutane and n-butane from the major propylene peak.

Column: GS-Alumina/KCl

50 m x 0.53 mm I.D.

J&W P/N: 115-3352

Carrier: Helium at 8 mL/min, measured at 35°C

Oven: 35°C for 2 min
35-190°C at 10°/min

Injector: Split 1:40, 200°C

0.2 mL propylene

Detector: FID, 200°C

Nitrogen makeup gas at 20 mL/min

1. Methane
2. Ethane
3. Ethylene
4. Propane
5. Cyclopropane
6. Propylene
7. Isobutane
8. Acetylene
9. n-Butane
10. Propadiene
11. trans-2-Butene
12. 1-Butene
13. Isobutylene
14. cis-2-Butene
15. Isopentane
16. n-Pentane
17. Propyne
18. 1,3-Butadiene
19. 1-Pentene

