


Impurities in Mixed Xylenes

Column: DB-WAXetr
60 m x 0.32 mm I.D., 0.25 μ m
J&W P/N: 123-7362
Carrier: Helium at 20 cm/sec,
measured at 145°C
Oven: 60°C for 10 min
60-150°C at 5°/min
150°C for 10 min
Injector: Split 1:150, 230°C
Detector: FID, 240°C



ASTM D2360

1. Benzene
2. Toluene
3. Ethylbenzene
4. *p*-Xylene
5. *m*-Xylene
6. *o*-Xylene
7. *n*-Butylbenzene (IS)

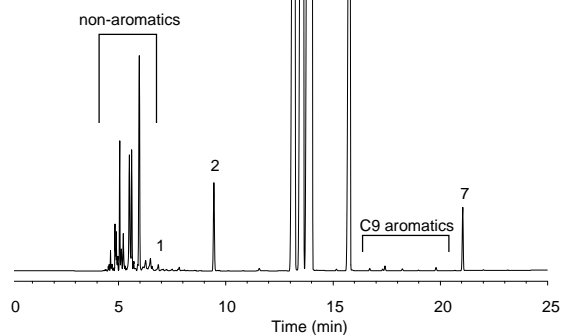

C590