

Acrylate Impurities on DB-200

Column: **DB-200**

30 m x 0.53 mm, 1.0 μ m

J&W P/N: **125-2032**

Carrier: Helium, 34.5 cm/sec at 35°C

Oven: 35°C for 5 minutes


10°/min to 200°C

Injector: Split, 1:10, 230°C

Detector: FID, 250°C

1. Methanol
2. Ethanol
3. Methyl acetate
4. Ethyl acetate
5. Methyl acrylate
6. Methyl propionate
7. Isobutanol
8. Butanol
9. Ethyl acrylate
10. Ethyl propionate
11. Methyl methacrylate
12. Isopropyl acrylate
13. Isobutyl acetate
14. Butyl acetate
15. Isobutyl acrylate
16. Dibutyl ether
17. Isobutyl propionate
18. Butyl acrylate
19. Butyl propionate

* Methyl acrylate impurities
• Ethyl acrylate impurities
◆ Butyl acrylate impurities

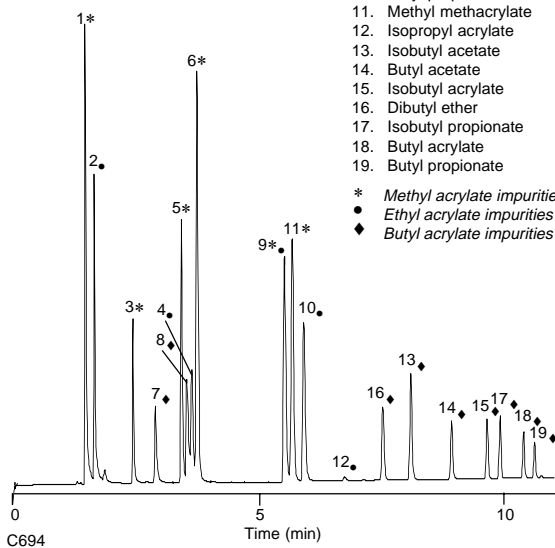

C694