

Agilent Lithography and Nanomanipulation Package, CAD-Based Software for the Agilent 5500, 5400, and 5100 AFM/SPM Data Sheet

Features and Benefits

- CAD-based drawing program offers user-friendly interface for easy operation
- Ability to design and save shapes for reuse helps conserve time and effort
- Ability to control parameters such as force or probe bias increases utility
- Ability to work “on-the-fly” or through scripts provides enhanced flexibility
- Move-tip tool provides real-time lithography
- Visual Basic® Scripting enables custom, programmable lithography
- Seamless integration with Agilent’s 5500, 5400, and 5100 AFM/SPM controller software offers simple upgrade path for AFM/SPM microscope users

Additional Capabilities

- Micro-machining via force variations
- Nano-patterning via probe bias and EC etching
- Nano-writing via surface potential controlled deposition and modification



Overview

Agilent’s lithography and nanomanipulation software is a CAD-based package that provides a nanolithography tool for scanning probe microscopy (SPM) researchers. It allows users to sketch various shapes on a canvas (including lines, poly-lines, circles, and boxes) that can then be mapped to a real sample surface. Individual shapes can be assigned a unique set of parameters. Each parameter set is associated with a color for easy visual reference.

Accessible parameters include force, probe bias, sample bias or potential, and probe speed. As shown in Figure 1, a user-friendly interface enables the user to edit parameters and segment properties, such as start and end points and the order of execution. The software offers users the ability to cut and paste predesigned shapes to assist in the building of more complex patterns. Once designed, shapes can be reused in future experiments and are easily resized.



Agilent Technologies

Move-Tip Tool Applications

The move-tip tool provides a complete set of functions for nanomanipulation. With numerous parameter settings, such as force set-point or tip bias, the user can inscribe a line, inscribe a pattern, or manipulate particles in real time. (See Figures 2 and 3.)

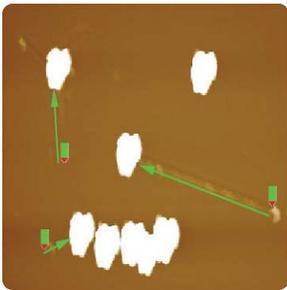


Figure 2. Making a smiling face with gold particles on a mica surface. Scan size: 2.5 μm x 2.5 μm .

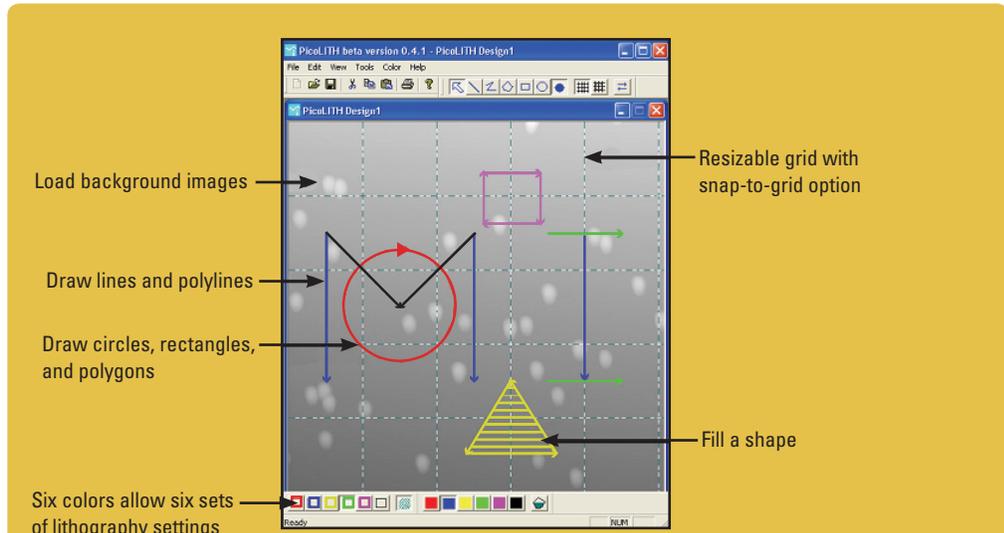


Figure 1a. Design canvas.



Figure 1b. Primitive property window.

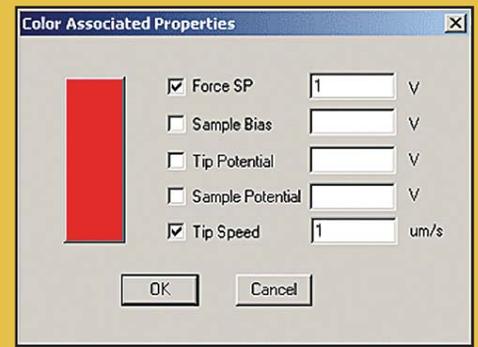


Figure 1c. Parameter settings window.

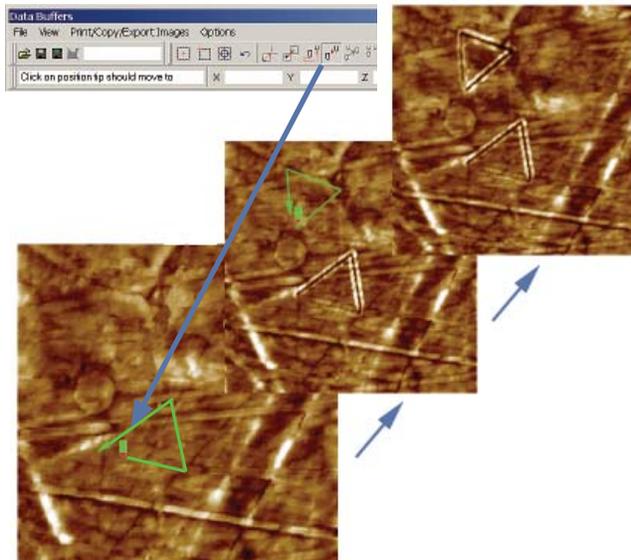


Figure 3. Sequence of drawing triangles on CD blank surface. Green lines indicate the path of the tip movement in contact mode. Images are taken in AAC mode. Scan size: 5.5 μm x 5.5 μm .

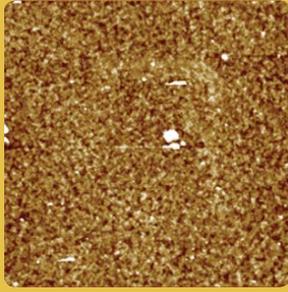


Figure 4a. AAC mode topography image of PZT film
Scan size: 14 μm x 14 μm

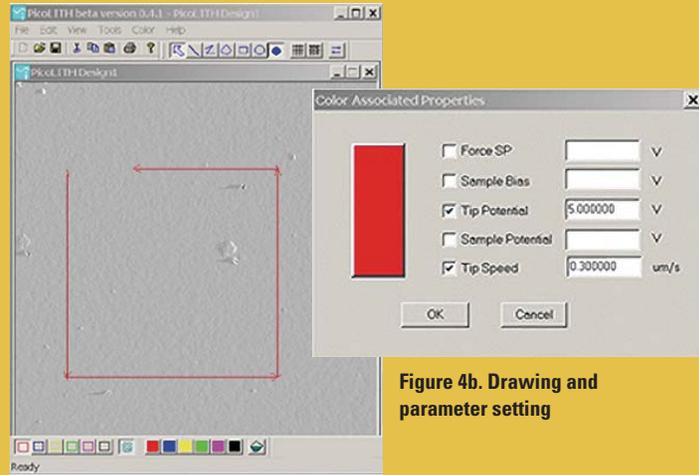


Figure 4b. Drawing and parameter setting

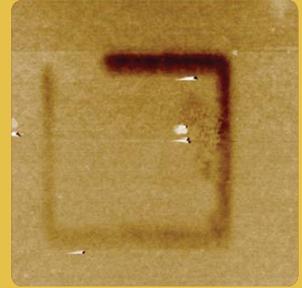


Figure 4c. EFM image after writing
Scan size: 14 μm x 14 μm .

Figure 4. Using Agilent's lithography and manipulation software to write with tip bias.

Visual Basic Scripting Applications

Agilent 5500, 5400, and 5100 AFM/SPM microscopes come with a 32-bit Windows® NT/2000/XP-based software package that incorporates Visual Basic Scripting capabilities. Visual Basic allows the user to program (via VBScript) most of the functions in the microscope software and define special sequences of desired action. By combining Agilent's closed-loop scanners and stages with VBScripting capability, the user is provided with a unique scripting tool for lithography applications. With predefined parameter settings (including STM/AFM tip bias, force set-point, and more), VBScripting makes drawing a line or a circle on the sample surface a straightforward procedure (Figure 4).

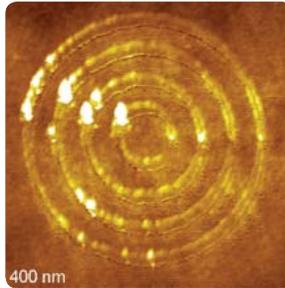


Figure 5. Drawing circles on PMMA film.
Scan size: 2.5 μm x 2.5 μm .

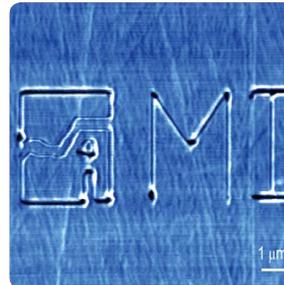


Figure 6. Simple writing on a blank compact disc surface.
Writing speed: 0.3 $\mu\text{m/s}$.
Force: 2 V (uncalibrated).
Scan size: 12 μm x 8 μm .

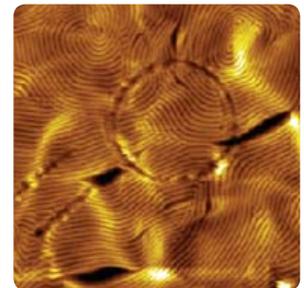


Figure 7. Inscribing a circle and lines on a polymer film by increasing the tip-sample interaction force.
Scan size: 5 μm x 5 μm .

AFM Instrumentation from Agilent Technologies

Agilent Technologies offers high-precision, modular AFM solutions for research, industry, and education. Exceptional worldwide support is provided by experienced application scientists and technical service personnel. Agilent's leading-edge R&D laboratories are dedicated to the timely introduction and optimization of innovative and easy-to-use AFM technologies.

www.agilent.com/find/afm

www.agilent.com

For more information on Agilent Technologies' products, applications or services, please contact your local Agilent office. The complete list is available at:

www.agilent.com/find/contactus

Phone or Fax

United States: (tel) 800 829 4444
(fax) 800 829 4433

Canada: (tel) 877 894 4414
(fax) 800 746 4866

China: (tel) 800 810 0189
(fax) 800 820 2816

Europe: (tel) 31 20 547 2111

Japan: (tel) (81) 426 56 7832
(fax) (81) 426 56 7840

Korea: (tel) (080) 769 0800
(fax) (080) 769 0900

Latin America: (tel) (305) 269 7500

Taiwan: (tel) 0800 047 866
(fax) 0800 286 331

Other Asia Pacific Countries:
tm_ap@agilent.com
(tel) (65) 6375 8100
(fax) (65) 6755 0042

Revised: 11/08/06

Product specifications and descriptions in this document subject to change without notice.

© Agilent Technologies, Inc. 2006
Printed in USA, December 20, 2006
5989-6004EN