

## MxPro™ QPCR Software 21 CFR Part 11 Compatible

- + Secure application login with unique username, encrypted password
- + Database file management for storing and tracking data
- + Audit trail tracks when files are created, modified, or archived
- + Report generation for audit trail, user accounts, and error logs



### MxPro™ Electronic Tracking (ET) QPCR Software

The MxPro™ ET software is built on Stratagene's existing MxPro software for the Mx3000P® and Mx3005P™ QPCR systems. The MxPro software is the graphical user interface for setting up, running, and analyzing data on the Mx™ QPCR Systems. This newest version of MxPro software uses an activation code to enable 21 CFR Part 11 compatible features. Four key features have been added to the ET version; secure application login, database file management, electronic audit trail, and report generation.

### Application Security

Access to the software is controlled through unique secure logins. Only the administrator has the ability to create, modify, and deactivate user accounts. This provides you access control to the application and to the database. In addition, the audit trail will track the user actions when creating or modifying a file. Each secure login is unique to ensure accurate electronic tracking in the application.

### Database File Management & Flexibility

The MxPro ET version includes a 2 GB Microsoft SQL® Server 2000 Desktop Engine (MSDE) database for storing and tracking data. The administrator has the option to use this database, create new databases, or use existing databases. Files can be easily imported and exported from the database and can be archived to other databases. This functionality is flexible and improves data organization, storage, and tracking.

### Audit Trail

Application security and database structure support the audit trail functionality. Each file has a date and time stamp with the associated action(s) performed on the file. All changes to a file are tracked to the user and only the administrator can view and print the audit trail.

### Report Generation

The administrator can generate reports for the audit trails, user accounts logs, and error logs. This creates a hard copy record of all the changes to a particular file or group of files.

For more information visit [www.stratagene.com/mxpro21cfrpart11](http://www.stratagene.com/mxpro21cfrpart11) or contact our Technical Service department at [qpcr@stratagene.com](mailto:qpcr@stratagene.com).

MxPro™ QPCR Software - 21 CFR Part 11 Compatible		Catalog No.
MxPro™ ET Software Activation*	New and existing Mx™ QPCR System users	401467
MxPro™ ET Software Multi-User** License Agreement	6 activations total	401468

#### Need More Information? Give Us A Call:

##### Stratagene US and Canada

Order: 800-424-5444 x3  
Technical Service: 800-894-1304 x2

##### Stratagene Europe

Order: 00800-7000-7000  
Technical Service: 00800-7400-7400

##### Stratagene Japan K.K.

Order: 3-5821-8077  
Technical Service: 3-5821-8076

[www.stratagene.com](http://www.stratagene.com)

\* Valid only for owners of Mx3000P® or Mx3005P™ QPCR Systems.

\*\* Valid only with purchase of Mx3000P® or Mx3005P™ QPCR System.

Mx3000P® is a registered trademark of Stratagene in the United States.  
MxPro™ and Mx3005P™ are trademarks of Stratagene in the United States.

Microsoft® and SQL® are registered trademarks of Microsoft.

For Research Use Only. Not for use in diagnostic procedures.

**STRATAGENE**

An Agilent Technologies Company

© Stratagene, an Agilent Technologies company, 2008  
Printed in USA, April 2008  
5989-8240ENUS