


Agilent ZORBAX Rapid Resolution High Definition Extend-C18 Column

Data Sheet

General Description

Agilent ZORBAX Rapid Resolution High Definition (RRHD) Extend-C18 threaded columns are specially designed for use with ultra high performance liquid chromatographs (UHPLCs) such as the Agilent 1290 Infinity LC and can be used up to an operating pressure of 1200 bar. Agilent ZORBAX Extend-C18 incorporates a patented bidentate organosilane combined with double endcapping to protect its ultra-pure (Type B) silica support from dissolution at high pH. Extend-C18 is specially designed for stable use with high pH mobile phases and is particularly well-suited for separating highly basic compounds as free bases. In addition, the structure of the unique bidentate-C18 bonded stationary phase also is useful for separations at low and intermediate pH with excellent stability and chromatographic separation properties. Extend-C18 packing is made by first chemically bonding a dense monolayer of propylene-bridged bidentate-C18 silane stationary phase to a specially prepared, ultra-high purity (99.995% SiO₂) ZORBAX Rx-SIL porous silica microsphere support. A schematic of the structure of this attached bidentate silane to the silica support is shown below.

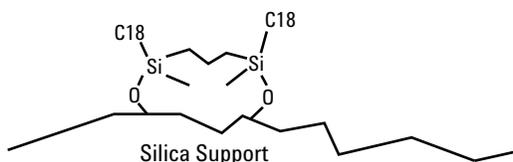

The bidentate-C18 bonded phase is double endcapped using proprietary reagents and procedures to obtain maximum deactivation of the silica support surface. The uniform, spherical Extend-C18 particles are based on ultra-high purity ZORBAX Rx-SIL that has a nominal surface area of 180 m²/g and a narrow controlled pore size of 80Å. This special ZORBAX silica support (Type B) is designed to eliminate or reduce strong adsorption of basic and highly polar compounds. Columns are loaded to a stable, uniform bed density using a proprietary high-pressure slurry-loading technique to give maximum column efficiency.

Column Characteristics

A typical quality control test chromatogram for a ZORBAX RRHD Extend-C18 2.1 mm id × 50 mm, 1.8-µm threaded column is shown in

Figure 1. The actual QC test and performance of your column is described on the Column Performance Report enclosed with your column. The efficiency found on the QC Performance Report may be higher than the efficiency found in your laboratory. The QC test system may vary from the LC used in your lab and has been modified from a standard system to minimize system volume. This allows a better evaluation of the column and assures a more consistent product for the chromatographer.

Safety Conditions

- All points of connection in liquid chromatographic systems are potential sources of leaks. Users of liquid chromatographic equipment should be aware of the toxicity or flammability of their mobile phases.
- These RRHD assembled columns are mechanically stable and have been tested to very high pressures to assure safe lab operation on a variety of LC instruments. The 2.1 and 3.0-mm ID columns will support 20,000 psi (1300 bar) operation. Opening columns will compromise these pressure limits.
- Because of the small particle size, dry ZORBAX packings are respirable. Columns should only be opened in a well ventilated area.

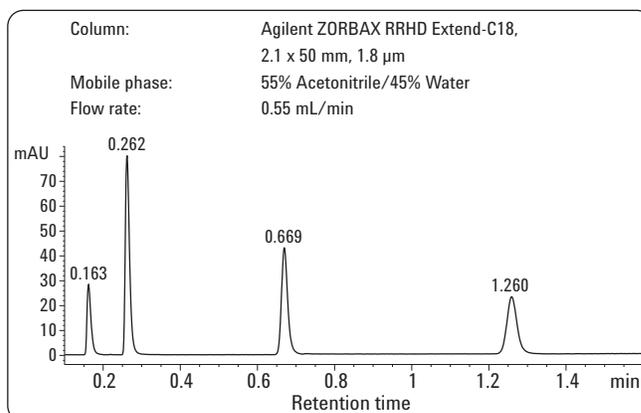


Figure 1. Agilent ZORBAX RRHD Extend-C18 QC chromatogram.



Agilent Technologies

pH Guidelines

The bidentate bonding to the silica support strongly contributes to the stability of the stationary phase at all pH values; hydrolysis at low pH is inhibited and degradation at intermediate and high pH by silica dissolution is moderated. When using organic buffers, it has been documented that the longevity of such a column at pH 11.0 can exceed 30,000 column volumes without significant separation degradation. This stability is similar to traditional silica-based columns, but only when they are operated at low or intermediate pH. Although this column is primarily proposed for high pH separations, it is highly competitive for low and intermediate pH applications. The Extend-C18 column has been used up to pH 11.5 with good results, providing certain operating parameters are followed. For high pH separations, organic buffers should be used and the operating temperature should not exceed 40 °C. Highest column lifetime is obtained at lower temperatures, such as 25 °C. Useful organic buffers and their properties are summarized below:

Organic base	pKa	Effective pH use range
Pyrrolidine	11.3	10.3–12.3
Triethylamine	10.7	9.7–11.7
1-Methyl-piperidine	10.3	9.3–11.3
Glycine	9.8	8.8–10.8
TRIS	8.1	7.1–9.1
Ammonia	9.2	8.2–10.2

All of these buffers can be made by titrating with HCl to the desired pH. Borate buffers apparently also can be used, but less experience is available with these materials. Whenever possible, carbonate and phosphate buffer should be avoided at both intermediate and high pH, since these materials enhance the solubility of the silica support to cause column failure, compared to the organic buffers listed above.

Studies indicate that methanol is a preferred organic modifier for high pH applications, since the rate of silica support dissolution and subsequent column degradation is slower with this solvent than with acetonitrile. Methanol also can produce better peak shapes when the column is used at intermediate pH. No preference for organic modifier was noted for low pH applications.

1-Methyl-piperidine as supplied by Aldrich Chemical Co., Inc. (M7,260-9) has been successfully used as a buffering material with the Extend-C18 column (for example, see *J. Chromatogr. A*, 797 (1998) 111). While this chemical is reputed to have 99% purity, prolonged use can result in some impurities building up on the inlet, changing column characteristics. Should this occur, a thorough purge of the column with 50% methanol/50% 0.05% trifluoroacetic acid generally restores the column to original performance. The operating pH of buffers used for separating basic compounds should be at least one pH unit (preferably 1.5 pH units) above the pKa of the basic compound of interest. (See J. J. Kirkland, *Current Issues in HPLC Technology, LC-GC Supplement*, May 1997, S46, Figure 13).

Best stability at pH >6 is obtained using organic buffers, and temperatures not exceeding 40 °C, as suggested in references on higher pH separations with silica-based columns [*J. Chromatogr. A* 797 (1998) 111; *J. Chromatogr. A* 762 (1997) 97; *J. Chromatogr. A* 728 (1996) 259; *J. Chromatogr. A* 691 (1995) 3]. Phosphate and carbonate buffers should be avoided for best stability of all silica-based columns at intermediate and high pH. Successful operation and good column lifetime at pH 11.5 and 40 °C with a buffer made from Aldrich pyrrolidine also has been demonstrated. Good column stability also has been found with a pH 10.5 mobile phase of ammonia/methanol, which is a useful medium for high pH separations when using a mass spectrometer as a detector. An overview of bidentate chemistry and chromatography was published in *Anal. Chem.* 70 (1998) 4344-4352.

Instrument Guidelines

Many HPLC instruments, including Agilent instruments use Vespel rotor seals in their injection valves (both manual and autosampler injection valves). Vespel is recommended for use up to pH 9.5. At pH's above 9.5, Vespel rotor seals will start to degrade, possibly causing plugging of downstream components in the flow path including the HPLC column. Therefore, Vespel rotor seals should be replaced with Tefzel rotor seals, which are stable up to pH 12.5, for applications using pH's above 9.5. The appropriate part numbers are listed below for Agilent instruments:

Agilent autosampler	Tefzel rotor seal p/n
Model 1090	1535-4900
Model 1050	0101-1849
Model 1100	0100-1849
Agilent manual injector	Tefzel rotor seal p/n
Model 1090 (Rheodyne 7010)	1535-4900
Model 1090 (Rheodyne 7125)	0101-0620
Model 1050 (Rheodyne 7125)	0101-0620
Model 1100 (Rheodyne 7125)	0101-0620

Operational Guidelines

- The direction of flow is marked on the column.
- These columns are packed and assembled for high pressure (up to 1200 bar) use. Disassembling the column will degrade column performance.
- Extend-C18 is compatible with water and all common organic solvents.
- Avoid use of this column below pH 2 or above pH 11.5; optimum operating range is pH 2 - 11.5 (see "pH Guidelines" section).
- Maximum operating pressure is 1200 bar (17000 psi).
- Maximum operating temperature is 60 °C.

NOTE: Extend-C18 columns are designed for stability over a wide pH range. When using these silica-based columns under conditions of pH > 6, maximum column lifetime is obtained by operation at low temperatures (< 40 °C) using low buffer concentrations in the range of 0.01 to 0.02 M. Column stability at pH > 6 is also enhanced by avoiding phosphate and carbonate buffers (see H. A. Claessens, M. A. van Straten, and J. J. Kirkland, *J. Chromatogr. (A)*, 728 (1996) 259).

Mobile Phase Selection

The bonded stationary phase is nonpolar in nature and is best used with mobile phases such as methanol/water or acetonitrile/water mixtures. Increasing the amount of organic component usually reduces the retention time of the sample. Due to the relatively high viscosity of recommended mobile phases, increased efficiency can be achieved with the use of column temperatures in the range of 40–60 °C; however, best column lifetime is achieved with operation at ≤ 40 °C. Gradient-elution techniques for this packing often use 5% methanol or acetonitrile as the initial solvent and 100% methanol or acetonitrile as the final solvent.

Column Care

Samples that contain particulate matter may plug the column inlet frit and should be filtered before injection into the column. The column inlet frit is nominally 0.5 µm and samples should be filtered through a 0.2-µm sample filter. To remove strongly retained materials from the column, flush the column with stronger (less polar) solvents. Solvents such as methanol, acetonitrile, or a 95/5 mixture of dichloromethane and methanol should remove most highly retained compounds. In extreme cases, dimethyl sulfoxide or dimethylformamide at low flow rates may also be used for this purpose. When switching between solvents with vastly different polarities, it may be necessary to first purge the column with a mutually miscible solvent such as isopropanol.

Since columns have 3/8-inch end nuts, a short 3/8-inch wrench should be used to attach the columns to the instrument to avoid any additional tightening of the end fittings. Over tightening the end fittings will cause damage and require column replacement

Storage Recommendations

Long term storage of silica-based, bonded phase columns should be in a pure organic solvent, preferably an aprotic liquid such as 100% acetonitrile. If the column was previously used with a buffered mobile phase, the buffer should first be removed by purging the column with 20–30 column volumes of a 50/50 mixture of methanol or acetonitrile and water, followed by 20–30 column volumes of the pure solvent. Before storing the column, the endfittings should be tightly capped with end-plugs to prevent the packing from drying out. Columns may be safely stored for short periods in most mobile phases. However, to protect equipment, it is desirable to remove salts from the instrument and column by purging the column with the same mobile phase without the buffer (for example, using 60/40 ACN/H₂O to remove a 60/40 ACN/0.02 M phosphate buffered mobile phase). Re-equilibration is rapid with the original mobile phase when using this approach, and any danger of corrosion from the salts is eliminated.

Agilent Ordering Information

For more information on our products, visit our Agilent Technologies home page on the World Wide Web at:
<http://www.agilent.com/chem/supplies>
For Technical support in the US and Canada, call 1-800-227-9770 or call your local Agilent sales office.

www.agilent.com/chem

Agilent shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.

Information, descriptions, and specifications in this publication are subject to change without notice.

© Agilent Technologies, Inc., 2010
Printed in the USA
December 2, 2010
Part No. 820214-002



Agilent Technologies