

SureSelect Human All Exon V6

Bridging the Gap



Key Benefits

Most comprehensive content, an exome solution for every application

- Up-to-date core content from relevant databases, targets more exons including hard-to-capture regions
- Easily add UTRs for translational research, COSMIC for cancer research, or custom content for specific applications

Deep and complete target coverage for definitive answers

- Highly optimized design for superior coverage and uniformity
- Increase breadth of coverage of targeted regions


Complete and flexible solutions from sample to data

- Compatible with SureSelect workflows for faster time to results
- Solutions from design, library prep, sample QC, automation and data analysis

Overview

Whole exome sequencing is a powerful technology used to analyze protein coding regions of the genome. This massively parallel analysis, coupled with the single molecule resolution of next-generation sequencing, has played a major role in accelerating disease gene discovery by providing a cost-effective, single reaction method for analysis of the exons that constitutes the 20,000 or so genes in the genome. The power of this technology extends to not only translational research, but also clinical research. For translational research, exome sequencing allows for the analysis of all protein coding regions to study the effect of DNA variants on normal biology and disease pathogenesis. In the clinical research setting, specifically in the study of rare disorders, the simultaneous analysis of these targets enable the facilitated detection of disease-associated variants out of the 20,000 to 50,000 that can be found in the human exome. Today, with the higher standards set for target coverage to increase confidence in variant detection, the challenge is to provide comprehensive designs that deliver complete target coverage.

Stringent Capture for High Performance and Deep Coverage of Targets

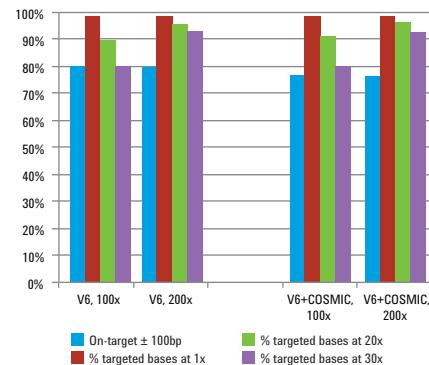

	SureSelect Human All Exon V6	SureSelect Human All Exon V6+COSMIC	Competitor I	Competitor N
Target Size	60Mb	66Mb	37Mb	64Mb
% DATABASES COVERED				
RefSeq	99%	99%	98%	99%
CCDS	99%	99%	99%	99%
GENCODE	99%	99%	98%	99%
HGMD_cds	99%	99%		
OMIM_cds	99%	99%		
COSMIC		96%		
Customization	YES	YES	NO	YES

Figure 1. SureSelect Human All Exon V6's highly optimized bait design coupled with stringent capture workflows enables a high on-target metric that results in a specific mapping of reads to targets for deep coverage. This exome targets updated content from relevant databases, including hard-to-capture regions, for comprehensive analysis of protein-coding regions.




Agilent Technologies

SureSelect Human All Exon V6

Deep and Complete Target Coverage, Comprehensive and Accurate Variant Calling

The SureSelect Human All Exon V6 is a high performing design that targets updated content from databases relevant to both the clinical and translational researchers. This content, which includes targets that in previous designs and other commercial designs have proven to be a challenge to capture, enables comprehensive coverage of protein coding regions (Figure 1).

In addition, with optimized bait design and boosting strategies that improve on coverage uniformity, the SureSelect Human All Exon V6 is able to provide high 20x end-to-end coverage of targeted exonic regions (Figure 2). To demonstrate, the coverage of genes important in constitutional research are shown. This improvement helps address the current observation of the inverse correlation between depth of coverage and breadth of coverage providing balanced coverage, maximizing variant information output for per exome sequenced.

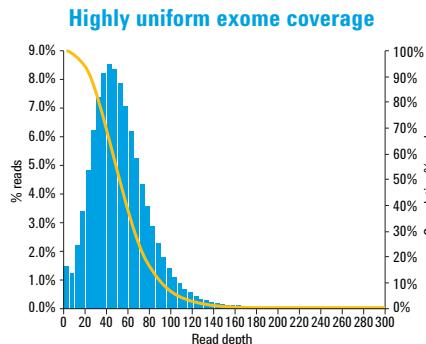

Deep coverage of 56 key regions in constitutional research for complete and accurate variant calling

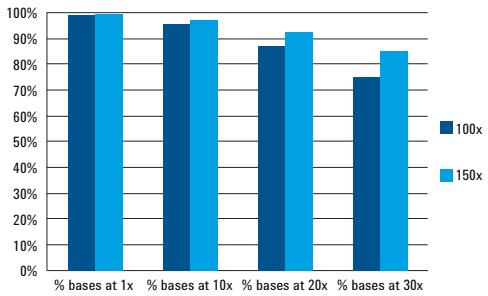


Figure 2. Optimized bait selection and boosting enables highly uniform coverage of the exome and each individual targeted exon as well. With this optimized design, targeted regions demonstrate excellent coverage which greatly facilitates variant identification. To demonstrate, coverage of 56 genes important in constitutional research is shown.

From Translational to Clinical Research, a Complete Exome Solution Tailored for Every Application

The utility of exome sequencing is evaluated as a balance between relevant information gained and cost of sequencing. In this respect, content definition becomes an application-specific question. To address the needs of the researchers in the constitutional space, typical reference databases used to curate and annotate new genes and transcripts were mined for SureSelect Human All Exon V6 ± UTR targets to facilitate detection of disease-associated variants. For cancer researchers on the other hand, the SureSelect Human All Exon V6 also provides an option that includes targets from both TCGA and COSMIC. These databases catalog variants that play roles in the initiation and progression of cancer, providing a better understanding of disease pathogenesis. Addition of custom content using SureDesign further enables this exome to be tailored to specific applications.

Accelerated workflows are critical to providing reduced turn-around time from sample to data. The SureSelect Human All Exon V6 is compatible with SureSelect workflows using the shortest hybridization times, as little as 90 minutes, providing faster sample to sequencing.

Data analysis using SureCall enables simplified workflows for single sample or trio analysis, suited for constitutional research or analysis of paired tumor and normal samples, typical in cancer research, addressing another critical bottleneck in translational or clinical research (Figure 3).

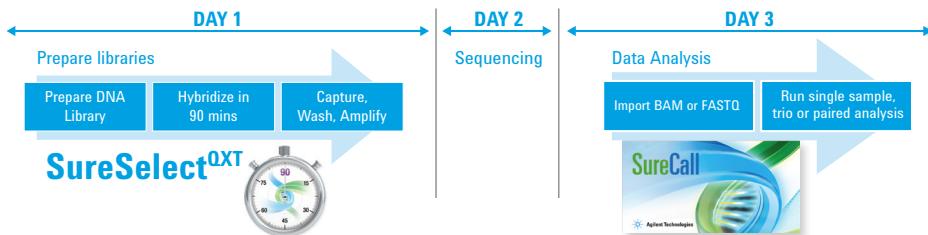


Figure 3. Faster workflows from sample to data are enabled by using SureSelect^{QXT}, the only single-day capture solution, coupled with SureCall, a guided analysis software for single sample, paired, or trio analysis.

Ordering Information:

	16rxn	96rxn	96rxn auto
SureSelect ^{XT} Human All Exon V6	5190-8863	5190-8864	5190-8865
SureSelect ^{XT} Human All Exon V6 +UTR	5190-8881	5190-8882	5190-8883
SureSelect ^{XT} Human All Exon V6 +COSMIC	5190-9307	5190-9308	5190-9309
SureSelect ^{XT} Human All Exon V6 Plus 1	5190-8866	5190-8867	5190-8868
SureSelect ^{XT} Human All Exon V6 Plus 2	5190-8869	5190-8870	5190-8871
SureSelect ^{XT2} Human All Exon V6	5190-8872	5190-8873	5190-8874
SureSelect ^{XT2} Human All Exon V6 +UTR	5190-8884	5190-8885	5190-9306
SureSelect ^{XT2} Human All Exon V6 +COSMIC	5190-9310	5190-9311	5190-9312
SureSelect ^{XT2} Human All Exon V6 Plus 1	5190-8875	5190-8876	5190-8877
SureSelect ^{XT2} Human All Exon V6 Plus 2	5190-8878	5190-8879	5190-8880

Find an Agilent customer center in your country:

www.agilent.com/genomics/V6

U.S. and Canada: 1-800-227-9770

cag_sales-na@agilent.com

For Research Use Only. Not for Use in Diagnostic Procedures.

PR7000-0152

© Agilent Technologies, Inc. 2015

Printed in USA, May 14, 2015

5991-5572EN