



Agilent CytoGenomics Software

Learn how to analyze chromosomal microarray data from blood, single cell, hematological and solid tumor samples

Anniek De Witte

Sr. Product Manager, CGH and NGS Software

Oct 1st 2013



Agilent Technologies

Agenda

1. Introduction

2. Live analysis of microarray data in CytoGenomics software:

- Blood
- Hematological tumor sample
- FFPE tumor sample
- Single cell

3. Q&A


What is CytoGenomics?

INTRODUCTION


Agilent CytoGenomics Software

Analyze, monitor, review all your samples in a few steps

- Workflow-based analysis tool for CGH and CGH+SNP microarray analysis
- Powerful algorithms: Accurate copy-number & LOH calls
- Designed specifically for cytogenetic research
- Available at no charge to all Agilent microarray customers
www.agilent.com/cytogenomics

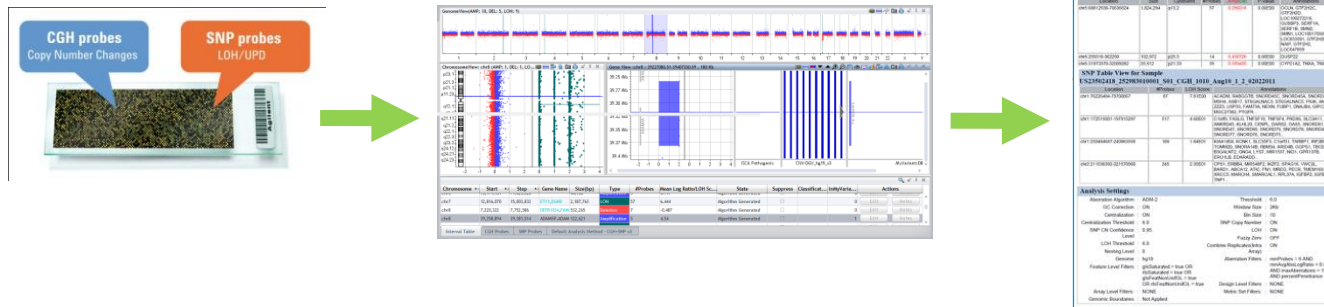


CGH+SNP wet lab and analysis workflow

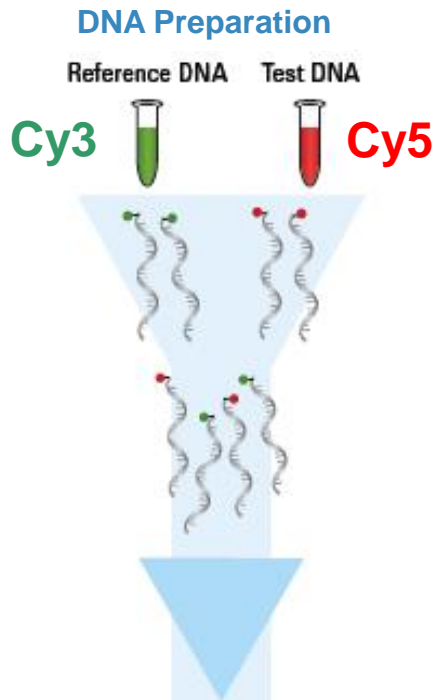


SNP data:
LOH/UPD


CGH data:
Copy number
changes

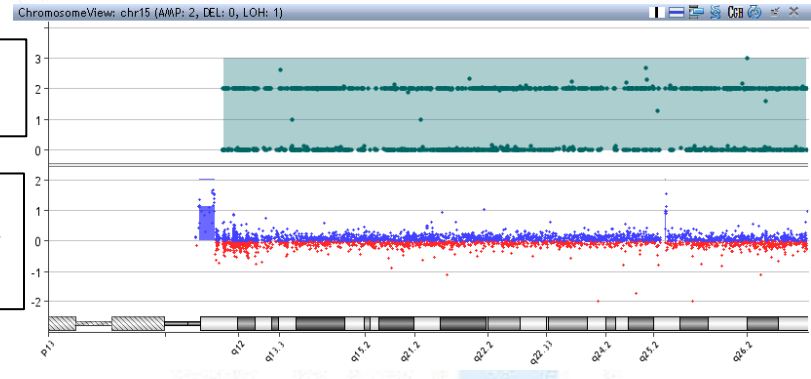

Hybridization

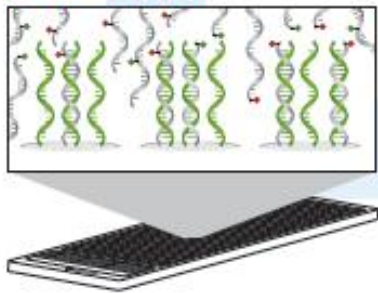


Data Analysis

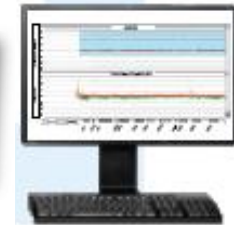




Scanning

Software

Schematic view of CGH data output

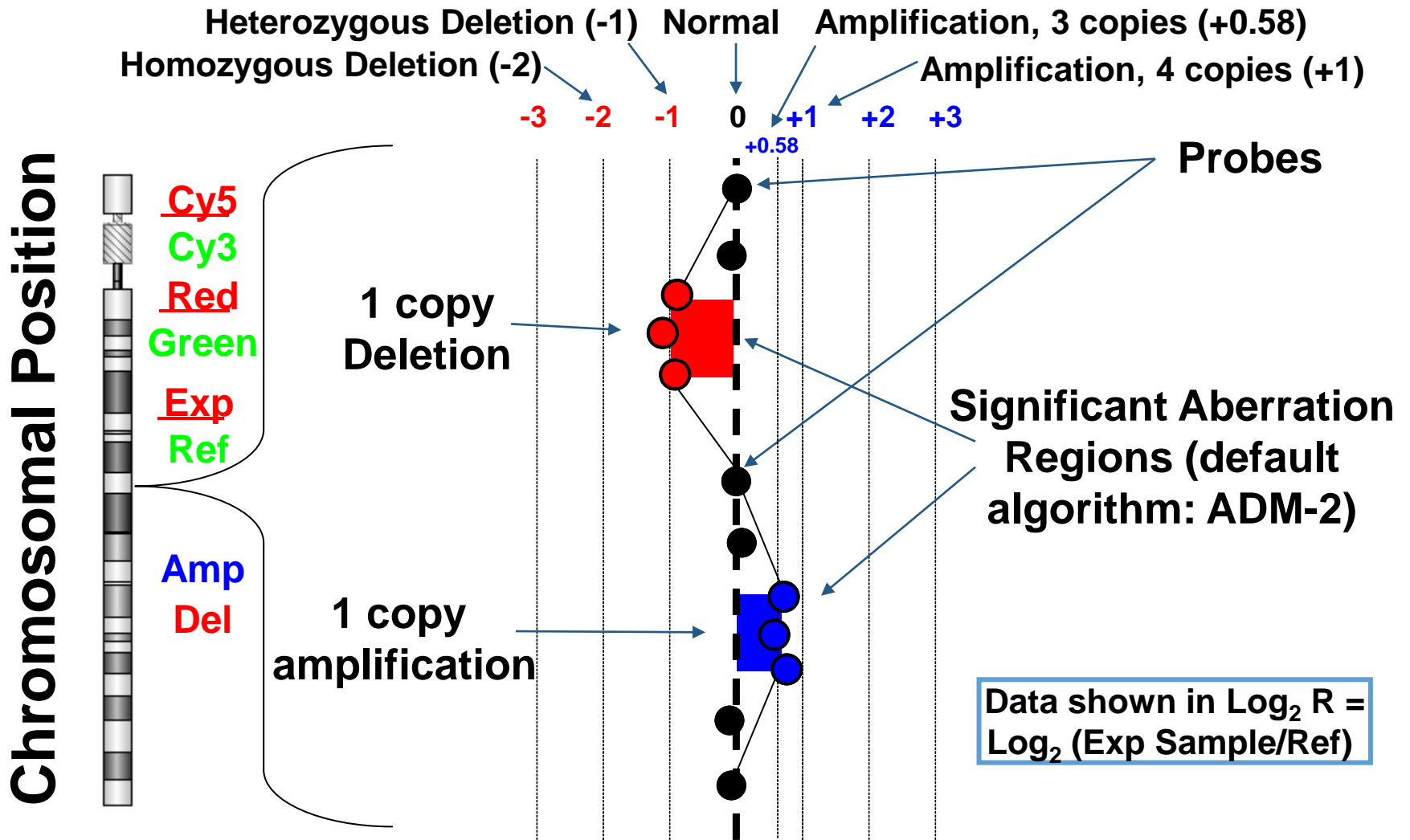


Mechanism of Agilent SNP Probes

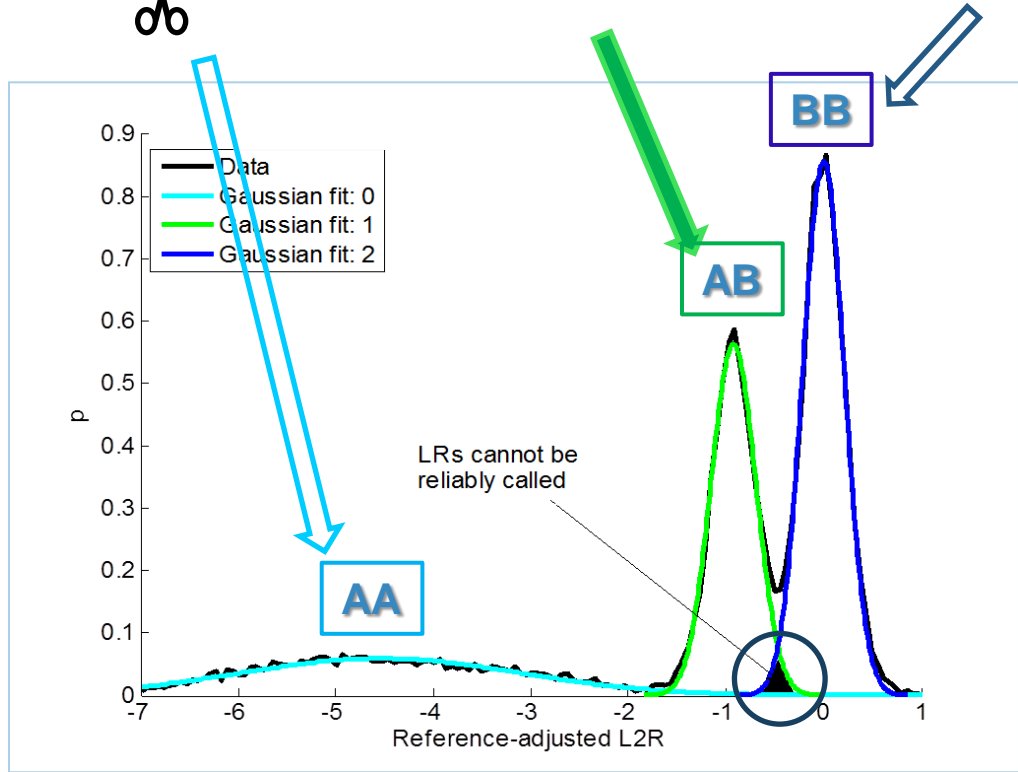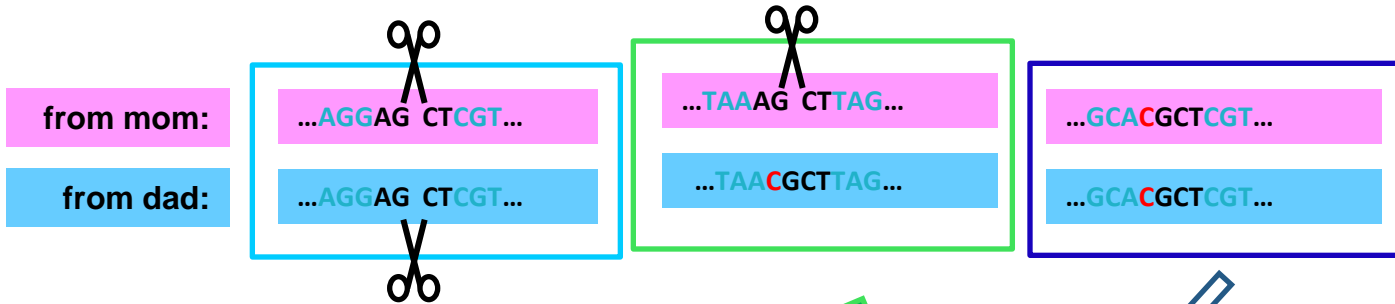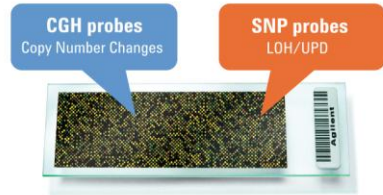


- Signal measures copies of uncut allele
- Raw SNP data: black curve
- Colored Gaussians: fits to distributions

Schematic view of SNP data output

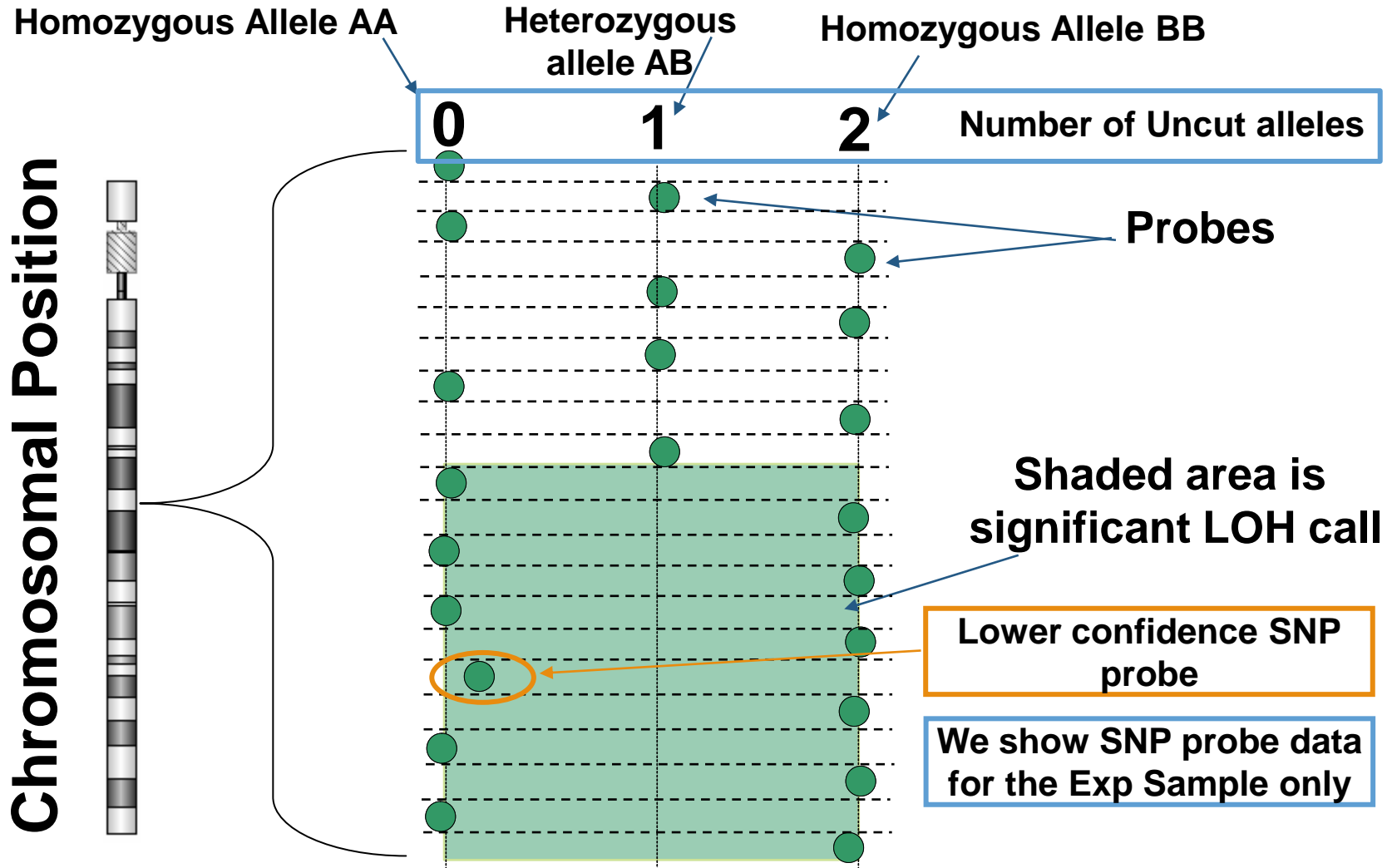

Clinical research on different sample types using SurePrint CGH microarrays and CytoGenomics

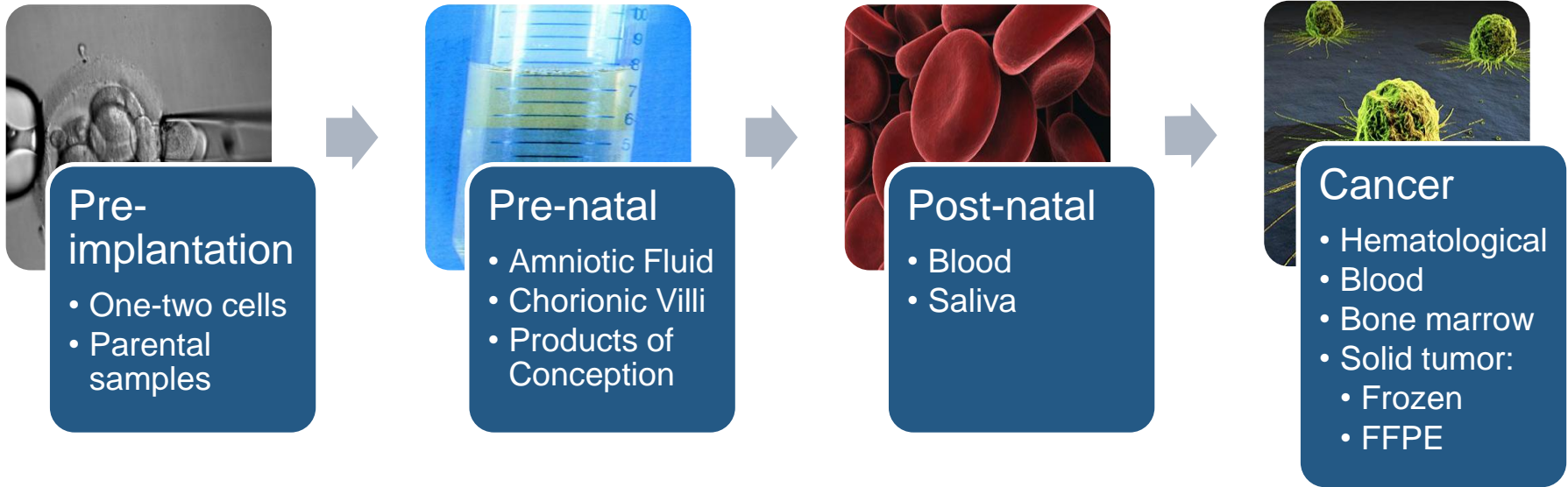

CYTOGENOMICS LIVE DEMO

CGH math: We measure the difference between 1 sample and 1 reference DNA, usually expressed as \log_2 (Sample/reference). Actual data are slightly compressed from ideal values.

| Description | Average Sample CN | Ref CN | Ratio (S/R) | Ideal \log_2 (Ratio) | Actual data |
|---------------------|-------------------|--------|-------------|------------------------|-------------|
| Diploid | 2 | 2 | 1 | 0 | 0 |
| Deletion | 1 | 2 | 0.5 | -1 | -0.9 |
| Trisomy | 3 | 2 | 1.5 | +0.58 | +0.53 |
| 50% mosaic deletion | 1.5 | 2 | 0.75 | -0.41 | -0.37 |
| 50% mosaic trisomy | 2.5 | 2 | 1.25 | +0.32 | +0.29 |
| 20% mosaic deletion | 1.8 | 2 | 0.9 | -0.15 | -0.13 |

Questions

