




## Overview


This instruction sheet could help you keep your firmware update utility current without downloading the entire 35 Megabyte package for version A.01.04. The only difference between versions A.01.02 and A.01.04 are the firmware files. These are stored in a set of subfolders. Each subfolder is associated with a different GC or Auto-sampler module. This instruction assumes that you have downloaded one or all of the updated firmware. Please read the entire sheet before beginning.

## Procedure


1. Launch Windows Explorer. Find the file that you downloaded from the internet.
2. Double-click on the download.exe file and save the extraction to a location on your PC. The default location is C:\extract.
3. Navigate to the extracted files. This picture shows the 5 files from the August 12<sup>th</sup> list.

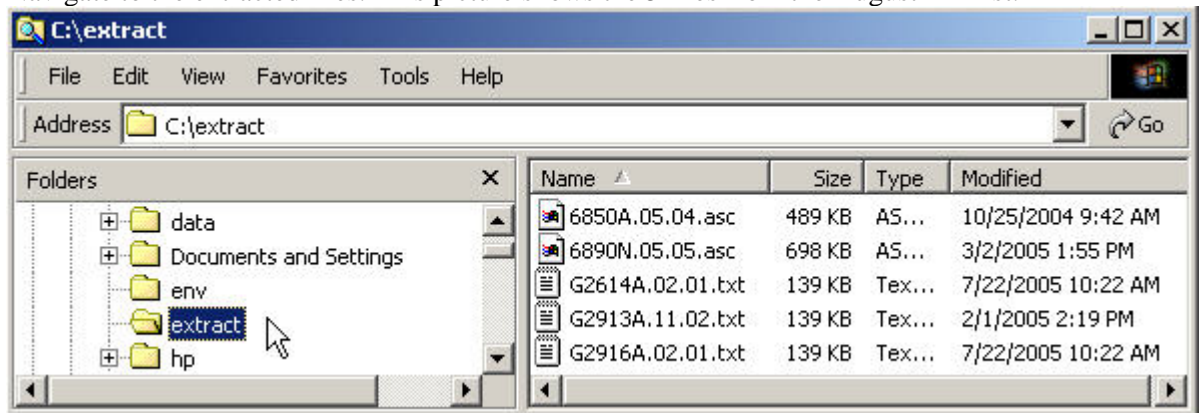

4. Navigate to your main drive and expand the Program Files folder by clicking on the + to the left of the folder. Expand the Agilent subfolder. Expand the Agilent GC Firmware Update Utility folder. Expand the Firmware folder.

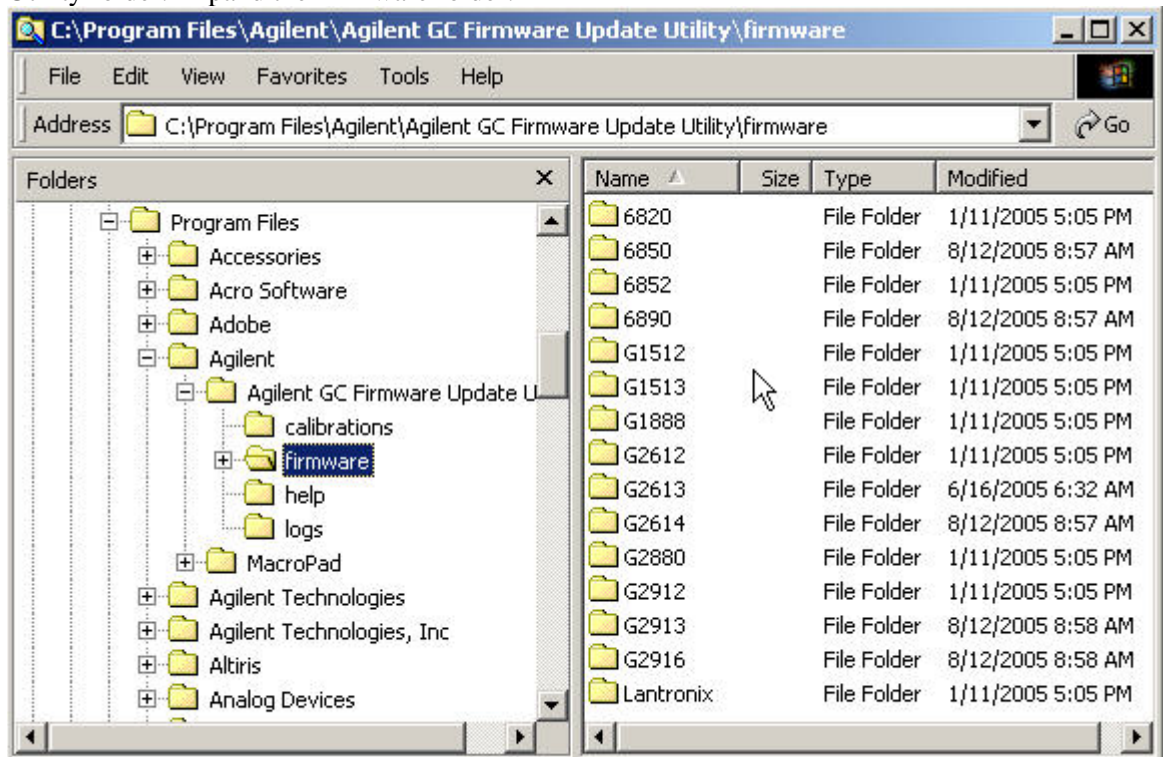

- Drag and drop each file into the appropriate firmware subfolder. For example 6890N.05.05.asc goes into the 6890 subfolder. Another example, the G2913A.11.02.txt goes into the G2913 subfolder.

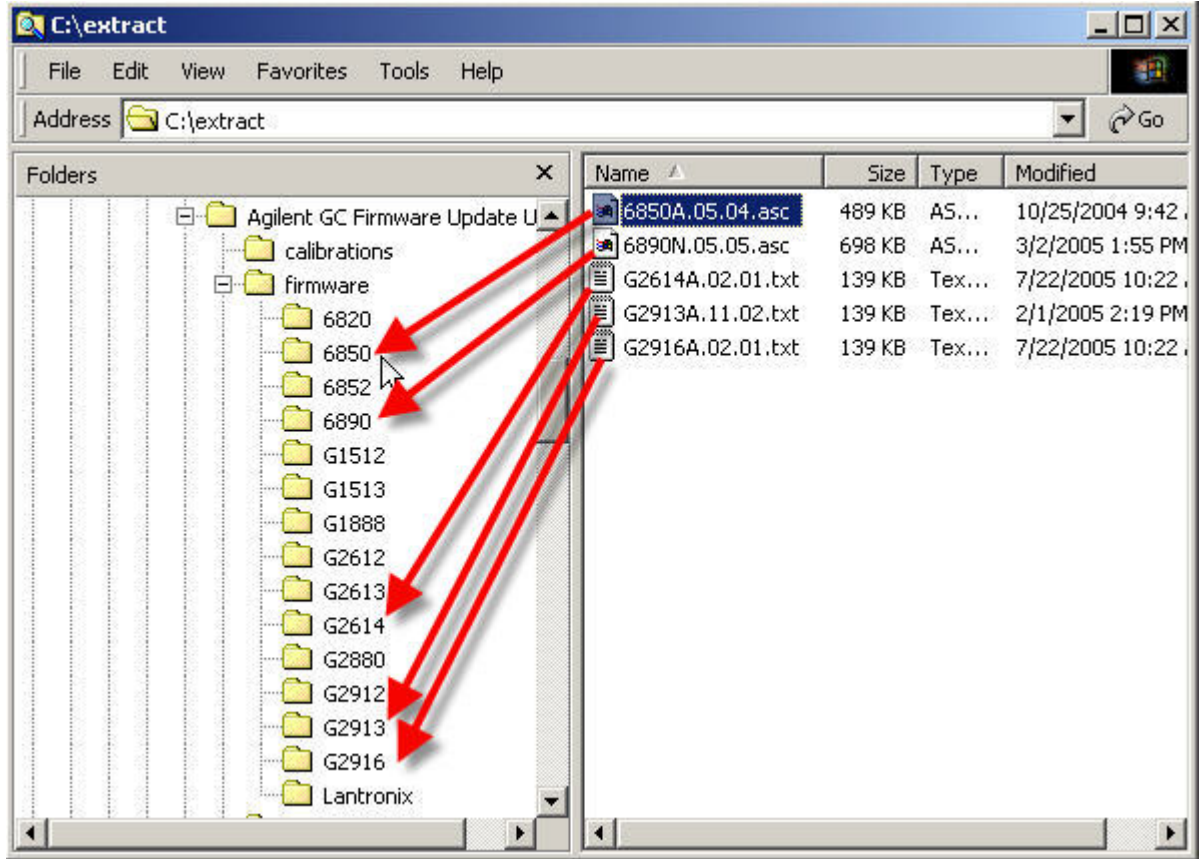

- Because these are read-only files, your PC may display a confirmation message. Click Yes to confirm and continue.



NOTE: The software will not recognize the new file if you put it in the wrong subfolder. If you attempt to update a module and the new firmware revision does not appear in the drop down list, you can either download the file again or move the file into the correct subfolder.

