

## Spraychamber Stand Instructions

The spraychamber stand is supplied as a flat plate. It is perforated for ease of folding and can therefore be readily folded manually – no tools are required and no special skill is needed.

### To assemble the stand:


- 1 Refer to Figures 1 and 2.

Lay the 'unformed' spraychamber stand on a suitable bench. Grasp one outer panel firmly as shown in Figure 1, and fold the panel upward so that it is perpendicular as shown in Figure 2.

- 2 Refer to Figures 2 and 3.


Grasp the opposite panel firmly as shown in Figure 2, and fold the panel upward so that it is perpendicular as shown in Figure 3.

- 3 Place the completed stand on the bench as shown in Figure 4. You may now use the stand to accommodate your complete spraychamber assembly including the liquid trap.

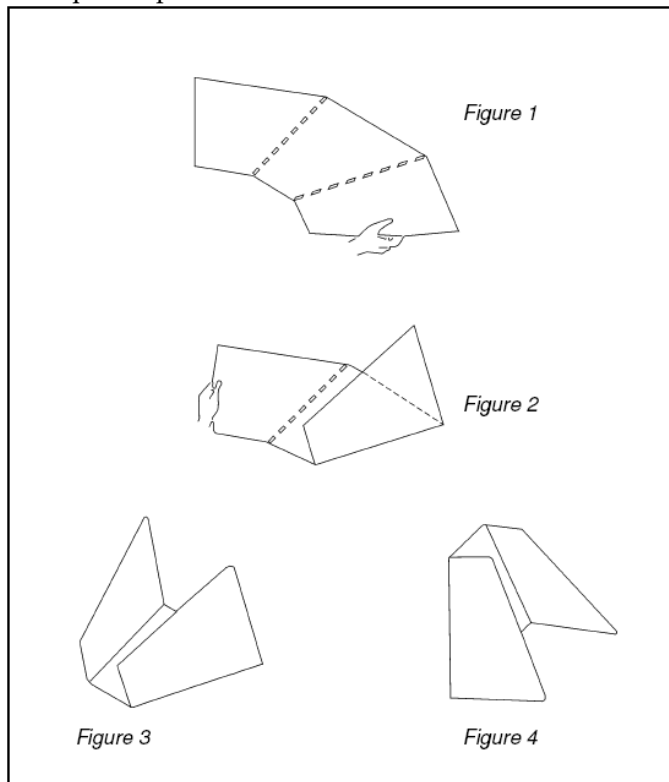

## Spraychamber Stand Instructions

This information is subject to change without notice.

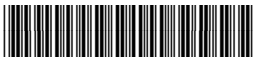

8510065700

Part Number: 8510065700

Edition 10/12

Issue 3

© **Agilent Technologies, Inc. 1999, 2010, 2012**

Agilent Technologies  
679 Springvale Road  
Mulgrave, VIC 3170