



Cary 50 Accessory Cable Kit

Installation category I
Pollution degree 2
Equipment class III

Introduction

The Accessory Cable kit allows you to use accessories such as the Temperature Probe on the Cary 50.

Unpacking

When you receive your Accessory Cable kit, unpack it and check that you have received all of these necessary parts:

- Accessory cable with bracket
- Accessory connector (ribbon cable attached)
- Screw (4 x 8 mm)

CAUTION

Inspect all parts for damage in transit. Any damage should be reported immediately. Refer to the accessory by model and part number.

Safety information

Before using this accessory, you must read the Safety Practices and Hazards section in your Cary operating instructions.

If the equipment is used in a manner not specified by the manufacturer, the protection provided by the equipment may be impaired.



Specifications

For environmental and power supply requirements, refer to your Cary operating instructions.

Connections

Computer internal: 40-way flat ribbon cable with 37-pin D-range connector (plugs into Cary 50 PC card)

Cary-computer accessory cable: 37-pin D-range connector from computer and 25-pin D-range connector mounted inside sample compartment

Weights

Packed: 1.0 kg (2.2 lb)
Unpacked: 0.82 kg (1.8 lb)

This information is subject to change without notice.



8510157200

Part Number: 8510157200

Edition 04/12

Issue 5

© Agilent Technologies, Inc. 1998, 1999,
2011 and 2012

Agilent Technologies
679 Springvale Road
Mulgrave, VIC 3170