



## Cary 50 Accessory Cable Kit

Installation category I  
Pollution degree 2  
Equipment class III

### Introduction

The Accessory Cable kit allows you to use accessories such as the Temperature Probe on the Cary 50.

### Unpacking

When you receive your Accessory Cable kit, unpack it and check that you have received all of these necessary parts:

- Accessory cable with bracket
- Accessory connector (ribbon cable attached)
- Screw (4 x 8 mm)

#### CAUTION

Inspect all parts for damage in transit. Any damage should be reported immediately. Refer to the accessory by model and part number.

---

### Safety information

Before using this accessory, you must read the Safety Practices and Hazards section in your Cary operating instructions.

If the equipment is used in a manner not specified by the manufacturer, the protection provided by the equipment may be impaired.



## Specifications

For environmental and power supply requirements, refer to your Cary operating instructions.

### Connections

Computer internal: 40-way flat ribbon cable with 37-pin D-range connector (plugs into Cary 50 PC card)

Cary-computer accessory cable: 37-pin D-range connector from computer and 25-pin D-range connector mounted inside sample compartment

### Weights

Packed: 1.0 kg (2.2 lb)  
Unpacked: 0.82 kg (1.8 lb)

This information is subject to change without notice.



8510157200

Part Number: 8510157200

Edition 04/12

Issue 5

© Agilent Technologies, Inc. 1998, 1999,  
2011 and 2012

Agilent Technologies  
679 Springvale Road  
Mulgrave, VIC 3170