



## Cary Temperature Probe Holder Instructions


Enclosed are probe holders for use with your Multicell Holder.

### CAUTION

Temperature probes are precision devices. To prevent damage, take care when inserting or removing probes and do not flex the probe cable more than necessary.

#### To install the probe holder on the temperature probe:

- 1 Separate the holder by unscrewing the two sections. A small o-ring will be found inside.
- 2 Push the upper section of the holder onto one of the temperature probes.
- 3 Push the black o-ring onto the probe.
- 4 Push the cell cap onto the probe and screw the two sections of the holder together.
- 5 You should now fit the probe and cell cap into the top of your cuvette in the Multicell Holder.

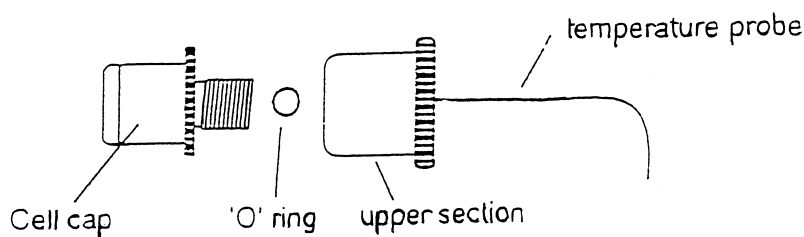

**Figure 1.** Assembly of the temperature probe holders



This information is subject to change without notice.



8510157900

Part Number: 8510157900

Edition 07/12

Issue 3

© Agilent Technologies, Inc. 2001, 2011,  
2012

Agilent Technologies  
679 Springvale Road  
Mulgrave, VIC 3170