

Peltier 4 Position Multicell Holder Accessory

Installation category I

Pollution degree 2

Equipment class III

Safety information

Before using this accessory, you must read the Safety Practices and Hazards section in your Agilent Cary Eclipse User's Guide.

A link to the Safety section can also be found on the main page of the Agilent Cary Eclipse Help.

If the equipment is used in a manner not specified by the manufacturer, the protection provided by the equipment may be impaired.

Introduction

The Peltier-thermostatted 4 Position Multicell Holder for the Agilent Cary Eclipse is an accessory that permits fluorescence measurements on up to four samples, while precisely controlling the temperature and rate of stirring.

Unpacking notes

Make sure that all of the adhesive tape, plastic and cardboard is removed before operating the Peltier Multicell Holder.

Installation

- 1 Turn the Agilent Cary Eclipse instrument power off using the switch on the front right side of the instrument. You may leave the PC turned on.
- 2 Slide the top panel of the sample compartment all the way back and completely remove the front panel by sliding it upwards. Ensure that the sample compartment is empty.
- 3 Lift the Multicell Holder by the handle and hold the end of the hose in your other hand. Rest the accessory at the front of the sample compartment in order to allow access to the D-shaped, 25 pin port at the left side of the rear sample compartment wall. Connect the cable plug to the 25 pin port.
- 4 Ensure that the fastening lever is in the open position (i.e., to the left). Hold the extended nose (the black metal flange which contains the stirring controls, see Figure 1) and align the two holes on the base of the accessory with the locating pins on the floor at the rear of the sample compartment. Move the fastening lever from left to right to lock the accessory into position.



Agilent Technologies

Peltier 4 Position Multicell Holder Accessory

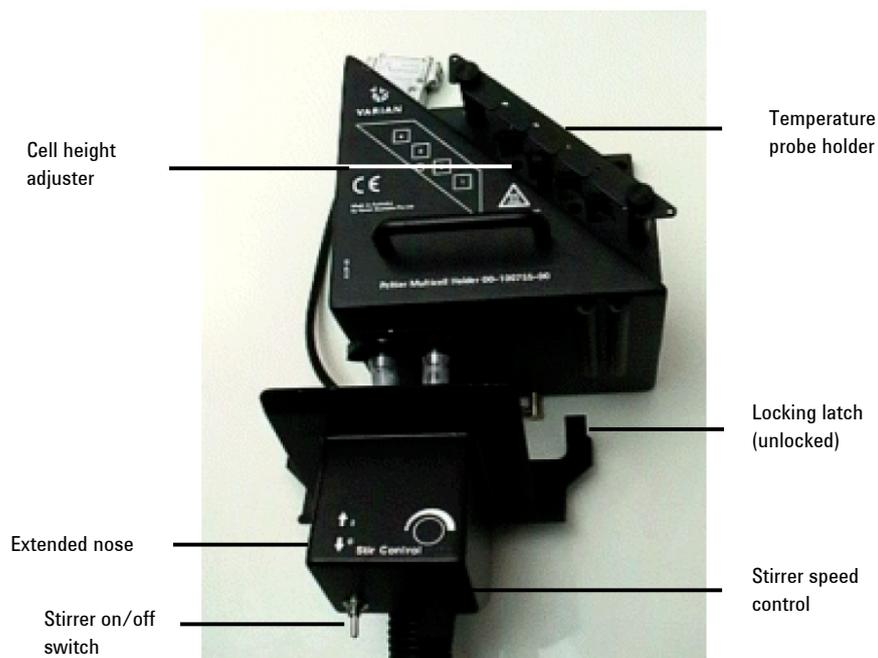


Figure 1. Peltier 4 Position Multicell Holder accessory

- 5 Ensure the two black locking latches at the back of the extended nose are pointing upwards (unlocked, see Figure 1). Position the extended nose into place by aligning its locating pins with the holes in the front of the sample compartment floor. Lock the extended nose into position by pushing the locking latches downwards.
- 6 Slide the special door for thermostatted accessories into position ensuring the electrical cable is tucked neatly into the sample compartment.
- 7 Connect the hose ends to the Temperature Controller inlet and outlet connections.

NOTE

The Peltier Multicell Holder must only be used with an Agilent Temperature Controller.

- 8 Connect the circular communications plug to the 14-pin socket on the Temperature Controller.

NOTE

You may need to rotate the communications plug before it will slot into the communications socket, as it has been designed to be inserted in one position only.

- 9 Ensure the Temperature Controller has sufficient liquid for optimal operation.

NOTE

The temperature probe holder (see Figure 1) that spans the two black posts on the multicell holder should be removed when inserting or removing cuvettes.

Operation

Operation of the accessory is through the Agilent Cary Eclipse software, with the exception of the stirrer control which is located on the front of the extended nose.

To use the multicell holder

- 1 From the 'Setup' dialog, select the 'Accessories' tab.
- 2 Select 'Multicell holder' and from the cell diagram, select the number of cells to use.
- 3 To enable temperature control (optional):
 - a Select **Temperature Control**. From the 'Device' option, select "Temperature controller".
 - b Enter the required temperature and select the monitor. This monitor will allow data collection to begin when the temperature of the monitoring device is within ± 0.5 °C.
- 4 To display the temperature of other items in the 'Status Display' window, select these items from the 'Temperature display' option. Figure 2 (below) shows a typical setup of the multicell peltier accessory.
- 5 Select **OK** to exit the 'Setup' window.
- 6 Select **Start** to begin data collection.

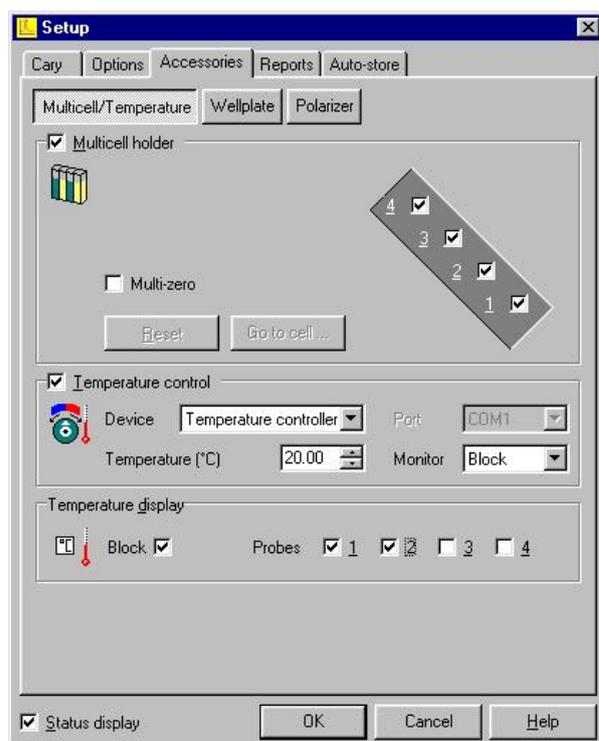


Figure 2. Accessories tab showing a typical setup of the multicell holder accessory

Specifications

For environmental and power supply requirements, refer to your Cary operating instructions. This product is suitable for indoor use only.

Connections

25-pin D-range connector on flying lead

14-pin circular lead (connect to Agilent Temperature Controller)

Two water circulation tubes (connect to Agilent Temperature Controller)

Weight

Packed	10.4 kg (22.9 lb)
Unpacked	4.2 kg (9.3 lb)

Dimensions

Packed	765 x 360 x 325 mm (30.1 x 14.2 x 12.8 in)
Unpacked	180 x 320 x 180 mm (7.1 x 12.6 x 7.1 in)

Maintenance and cleaning

All cleaning should be done with a soft cloth. If necessary, this cloth can be dampened with water or a mild detergent to clean black painted surfaces. Do not use organic solvents or abrasive cleaning agents.

This information is subject to change without notice.



8510183100

Part Number: 8510183100

Edition 04/12

Issue 5

© Agilent Technologies, Inc. 2000, 2011
and 2012

Agilent Technologies
679 Springvale Road
Mulgrave, VIC 3170