# Agilent Cary UV Workstation

## Software Registration Instructions

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### Introduction

SubscribeNet allows users to manage software entitlements, order media or download software and we recommend users renew subscriptions annually to have full access to the newest updates.

After installation of Cary UV Workstation, you will have 60 days to complete software registration as described in these instructions. To connect to an instrument after the 60-day period, you must register your software.

#### You will need the following:

- Activation code(s) – found in the lavender envelope(s) shipped with the software
- An internet connection to register your software
  - Is your instrument computer not connected to the internet? Perform Steps 4-19 on another computer that can access the internet.
  - USB drive – copy the license file generated at the end of Step 19 onto the USB and then onto the computer the Cary UV Workstation software is installed on
- SubscribeNet login credentials if your company has a pre-existing account
  - If you aren't sure, first check with your IT department or email [subscribenet\\_support@agilent.com](mailto:subscribenet_support@agilent.com)

### Registration

The main steps to register your software are:


- 1 Get the computer name and MAC address for the computer the Cary UV Workstation software was installed on from the OpenLab Control Panel application.
- 2 Create a SubscribeNet account or login with existing SubscribeNet username and password.
- 3 Enter any additional activation codes, if required.
- 4 Create the License file.

5 Import the License file into OpenLab Control Panel.

To find the computer name and MAC address:

- 1 Double-click the OpenLab **Control Panel** icon on the computer desktop or click the Windows **Start** button > **Agilent Technologies** > **Control Panel**.
- 2 Click **Licenses** from the **Administration** tab.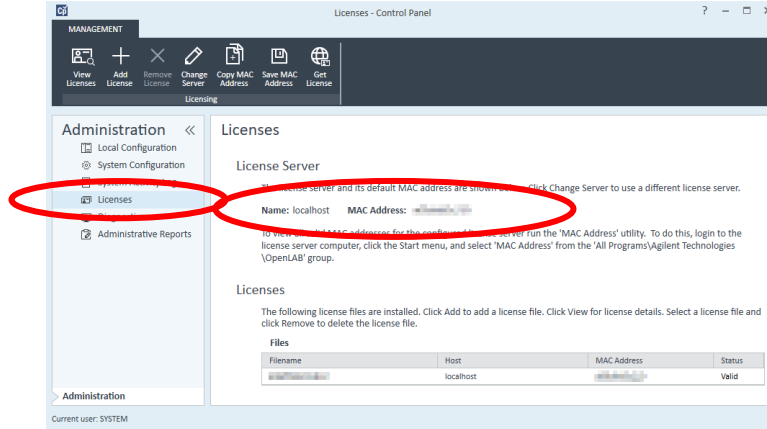

3 Take note of the server **name** and the **MAC Address**. You will need them in Step 15.

To create a **SubscribeNet** account or login with existing **SubscribeNet** credentials:

4 Press CTRL on your keyboard and then click <https://agilent.subscribe.net.com>.

**NOTE**

If your computer is not connected to the internet, print these instructions out and take them, and the USB, to a computer that will be used to continue your software registration.

5 If you already have a **SubscribeNet** login and password from previously registered software, enter in that information, click **Login** and then skip to **Step 11**. Otherwise, go to **Step 6**.

**NOTE**

If you are generating additional **Academic Licenses** from an already registered **Academic License** activation code skip to **Step 14**.

6 Click '**CLICK HERE**' to create a Login ID for a new account.

**Electronic Software and License Delivery**

Please login. Your Login ID is your Email address.

**Login ID**

**Password**

Remember my password until I logout

Login

If you have forgotten your login ID, password, or are not sure whether you have an account use our [Password Finder](#).

**SubscribeNet new account registration.**

Customers who have an authorization code from their Agilent product purchase may [CLICK HERE](#) to register and create a new **SubscribeNet** Account and Login ID.

7 Enter in the required details. The **Authorization Code(s)** are printed on stickers in the lavender envelope(s) provided in the software box.

**Electronic Software and License Delivery**

**Registration Information**

Registration information is used to create a new SubscribeNet Account and Login ID (email address) for the Account. An authorization code you received from your product purchase is required.

**Important Email Notes**

- Please ensure the email address is entered correctly or you will not receive a welcome email with your Login ID and Password.
- If the email entered has already been used for an account Login ID, you will not receive a welcome email. The new account will display as an additional selection after you log in to SubscribeNet.

\* Indicates a required field.

Authorization Code\*

Email\*

Re-enter Email to Verify\*

Company\*

Department\*

First Name\*

Last Name\*

Job Title

Phone\*

Address 1\*

Address 2

Address 3

City\*

State/Province\*

Postal Code

Country\*

- 8 Press **Submit**. You will receive an email with your username and password. If you do not receive an email within 15 minutes first check your email Spam or Junk folder. Otherwise, contact [subscribenet\\_support@agilent.com](mailto:subscribenet_support@agilent.com).
- 9 Open the email and click the SubscribeNet link. Enter your Login ID (the email address you registered with) and password and log in.
- 10 Follow the prompts to change the password.

To enter any additional activation codes, if required:

**NOTE**

When registering Academic License authorization codes, perform the following steps in the registration process once on *each* computer the Cary UV Workstation software is installed on. You can use the same SubscribeNet credentials for each registration, but the MAC address found in Step 2 and the license file generated in Step 14 must be specific to each computer.

- 11 If you used an existing username and password to log into SubscribeNet, click **Register Authorization Code**. Otherwise, go to Step 14.

Agilent Technologies | Our measure is your success. | Helping our customers achieve better results, faster than ever.

Your Login is: \_\_\_\_\_ Account: \_\_\_\_\_

**Electronic Software and License Delivery**

**Register Authorization Code**

Authorization Codes represent license or subscription products that can be added to your Account Lic using this page.


You may enter up to ten Authorization Codes, separated by semicolons.

Please enter an Authorization Code:


- 12 Enter in the Authorization Code included in the lavender envelope provided in the Cary UV Workstation software box.
- 13 Press **Submit**.

To create the license file:

- 14 Click **Generate or View Licenses**.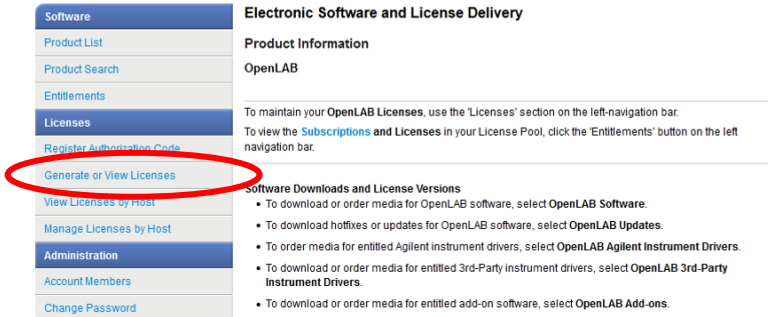

15 Select the tick box of the license to generate. Click **Generate**.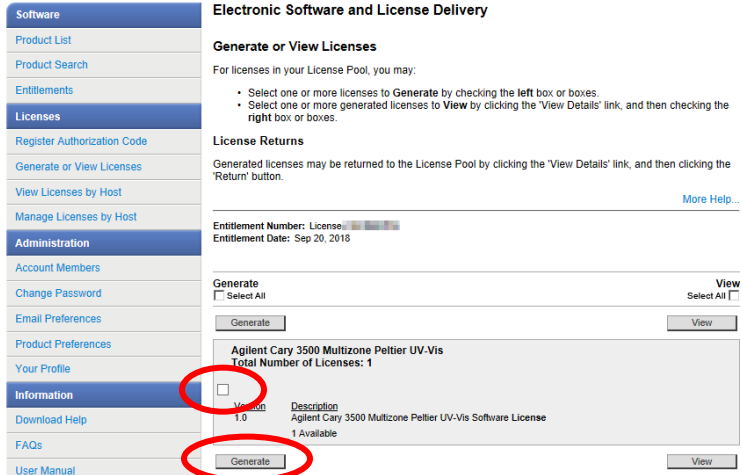


16 On the SubscribeNet window, enter the MAC Address and Name you noted in Step 3. Enter the **MAC Address** into the **HostID** field and the **Name** into the **Hostname** field. Enter a Description of your computer.

17 Leave the 'Number of Licenses to Generate' box set as the default value of 1.


**NOTE**

When registering Academic License authorization codes, enter 1 in the 'Number of Licenses to Generate' box.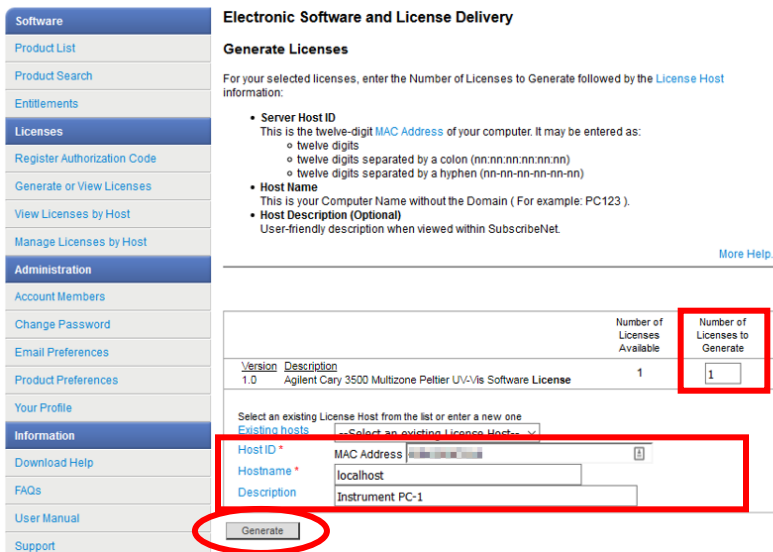

18 Click **Generate** to generate the license file.

- Click **Download License File**. A dialog box may appear. Follow the prompts to save the file or note where the file was automatically saved to, typically in your C:\Users\Username\Downloads folder.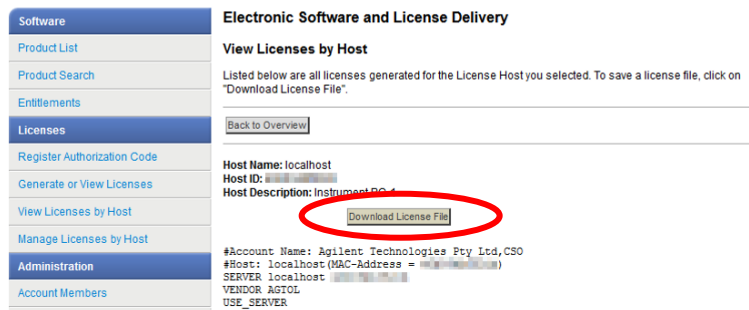


**NOTE**


If you are performing this registration on a computer other than the one the Cary UV Workstation software was installed on, navigate to the saved license file and then save it to a USB drive. Insert the USB drive into the computer that the Cary UV Workstation software is installed on before proceeding.

**To import the License file into OpenLab Control Panel:**

- In the OpenLab Control Panel, click **Add License**.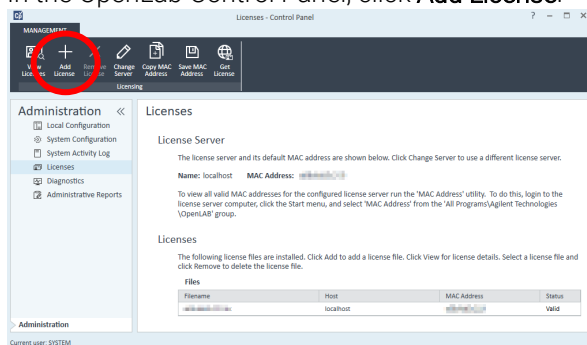

- Click **Select a license file** and then click **Browse**.

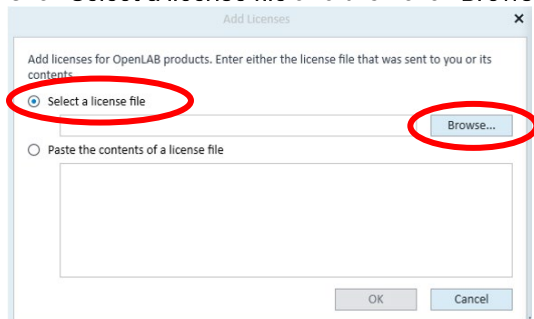- Navigate to and then select your previously saved license file on the computer or USB. Click **Open**.
- Select **OK** to Add your license.
- Your license will now appear in the list of Licenses in the OpenLab Control Panel and in SubscribeNet when you click 'Generate or View Licenses'.
- Restart the computer to complete the process.

**NOTE**

When registering Academic License authorization codes, repeat Steps 2-5 and then Steps 14-25 on each computer the Cary UV Workstation software is installed on. You will need to log in to SubscribeNet using the same credentials as the first registration before continuing with Step 14.

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© Agilent Technologies, Inc. 2018
Agilent Technologies Australia [M] Pty Ltd