

Agilent CrossLab Start Up Services

Agilent 7650 ALS Site Preparation Checklist

Thank you for purchasing an instrument from **Agilent Technologies**. CrossLab Start Up is focused on helping customers shorten the time it takes to start realizing the full value of their instrument investment.

Correct site preparation is the key first step in ensuring that your instruments and software systems operate reliably over an extended lifetime. This document is an **information guide and checklist** prepared for you that outlines the supplies, space, and utility requirements for the system set up in your lab.

Introduction

Customer Information

- If you have questions or problems in providing anything described as part of *Customer Responsibilities* below, please contact your local Agilent or partner support / service organization for assistance prior to delivery. In addition, Agilent and/or its partners reserve the right to reschedule the installation dependent upon the readiness of your laboratory.
- Should your site not be ready for whatever reasons, please contact Agilent as soon as possible to re-schedule any services that have been purchased.
- Other optional services such as additional training, operational qualification (OQ) and consultation for user-specific applications may also be provided at the time of installation when ordered with the system but should be contracted separately.
- Please refer to the other peripheral products (ie, sampling devices, etc.) for site preparation requirements.

Customer Responsibilities

Ensure that your site meets the following specifications before the installation date. For details, see specific sections within this checklist, including:

- The necessary laboratory or bench space is available.
- The required **environmental conditions for the lab** as well as laboratory gases, tubing.
- The **power requirements** related to the product (e.g. **number & location** of electrical outlets).
- The **required operating supplies** necessary for the product and installation.
- While Agilent is delivering **Installation and Introduction** services, users of the instrument should be present throughout these services; otherwise, they will miss important operational, maintenance and safety information.
- Please consult the **Special Requirements and Other Considerations** section below for other product-specific information
- For more details, please consult the product-specific site preparation or pre-installation manual.

Important Customer Web Links

- To access **Agilent training and education**, visit <http://www.agilent.com/chem/training> to learn about training options, which include online, classroom and onsite delivery. A training specialist can work directly with you to help determine your best options.
- To access the **Agilent Resource Center** web page, visit <https://www.agilent.com/en-us/agilentresources>. The following information topics are available:
 - Sample Prep and Containment
 - Chemical Standards
 - Analysis
 - Service and Support
 - Application Workflows
- The **Agilent Community** is an excellent place to get answers, collaborate with others about applications and Agilent products, and find in-depth documents and videos relevant to Agilent technologies. Visit <https://community.agilent.com/welcome>
- Videos about specific preparation requirements for your instrument can be found by searching the **Agilent YouTube** channel at <https://www.youtube.com/user/agilent>
- **Need to place a service call?** [Flexible Repair Options | Agilent](#)

Site Preparation

Dimensions and Weight

Identify the laboratory bench space before your system arrives based on the table below. Pay special attention to the total height and total weight requirements for all system components you have ordered and avoid bench space with overhanging shelves. Also pay special attention to the total weight of the modules you have ordered to ensure your laboratory bench can support this weight.

Special notes

The following table provides dimensions and weight requirements.

Instrument Description	Weight		Height*		Depth**		Width***	
	Kg	lbs.	cm	in	cm	in	cm	in
G4567A, 7650 ALS (mounted on 7890 GC)	4.5	9.9	94	37	50	19.68	83	32.67
G4567, 7650 ALS (mounted on 7820 GC)	4.5	9.9	95	37.4	79	31.10	85	33.46

*Above bench surface of top of a 7650 injector as mounted on the noted GC.

** Extension of the 7650 injector past front of the noted GC.

*** Extension of the 7650 injector past left side of the noted GC.

Environmental Conditions

Operating your instrument within the recommended temperature ranges ensures optimum instrument performance and lifetime.

Special notes

- Performance can be affected by sources of heat & cold, e.g. direct sunlight, heating/cooling from air conditioning outlets, drafts and/or vibrations.
- The bench or supporting surface must be vibration free.
- The site’s ambient temperature conditions must be stable for optimum performance.
- Pollution degree 2, Installation Cat II

The following table may help you calculate the additional BTUs of heat dissipation from this new equipment. Maximums represent the heat given off when heated zones are set for maximum temperatures.

Instrument Description	Operating Temperature Range °C (F)	Operating Humidity Range %
G4567A, 7650 ALS	15 to 35°C (59 to 95°C)	5 to 95%

Power Consumption

Special notes

- If a computer system is supplied with your instrument, be sure to account for those electrical outlets.
- Mains supply voltage fluctuations up to \pm of the nominal voltage

Instrument Description	Line Voltage and Frequency V, Hz	Maximum Power Consumption VA	Maximum Power Consumption W
G4567A, 7650 ALS	Americas: 120 single phase (-10% / +10%) 50/60 Hz		
G4567A, 7650 ALS	200/220/230/240 Single/split phase (- 10% / +10%) 50/60 Hz		

Required Operating Supplies by Customer for Installation

Special notes

- For information on Agilent consumables, accessories, and laboratory operating supplies, please visit: <https://www.agilent.com/en-us/agilentresources>

Special Requirements and Other Considerations

- Only one 7650 ALS can be operated on a compatible GC at a time, in either the front or back injection port.
- When used on an Agilent 7890 GC, an Agilent 7693A 16 vial injector can also be operated on the back injection port with a 7650 ALS operated on the front injection port.
- The injector will not operate if it is not mounted on a GC, if the turret is not installed, or if the door is open.

Service Engineer Review (Optional)

Service Engineer Comments

If the Service Engineer completed a review of the Site Preparation requirements with the customer, the Service Engineer should complete the following Comments section. Both the Service Engineer and the customer should complete the Site Verification section below.

If there are any specific points that should be noted as part of performing the site preparation review or other items of interest for the customer, please write in this box.

Site Preparation Verification

Service Request Number:

Date of Review:

Service Engineer Name:

Customer Name:

Service Engineer Signature:

Customer Signature:

Total number of pages in this document: