

## Agilent Bio/SEC Software G7857AA \_ Site Preparation Checklist

Thank you for purchasing an Agilent instrument. To get you started and to assure a successful and timely installation of your Agilent **software**, please refer to this specification or set of requirements.

Correct site preparation is the key first step in ensuring that your instruments and software systems operate reliably over an extended lifetime. This document is an **information guide and checklist** prepared for you that outlines the supplies, consumables, space and utility requirements for your equipment for your site.

### Customer Responsibilities

**Make sure your site meets the following prior specifications before the installation date. For details, see specific sections within this checklist, including:**

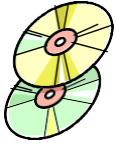
- The computing environment and the necessary space is made available
- The number & location of electrical outlets for your computer systems and peripherals are planned
- That your site meets the software, hardware and networking specifications below
- Locate your sales order information, software authorization codes and/or software licenses/certificates
- The necessary software media, disks etc are available including upgrade/update disks
- That a suitable backup solution is identified for your software
- Availability of a system/network administrator as needed to connect to your intranet
- Please consult Other Requirements section below for other product-specific information.
- Complete Final Check: Software Site Preparation Tool. Please visit the following Agilent website to download the Software Installation Site Preparation Tool.  
<http://www.chem.agilent.com/en-US/Support/Downloads/Utilities/SWSitePrepTool/Pages/Default.aspx>  
This easy-to-use tool verifies that your PC meets the necessary hardware and software pre-requisites before installing the Agilent software.

**If Agilent is delivering installation and familiarization services, users of the instrument should be present throughout these services; otherwise, they will miss important operational, maintenance and safety information.**

- Author to add specific customer activities or additional help, information, or guidance for the customer to follow prior to installation>.

### Important Customer Information

1. If you have questions or problems in providing anything described as a Customer Responsibilities above, please contact your local Agilent or partner support/service organization for assistance prior to delivery. In addition, Agilent and/or its partners reserve the right to reschedule the installation dependent upon the readiness of your laboratory.
2. Should your site not be ready for whatever reasons, please contact Agilent as soon as possible to re-arrange any services that have been purchased.
3. Other optional services such as additional training, operational qualification (OQ) and consultation for user-specific applications may also be provided at the time of installation when ordered with the system, but should be contracted separately.

**Agilent Bio/SEC Software G7857AA \_ Site  
Preparation Checklist**

**Software Specifications for  
Workstations, Clients, and Servers**
**Special Notes:**

1. Upgrade system to the latest MS Windows Service Pack

Specification Description	Minimum	Recommended (if applicable)	Comments (if applicable)
Operating system type(s), versions	* Microsoft® Windows® 7 SP1 Professional & Enterprise (32 bit and 64 bit). updated to latest Windows patches  * Microsoft® Windows® 8.1 Professional & Enterprise (64 bit) updated to latest Windows patches	* Microsoft® Windows® 10 Professional (64 bit). updated to latest Windows patches	* The USB driver supplied for the Agilent 1260 Infinity MDS is compatible with Microsoft® Windows® 10, A Microsoft® Windows® 8.1 and Microsoft® Windows® 7.
O/S .NET & other add-ons	.NET 3.5. 4.0 and Visual C++ ® tools if required		Installed as part of Bio-SEC
Language settings/compatibility	English US, UK	English US, UK	
Regional settings/compatibility	N/A		
Account settings/privileges	Full admin rights		
Specific drivers	N/A		


**Computer Hardware Specifications for  
Workstations, Clients, and Servers**
**Special Notes**

1. Author to add special considerations or notes.

Specification Description	Minimum	Recommended (if applicable)	Comments (if applicable)
Processor type & speed	Pentium 4 operating at 3.2 GHz or greater microprocessor	Pentium 4 operating at 3.2 GHz or greater microprocessor	

**Agilent Bio/SEC Software G7857AA \_ Site  
Preparation Checklist**

Memory	4 GB (RECOMMENDED 8 GB)	8GB	
Internal Storage/devices/media	>250GB	250GB+	
External Storage/devices/media	N/A		
Video devices	17" 1280x1024 resolution (SXGA)	19", 1440x900)	
Audio devices	N/A		
Printing devices	N/A		
Pointing devices	Microsoft compatible mouse		
Additional devices/ports	Minimum 2 USB ports.	2 USB ports +	


**Networking Specifications**
**Special Notes**

1. Author to add special considerations or notes.

Specification Description	Minimum	Recommended (if applicable)	Comments (if applicable)
Network type, bandwidth, speed, protocol etc	N/A		
Additional network or instrument devices/cards requirements	N/A		

**Important Customer Web Links**

- For additional information about our solutions, please visit our web site at <http://www.chem.agilent.com/en-US/Pages/HomePage.aspx>
- Need to get information on your product?  
Literature Library - <http://www.agilent.com/chem/library>
- Need to know more?  
Customer Education – <http://www.agilent.com/chem/education>
- Need technical support, FAQs? – <http://www.agilent.com/chem/techsupp>
- Need supplies? – <http://www.agilent.com/chem/supplies>