



Agilent InfinityLab Ultivo LC/TQ

Site Preparation Guide







Unpacking

Delivery and Installation

Do not open any shipping containers until an Agilent representative is present. Warranty claims for missing items will not be honored unless an Agilent representative is present to verify the contents of each container as it is unpacked. The actual shipping containers become your property and should not be returned to Agilent Technologies.

If the product is coming from a cold storage environment, allow time for thermal equilibration to room temperature before opening the shipping containers. This will avoid exposing the instrument to condensing humidity.

1	If the product is coming from a cold storage environment, allow time for thermal equilibration to room temperature before opening the shipping containers. This will avoid exposing the instrument to condensing humidity.	 A grayscale photograph of a laboratory interior. In the foreground, a person is operating a forklift, moving a large, light-colored shipping container. The container has the Agilent Technologies logo and some handling symbols on its side. In the background, there are laboratory benches with various pieces of equipment and glassware.
2	If the product is coming from a cold storage environment, allow time for thermal equilibration to room temperature before opening the shipping containers. This will avoid exposing the instrument to condensing humidity.	 A grayscale photograph of a laboratory interior. In the foreground, a large, light-colored shipping container sits on a wooden pallet. The container has the Agilent Technologies logo on its side. To the right of the container, there is a circular clock icon with a red hand pointing to the 12 o'clock position, indicating a time period for thermal equilibration.

3	Do not open any shipping containers until an Agilent representative is present. Warranty claims for missing items will not be honored unless an Agilent representative is present to verify the contents of each container as it is unpacked.	
4	Do not open any shipping containers until an Agilent representative is present. Warranty claims for missing items will not be honored unless an Agilent representative is present to verify the contents of each container as it is unpacked.	
5	The actual shipping containers become your property and should not be returned to Agilent Technologies.	