



# Agilent InfinityLab Ultivo LC/TQ

## Site Preparation Guide

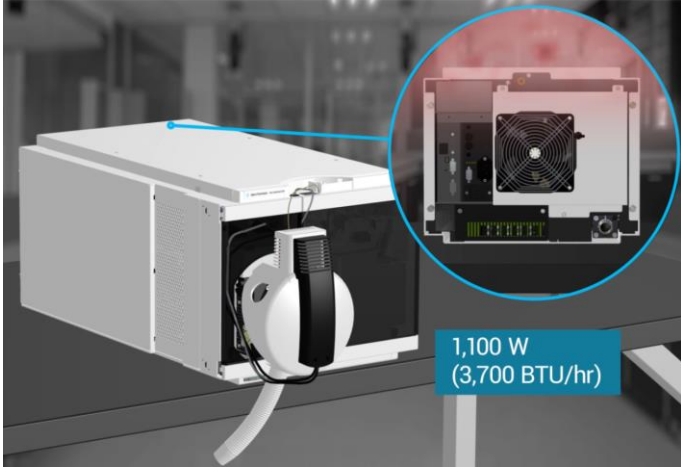
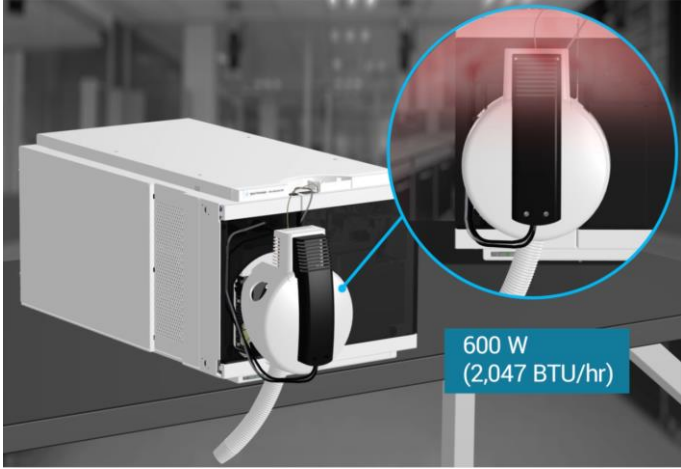


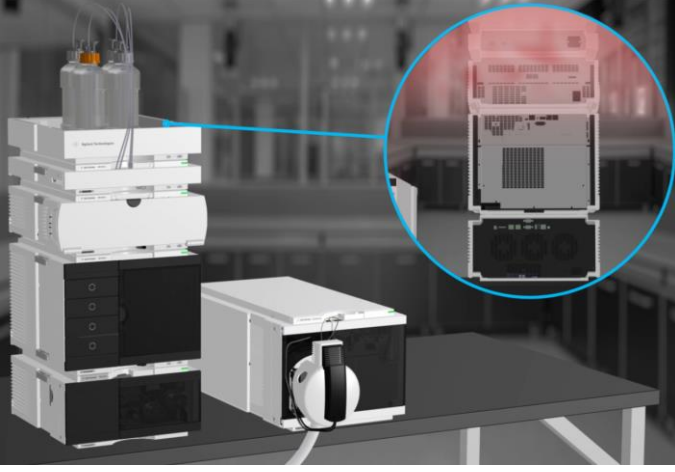
## **Temperature and Humidity**

The Ultivo LC/TQ is specified for operation under the following conditions:

- 15 to 35°C (59 to 95°F)
- Constant temperature (< +/-3°C from calibration temperature)
- < 85% relative humidity at 35°C
- Non-condensing, non-corrosive atmosphere

Environmental control systems must maintain these temperatures and humidity ranges.

1	The Ultivo LC/TQ system dissipates up to 1,100 Watts (3,700 BTU/hr).	
2	Approximately 600 Watts (2,047 BTU/hr) are removed with the source exhaust.	

3	<p>The LC and data system also contribute significantly to the cooling load. The exact amounts will depend on their configurations.</p>	
4	<p>The LC and data system also contribute significantly to the cooling load. The exact amounts will depend on their configurations.</p>	