



# **Agilent 6500 Series Q-TOF LC/MS**



## **Site Preparation Guide**






## **Unpacking**

Do not open any shipping containers until an Agilent representative is present. Warranty claims for missing items will not be honored unless an Agilent representative is present to verify the contents of each container as it is unpacked.

The actual shipping containers become your property and should not be returned to Agilent Technologies. The flight tube shipping insert is also your property and does not need to be returned to Agilent Technologies.

1	<p>If the product is coming from a cold storage environment, allow time for thermal equilibration to room temperature before opening the shipping containers. This will avoid exposing the instrument to condensing humidity.</p>  A grayscale photograph showing a person operating a forklift in a warehouse setting. The forklift is carrying a large, rectangular shipping container. The background shows industrial shelving and equipment.
2	<p>If the product is coming from a cold storage environment, allow time for thermal equilibration to room temperature before opening the shipping containers. This will avoid exposing the instrument to condensing humidity.</p>  A grayscale photograph of a large, rectangular shipping container in a warehouse. The container has the Agilent Technologies logo on its side. A circular clock overlay is positioned to the right of the container, with a red hand pointing to the 12 o'clock position.

<p>3 Do not open any shipping containers until an Agilent representative is present. Warranty claims for missing items will not be honored unless an Agilent representative is present to verify the contents of each container as it is unpacked.</p>	
<p>4 Do not open any shipping containers until an Agilent representative is present. Warranty claims for missing items will not be honored unless an Agilent representative is present to verify the contents of each container as it is unpacked.</p>	
<p>5 The actual shipping containers become your property and should not be returned to Agilent Technologies.</p>	

- 6 The actual shipping containers become your property and should not be returned to Agilent Technologies.

