

Agilent CrossLab Start Up Services

Agilent OpenLab CDS v3.0 Workstation / Workstation Plus Software Site Preparation Checklist

Thank you for purchasing software from **Agilent Technologies**. CrossLab Start Up is focused on helping customers shorten the time it takes to start realizing the full value of their instrument system investment.

Correct site preparation is the key first step in ensuring that your instruments and software systems operate reliably over an extended lifetime. This document is an **information guide and checklist** prepared for you that outlines the software requirements for the system set up in your lab.

Introduction

Customer Information

- If you have questions or problems in providing anything described as part of Customer Responsibilities below, please contact your local Agilent or partner support / service organization for assistance prior to delivery. In addition, Agilent and/or its partners reserve the right to reschedule the installation dependent upon the readiness of your laboratory.
- Should your site not be ready for whatever reasons, please contact Agilent as soon as possible to re-schedule any services that have been purchased.
- Other optional services such as additional training, operational qualification (OQ) and consultation for user-specific applications may also be provided at the time of installation when ordered with the system but should be contracted separately.

Customer Responsibilities

Ensure that your site meets the following specifications before the installation date. For details, see specific sections within this checklist, including

- The **computing environment** and the necessary space is made available.
- The number and location of **electrical outlets** for your computer systems and peripherals are planned.
- Your site meets the **software, hardware, and networking specifications** described later in this document.
- Locate your **sales order information** and the **software entitlement email** sent to the person named as the software contact person during purchase.
- Check for Hotfix, Microsoft updates, and patch compatibility appropriate for the customer's system. **Critical Note: Current Microsoft updates and Windows version must be installed on the customer provided PC before installation can begin.**
- The necessary **software media** is available, including upgrade/update media. Note: physical software media is only delivered if requested at the time of purchase. Software may be downloaded from the customer's **Agilent Online Account** and should be downloaded prior to the installation.
- That a suitable **backup solution** is identified for your software.

- If Agilent is delivering **Installation and Introduction services**, users of the instrument system should be present throughout these services. Otherwise, they will miss important operational, maintenance, and safety information.
- Availability of a **system/network administrator** as needed to connect to your intranet.
- Please consult the **Special Requirements** section for other product-specific information.
- Complete Final Check: Run the System Preparation Tool (SPT). On the OpenLab CDS 3.0 installation media run the *SystemPreparationTool.exe* under \Setup\Tools\SPT to complete the final check.

The System Preparation Tool (SPT) checks and applies Windows settings on your machine. The majority of the steps are automated by the tool. Applying these settings helps avoid issues during installation and use of the product. Please read the System Preparation Report and manually make the mandatory and recommended adjustments that are outlined in the report.

Important Customer Web Links

- To access Agilent training and education, visit <http://www.agilent.com/chem/training> to learn about training options, which include online, classroom and onsite delivery. A training specialist can work directly with you to help determine your best options.
- To access the **Agilent Resource Center** web page, visit <https://www.agilent.com/en-us/agilentresources>. The following information topics are available:
 - Sample Prep and Containment
 - Chemical Standards
 - Analysis
 - Service and Support
 - Application Workflows
- The **Agilent Community** is an excellent place to get answers, collaborate with others about applications and Agilent products, and find in-depth documents and videos relevant to Agilent technologies. Visit <https://community.agilent.com/welcome>
- Videos about specific preparation requirements for your instrument can be found by searching the **Agilent YouTube** channel at <https://www.youtube.com/user/agilent>
- OpenLab CDS 3.0 uses **Agilent Electronic Licensing**, allowing all licensing entitlements to be viewed and managed through the Licenses & Agreements page of your **Agilent Online Account**. You can access your **Agilent Online Account** by visiting <https://www.agilent.com>, logging in with your account credentials, expanding your account dropdown, and selecting the "Dashboard" option. Software installation media and updates can be downloaded by following the "Download/Activate software" link for the associated Entitlement.
- Some OpenLab instrument drivers and add-ons use **Agilent SubscribeNet** to manage and administer their licenses and provide users with access to download the latest software/updates.
To access **Agilent SubscribeNet**, visit <https://agilent.flexnetoperations.com>.
- **Need to place a service call?** [Flexible Repair Options | Agilent](#)
- Links to the Software Status Bulletin (SSB) and Software Release Bulletin (SRB) are in the Software Bulletins.pdf located under Setup\Tools\Support\History on the installation media.
- The full OpenLab Help & Learning system with most current Help and Resources is available via <https://openlab.help.agilent.com>. The Online version features localized User Manuals.

Site Preparation

Software Specifications for OpenLab CDS standalone Workstation

Special notes

- If you have purchased a system including hardware instrumentation, refer to the instrument Site Preparation Checklist for requirements regarding laboratory bench space requirements.

Specification Description	Details
Operating system name, version	Windows 11, Enterprise or Professional, 64-bit (24H2 or greater) Windows 10, Enterprise or Professional, 64-bit (21H2 or greater)
.NET Framework (64-bit)	.NET 3.5.x (<i>must be enabled on Windows</i>) and .NET 4.8 or higher (<i>installed by OpenLab CDS v3.0 Installer</i>)
.NET (64-bit) Desktop Runtime	6.0 and 8.0 (<i>installed by OpenLab CDS v3.0 Installer</i>)
Asp.NET Runtime	6.0 and 8.0 (<i>installed by OpenLab CDS v3.0 Installer</i>)
Web browser	Google Chrome 98 or higher Microsoft Edge
Antivirus Software	Microsoft Defender Trellix MVISION Plus Symantec Endpoint Security CrowdStrike
Language settings/compatibility	English Western European Language (<i>CDS will always appear in English</i>) Chinese (<i>Non-localized instrument drivers are supported and will always appear in English</i>) Japanese (<i>Non-localized instrument drivers are supported and will always appear in English</i>) Brazilian Portuguese (<i>Non-localized instrument drivers are supported and will always appear in English</i>)
Account settings/privileges	Domain user with local administrator privilege required for installation and configuration

Computer Hardware Specifications for Workstations

Specification Description	Workstation	Workstation Plus
Processor type and speed	Intel® i5, i7, or Xeon E3 or equivalent 3.0 GHz or greater, 4 Core	
Physical memory (RAM)	1-2 instrument points: 8 GB 3 or more instrument points: 16 GB	16 GB
	Ensure that at least 4 GB is reserved for the Windows operating system.	
Hard disk	1 x 500 GB 7200 RPM SATA drive minimum or equivalent solid-state drive. SSD is recommended for better performance.	1 TB or 2x 500 GB 7200 RPM SATA drive (minimum). If the computer has a disc array controller we recommend 2 x 1 TB in RAID1, and SSD drive.
RS-232-port	1 serial port required for selected instruments that are still using RS-232 communication. See instrument specifications for details.	
USB Port	USB 3.0 might be required for installation, if USB media was ordered	
LAN Card	100 MB/1 GB LAN for instrument control 2nd LAN card required for house, to isolate the instruments data traffic from the lab intranet connection.	
Graphic resolution	1600 x 900 minimum 1920 x 1080 recommended	

Computer Hardware Specifications for Workstations with any LC/(Q-)TOF

NOTE: Standalone Workstation with Content Management (Workstation Plus) is not supported with LC/(Q-)TOF or MassHunter DA

Specification Description	Workstation
Processor type and speed	Intel® Xeon W3-2423 (4.2 GHz, 15 MB cache, 6 cores) or equivalent 4.2 GHz or greater, 6 Core or more
Physical memory (RAM)	32 GB minimum, 64 GB recommended (for one LC /(Q)TOF instrument)
Hard disk	1 TB M.2 NVMe SSD Boot Volume and 12 TB 7200 RPM SATA 6G Hard Drive
RS-232-port	1 serial port required for selected instruments that are still using RS-232 communication. See instrument specifications for details.
USB Port	USB 3.0 might be required for installation, if USB media was ordered
LAN Card	10 GbE LAN for instrument control of G6575A. 1 GbE LAN for G6549A or G6230C. 2nd LAN card required for lab intranet connection, to isolate the instrument's data traffic from the lab intranet connection.
Graphic resolution	1600 x 900 minimum; 1920 x 1080 recommended

Networking Specifications

Special notes

When using LAN communications to connect Workstation / Workstation Plus to an instrument, use one of these methods:

- Connect via an isolated switch using standard CAT-5E or CAT-6 network cabling.
- LAN communication hardware should be 100/1000 Mbps (or higher) speed capable. 10 GbE required for some LC/(Q-)TOF configurations. The J4100 Jet Direct Card is not supported. Use a G1369 LAN interface card instead.
- NIC teaming (See *Note 1*): LAN cards should not be teamed on Workstation / Workstation Plus.
- LAN communication must be on the same subnet as instruments, and preferably on the same segment.

Note 1: Network Interface Card (NIC) teaming is also known as Load Balancing and Failover (LBFO)

Specification Description	Supported
Network type, bandwidth, speed, protocol, etc.	Internet Protocol Version 4 (TCP/IPv4) only <i>Internet Protocol Version 6 (TCP/IPv6) is not supported</i>
IP Address	Static or DHCP Reservation

Special Requirements

- Please refer to the **CDS_v3.0_Requirements_en.pdf** on the installation media for supported hardware, software, instruments, and firmware requirements.
- If the customer uses a Firewall other than the Windows Firewall, ensure that all required ports can be used by OpenLab CDS. For details, see the *Firewall Settings* section of *Chapter 3 Network Requirements* in the *OpenLab CDS Workstations, Clients, and Instrument Controllers Requirements and Supported Instruments (CDS_v3.0_Requirements_en.pdf)*.

Service Engineer Review (Optional)

Service Engineer Comments

If the Service Engineer completed a review of the Site Preparation requirements with the customer, the Service Engineer should complete the following sections below.

If there are any specific points that should be noted as part of performing the service review or other items of interest for the customer, please write in this box.

Site Preparation Verification

Service Request Number:

Date of Review:

Service Engineer Name:

Customer Name:

Service Engineer Signature:

Total number of pages in this document: