

**ChemStore C/S G2181BA Standalone or client, revision B.04.02****ChemStation Plus Security Pack G2183AA, revision B.04.02****Site Preparation Specification****Purpose of Procedure**

To assure that the installation of Agilent instruments and systems can be completed successfully by careful preparation and evaluation of the installation site and by ensuring the availability of appropriate utilities, consumables and supplies.

**Customer Responsibilities**

Customers should ensure that all necessary operating supplies, consumables and usage dependent items such as columns, vials, syringes and solvents required for the successful installation of instruments and systems are available. In addition, customers must provide the Agilent Support Provider with appropriate administrator-level access to the system during the installation. Installation sites should be prepared in accordance with the following specifications.

**Important Information**

If you have problems in providing any of the following, please contact your local Agilent office for assistance. Assistance with user specific applications may be provided but should be contracted separately. Users of the instrument should be present throughout the installation and familiarization otherwise important operational, maintenance and safety information may be missed.

**Procedure Checklist****A) ChemStore C/S Requirements****PC Hardware Requirements**

- Pentium IV processor operating at 1.5 GHz MHz or higher
- Memory: At least 512 MB (XP) or 1 GB (Vista, Windows 7)
- Hard-disk with at least 4 GB free disk space
- Display: min. resolution 1280 x 1024, more than 256 colors
- For ChemStore clients: Network Interface Card and 10 BaseT network connection with a speed of 10 MB/s, at least.
- A backup device or network-based backup strategy

**Software Requirements Operating System**

- Windows XP Professional with Service Pack 3
- Or Windows Vista Business with Service Pack 2
- Or Windows 7 Business (32-bit)
- Or Windows 2003 server R2 with Service Pack 2
- Internet Explorer 7.0 or 8.0
- User logon with Administrative capabilities

**Software Requirements Chromatographic Data system**

- ChemStation B.04.02 or higher for
  - A/D (G2172BA)
  - CE or CE/MSD (G1601B/G2201BA)
  - GC (G2070BA)
  - LC or LC/MSD (G2170BA,G2710BA)

**B) Additional Requirements for Security Pack****The following requirements only apply if Security Pack G2183AA or G2187AA is installed on top of ChemStore C/S**

- At least one NTFS partition configured
- For PCs other than Agilent Bundle PCs: Setup folder permissions as outlined in the document 'Configure XP.pdf' on the ChemStation installation CD.
- At least one user logon as member of the user group "Administrators" and one user logon as member of the user group "Users" set up.
- Members of the user group "Users" must not be member of the "Administrators" group