



Title: Patch 2 for MassHunter for ICP-MS G7201D D.01.03

Patch Number: G7201D D.01.03.02

Supersedes: False

Applies to: MassHunter for ICP-MS G7201D

Revision: D.01.03

Summary: Patch 2 for ICP-MS MassHunter (G7201D, revision D.01.03) has been released.

To Be Installed By: Agilent-Qualified Personnel or Customer

Localized: Yes

Prerequisites:

Installation of MassHunter for ICP-MS G7201D D.01.03 Workstation.

Situation: Patch 2 for MassHunter for ICP-MS G7201D D.01.03 ("Patch.msp") resolves the following issues.

This Patch resolves following issue/s:

Patch.msp

KPR#	Issue ID	Description
1145523	ICPMH-12910	E-Signature information is created for each analysis result. However, when an analysis result is copied using "Save As", the copied result is linked to the E-Signature information of the original result.

New features:

n/a

**Solution/Action:**

Solution recommended for All customers (Yes/Specific Failure Only): **Yes**

The issues listed above are resolved with this patch for MassHunter G7201D D.01.03
This software patch should be applied to all MassHunter systems G7201D D.01.03.

Patch and Instruction Download

Agilent.com Home.
Then search with G7201D

<https://www.agilent.com/en/support/software-informatics/atomic-spectroscopy-data-systems/icp-ms-masshunter-software/g7201d-d0103-patch02>

Additional Information:

- "Patch.msp" includes all fixes for English, Chinese, and Japanese versions.

Compliance Information:**Validation Notes:**

Agilent Technologies confirms that the issues listed in the situation section are resolved and that no related functionality is adversely impacted after the application of this patch. This patch was developed, tested, and passed according to the Agilent Technologies Patch Release Process.

Risk Assessment:

This update patch has been developed to specifically address the issues as described above. Agilent has evaluated the possibility of adverse effects to other functions through our release processes and has found the risk to be very low.

Recommended Validation tasks:

Customers should verify the update patch is installed correctly by running the IQT tool on all machines where the software update was applied. Instructions on how to run this are listed in the installation procedure below. Any additional validation requirements would be in accordance with procedures specific to a customer site based on the changed noted in this document.

Installation:

This patch must be installed on MassHunter Workstation.

Installation Instructions:

Please refer to the Instruction Procedure document which also can be downloaded.

Files Updated with Patch 2 (Patch.msp):

Root Path\Path\File Name	Version
C:\Program Files\Agilent\ICPMH\Bin\IcpDaAppLogic.dll	17.1.748.12
C:\Program Files\Agilent\ICPMH\Bin\IcpMHServiceProvider.exe	17.1.748.12
C:\Program Files\Agilent\ICPMH\Bin\SIVersion200.dll	17.1.748.12

Files Added with Patch 2:

Root Path\Path\File Name	Version

Uninstall Supported (Y/N): **Y**

Patch Release Date:	16 Dec 2024	Document Version:	1.0
Disclaimer			
THIS DOCUMENT AND ACCOMPANYING SOFTWARE MODULES ARE PROVIDED "AS IS" WITH THE SAME WARRANTY CONDITIONS SPECIFIED IN THE SOFTWARE LICENSE AGREEMENT FOR THE LICENSED PRODUCT.			
©2024 Agilent Technologies Printed in Japan			