

**Removing and Re-Installing
ChemStore and Security Pack Software
on Microsoft Windows NT 4.0
G2181BA & G2183AA, Rev. B.01.03-B.02.01**

Warning:

USE EXTREME CAUTION. The following procedure is intended to aid in the reloading of software for the Agilent Technologies family of ChemStation Plus products and potentially involves the modification of Windows system files and renaming of the current ChemStation Plus application subdirectories.

AGILENT TECHNOLOGIES MAKES NO WARRANTY OF ANY KIND WITH REGARD TO THIS MATERIAL, INCLUDING, BUT NOT LIMITED TO, IMPLIED WARRANTIES OF MERCHANTABILITY AND FITNESS FOR A PARTICULAR PURPOSE. AGILENT TECHNOLOGIES SHALL NOT BE LIABLE FOR ERRORS CONTAINED HEREIN., OR FOR INCIDENTAL OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH THE FURNISHING, PERFORMANCE, OR USE OF THIS MATERIAL.

This procedure assumes the user has knowledge of using a mouse and Windows programs.

Please read all instructions, first. If you do not understand any sections, **DO NOT PERFORM THIS PROCEDURE!!**

NOTE: If you need to remove and re-install your ChemStation software, please refer to the following document:


[ChemStation Re-Installation Procedure](#) (A10572.pdf)

NOTE: If you want to remove ChemStore from your system, and do not intend to re-install it, please see "[Removing ChemStore and Security Pack completely from a system](#)" at the end of this document.

This document is believed to be accurate and up-to-date. However, Agilent Technologies, Inc. cannot assume responsibility for the use of this material. The information contained herein is intended for use by informed individuals who can and must determine its fitness for their purpose.


Remove the ChemStore and Security Pack software

1. Log into Windows as an administrator. If you are operating in a domain environment, log in as a domain administrator.
2. From the Windows® Taskbar, select the **Start** button.
3. From the Start menu, select **Settings**.
4. From the Settings menu, select **Control Panel**.
5. Once the Control Panel is displayed, double-click on **Add/Remove Programs**. You should now have a panel displayed that looks like this:

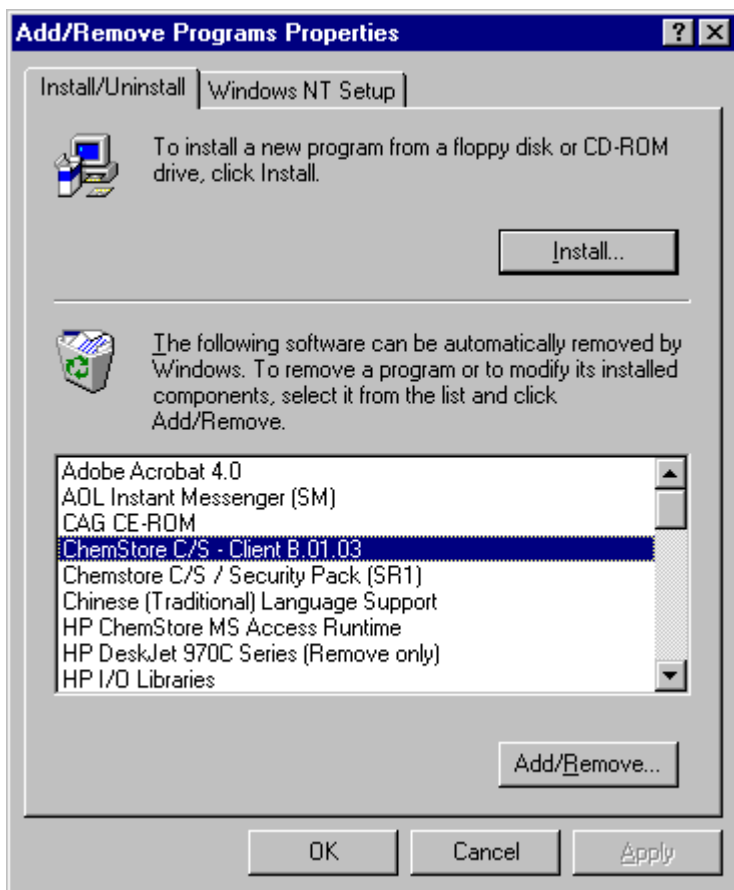

6. Highlight the ChemStore C/S – Client B.01.03 line (as in the above picture), and click on the **Add/Remove** button.
7. When the uninstallation program prompts you, asking if you are sure you want to uninstall ChemStore, click on the **Yes** button.

Please note: You may receive a panel that looks similar to the following:

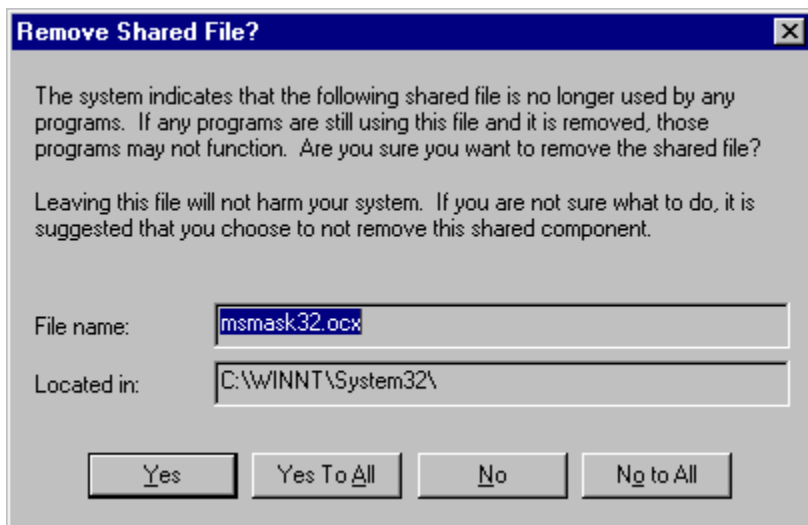


If so, click on the “**Yes to All**” button.

8. Once the software removal has completed, click on the **OK** button to exit.
9. If you are using ChemStore C/S Security Pack software, remove it from the system using the same process as above in steps 6 through 8.

Rename the ChemStore application subfolder

1. From the Windows Taskbar, select the **Start** button.
2. From the **Start** menu, select **Find**, and then **Files or Folders...**
3. Drop down the Look In selections, and select **My Computer**.
4. Enter “chemstor” in the Named: field (omitting the quotation marks). Note that there is no ‘e’ at the end of “chemstor”.
5. Click on the **Find Now** button.
6. In the results display table, find the ChemStor folder, and note the full path name for the folder (e.g. C:\HPCHEM\CHEMSTOR).
7. Right-click on the Chemstor folder, and select **Rename**.
8. Enter “chemstor.old” in the name field that is now active (omitting the quotation marks), and hit the **Enter** key. You may receive a warning message similar to the following:

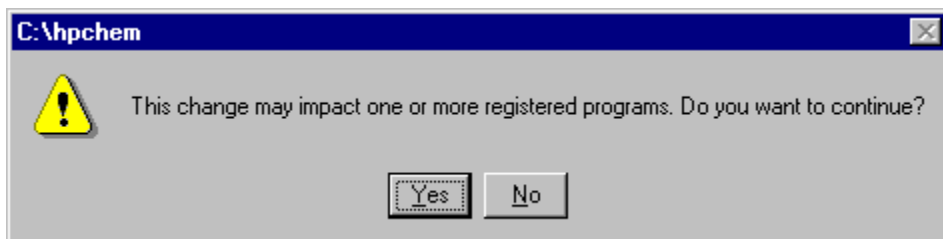


Click **Yes** to continue the renaming operation. You should now see the name displayed as chemstor.old on the folder.

9. Reboot your computer.

Re-Installing ChemStore and Security Pack

The instructions for installing ChemStore and Security Pack are contained in Agilent ChemStore C/S Installation Guide that was provided with your software. For more information, you may wish to refer to Chapters 1 or 2 for Standalone or Client/Server installation instructions, respectively. Here is a description of the installation process:

1. Find the appropriate installation CD. This should be labeled “ChemStation Plus Security Pack, Security Pack for UV-vis and ChemStore C/S”. It will also have a large 21 CFR Part 11 logo on it.
2. Insert the CD into the CD-ROM drive.
3. From the Windows Taskbar, click the **Start** button.
4. Select **Run...** from the menu.
5. Enter <drive>:\G2181_setup.exe in the Open: field (where <drive> is the letter that designates the CD-ROM drive), and click **OK**.
6. ChemStore software installation will begin. Click **Next** after the initial display screen, and **Yes** at the software license agreement. When prompted, enter your software license number and click **Next** again.
7. When prompted for the path, check to make sure that the path is the same as it was in Step 6 of “[Rename the ChemStore application subfolder](#)”, above (excluding the ChemStore portion of it), and click **Next**. The software will begin to copy files to the hard drive of your computer.
8. A panel will now display with two select boxes that must be checked in order for the correct software to be installed. It looks like this:

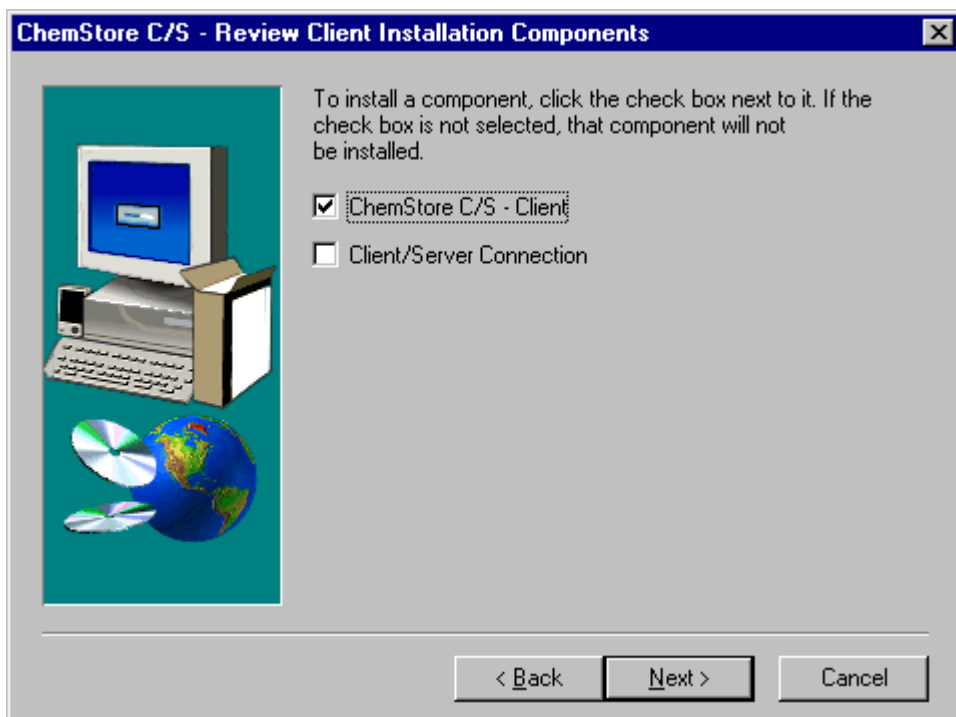

If you are re-installing a standalone ChemStore system, leave the boxes as they are when they come up. If you are re-installing a client/server ChemStore client, then both boxes should be checked. Click **Next** twice to continue. The system will begin to copy files.

9. You may receive a message indicating the time-zone variable TZ has been modified. Click **OK**
10. If you are performing a re-installation of client/server related system, you will be prompted for the name of the server and database instance being used. Please contact your system administrator for this information. Click **Next** twice to continue installation.
11. Reboot your computer when prompted. Installation is complete.

Removing ChemStore and Security Pack completely from a system

Follow the instructions in "[Remove the ChemStore and Security Pack software](#)" above. Complete through step 5 of "[Rename the ChemStore application subfolder](#)", and proceed with the following:

6. In the results display table, find the ChemStor folder, and right-click on it. Select **Delete** from the menu. Click **Yes** on the delete confirmation panel. The ChemStore program has now been removed from your system.
7. Back up your win.ini file according to the following procedure:

- a) Place a formatted floppy disk in the floppy drive of the computer you are working on.
 - b) From the Windows Taskbar, select the **Start** button
 - c) From the Start menu, select **Find**, and then **Files or Folders...**
 - d) Drop down the **Look In** selections and select **My Computer**
 - e) In the Named field, enter "win.ini" (excluding the quotation marks)
 - f) Click on the **Find Now** button
 - g) In the results display table, right-click on the win.ini file and select **Send To...**
 - h) From the **Send To...** menu, select **3½ Floppy (A)**.
 - i) After the file is copied, close the **Find: Files** panel.
8. From the Windows Taskbar, select the **Start** button.
 9. Select **Run...** from the menu.
 10. Enter "sysedit" (without the quotation marks) in the Open: box and hit the **enter** key.
 11. Find the panel for 'win.ini' and click on the top to bring it forward.
 12. Click on **Search...** and select **Find** from the menu.
 13. In the Find: field, enter "PCS" (excluding quotation marks) and click the **Next** button.
 14. Scroll down in the section below the [PCS] line, and find a line that begins with "ChemStore C/S=". Delete this entire line.
 15. Hit the **F3** key to find the next PCS section (usually labeled as [PCS,1]).
 16. Scroll down in the section below the [PCS,1] line, and find a line that begins with "Addons=". If the value that Addons is set to is 1 (the line reads Addons=1), then delete the entire line. If not, set the value to one less than its current value (set it to 1 if the value is currently 2).
 17. In the same section, find the line that begins with "Addon1=". If this line is set to a path value that includes the file name 'hpdbif00.mcx', delete the entire line.

If the "Addons=" line was set to a value of >1, then there is more than one "AddonX=" (where X is a number) line in this section. Delete the line containing the above file name, but be sure that the remaining "AddonX=" lines begin with X having the value of 1, and increasing as consecutive integers (e.g. Addon1=, Addon2=, etc.).
 18. There may be a line in this section beginning with "ChemStore C/S =". If so, delete the entire line.

19. Repeat this procedure (steps 15 through 18) until you find no more [PCS,X] sections (where X is a number from 1 to 4).
20. Select **File** from the menu at the top of the panel, and select **Save**.
21. Exit System Editor by selecting **File**, and then **Exit**.

NOTE: You may, instead of deleting the lines, add a semicolon (;) at the beginning of each line. This has the effect of causing these lines to be ignored.