



Agilent Technologies

Innovating the HP Way

G1600A CE Electrochromatography System Preparation

This document is believed to be accurate and up-to-date. However, Agilent Technologies, Inc. cannot assume responsibility for the use of this material.


The information contained herein is intended for use by informed individuals who can and must determine its fitness for their purpose.

Preparing the Agilent CE Instrument for High-Pressure Use

- 1 Ensure gas source is switched off prior to connection.
- 2 Connect push-fit connector and gas line to the receptor at the rear of the Agilent CE instrument.
- 3 Connect the gas line to the source.
- 4 Switch on gas supply.
- 5 Select CEC from the CE mode selection menu or by clicking on the option button next to CEC (Figure 60).

Figure 60

CE Mode Selection Menu

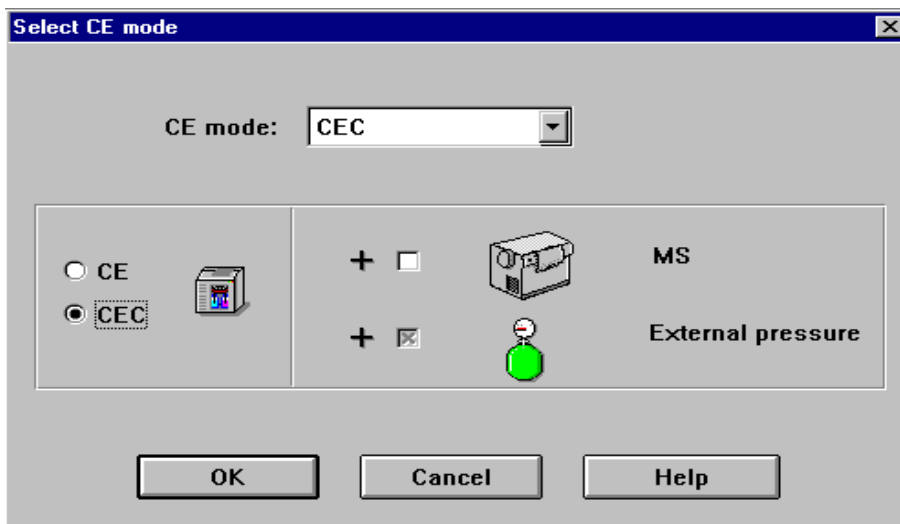

- 6 A new icon will appear on the GUI and the Mode field will indicate CEC (Figure 61).

Capillary Electrochromatography

Preparing the Agilent CE Instrument for High-Pressure Use

Figure 61

GUI

