

How to Install the Patch 003 File for ICP-MS MassHunter Rev.B.01.01

1 Install the Patch 3

- 1.1 Put the 7700 / 7500 in shutdown mode from MassHunter.
- 1.2 Close MassHunter.
- 1.3 Copy patch file “g7200b_0101_003_en.exe” to the MassHunter PC (for English version PC; for Japanese version, copy “g7200b_0101_003_ja.exe”).
- 1.4 Double click the patch file.
- 1.5 Follow the onscreen instructions.

Note:


Installation status bar pauses during the installation because the dialog box to install the factory data is waiting for an action. The dialog box may be hidden behind the installation status dialog box, clicking “**Cancel**” in the dialog box to install the factory data will allow the installation to continue.

2 Install Excel Add-in for LabQC

- 2.1 Click on **[Start]** on the Windows Task bar and select **[All Programs] > [ICP-MS MassHunter Workstation] > [Install Excel Add-in for LabQC]**.
- 2.2 Install the program following the instructions on the installer.

3 Confirm Build No.

- 3.1 Open **[OffLine Data Analysis]**.
- 3.2 Select **[Help] > [About]**, and the splash screen will appear indicating the Version, Build number and Patch number as follows.

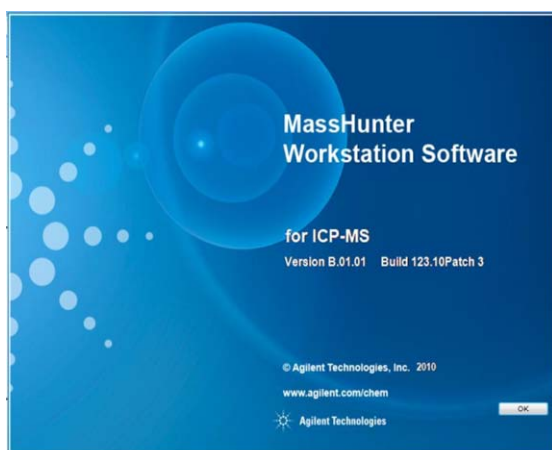

- 3.3 Click **[OK]** to exit the screen.

- 4 Switch the 7700/7500 power on.
- 5 7700 or 7500 status indication initially shows Red after turning on the power, then it is turned off after a few minutes. After indicator turns off, click ICP-MS Top.