

The Refractive Index Calibration Procedure

When required	Recommended after exchange of the optical unit or RIM board.
Tools required	Laboratory balance
Parts required	DAB/Ph Eur/BP/JP/NF/USP Grade Sucrose Syringe 9301-1446 Syringe needle 9301-0407 Sample filter 5061-3367 PEEK fitting 0100-1516

Steps

- 1** Preparation of the Sucrose calibration solution.
- 2** Preparing the pump.
- 3** Start Refractive Index Calibration.
- 4** Flushing the degasser and pump.
- 5** Purging the sample and reference cells.
- 6** Fill the sample cell with calibration solution.
- 7** Calibrate Refractive Index.

Preparation of the Sucrose sucrose Calibration Solution

- 1** To prepare 25 ml of the calibration solution 87.5 mg of the Sucrose sample is required.
- 2** Add the weighed amount of sample into a suitable volumetric flask.
- 3** Dispense 10 ml of LC grade water into the flask and shake to dissolve.
- 4** Dilute the contents of the flask to volume with LC grade water.
Wait five minutes and shake again. The solution is now ready for use.

Preparation of the Pump

- 1** Fill a suitable solvent bottle with LC grade water.
- 2** Connect this bottle to Channel A of the pump, A1 if a binary pump.



Agilent Technologies

Innovating the HP Way

Calibration Procedure for the 1100 Series Refractive Index Detector

This document is believed to be accurate and up-to-date. However, Agilent Technologies, Inc. cannot assume responsibility for the use of this material.
The information contained herein is intended for use by informed individuals who can and must determine its fitness for their purpose.

Windows® and Windows NT® are registered trademarks of Microsoft Corp
a25879c.doc

[http:// www.chem.agilent.com](http://www.chem.agilent.com)

Page 1 of 6


The Refractive Index Calibration Procedure

Start Refractive Index Calibration

- 1 From the user interface start the RID Calibration (see Figure 18).
 Agilent ChemStation: *Diagnosis - Maintenance - RID Calibration*
 Control Module: *System - Tests - RID - Calibrate.*
- 2 Follow the instructions with reference to the steps below.

Figure 18

Start Refractive Index Calibration

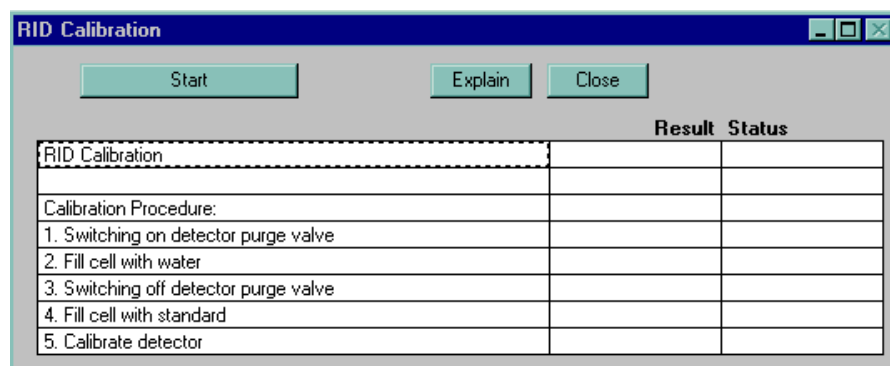

Purging the Sample and Reference Cells


- 1 The purge valve will automatically switch to the On position.
- 2 Using a syringe or LC pump flush the sample and reference cell with about 20 ml of LC grade water.
- 3 Click continue and the purge valve will automatically switch to the Off position.

Fill the Sample Cell with Calibration Solution

- 1 Remove the inlet capillary or flushing syringe from the in port.
- 2 Take the syringe and fix the needle to the syringe adapter.
- 3 Suck about 1.5 ml of the calibration sample into the syringe.
- 4 Keep the syringe in a horizontal position.
- 5 Remove the needle.
- 6 Add the filter to the syringe and fit the needle to filter.

Figure 19

Syringe with Sample Filter

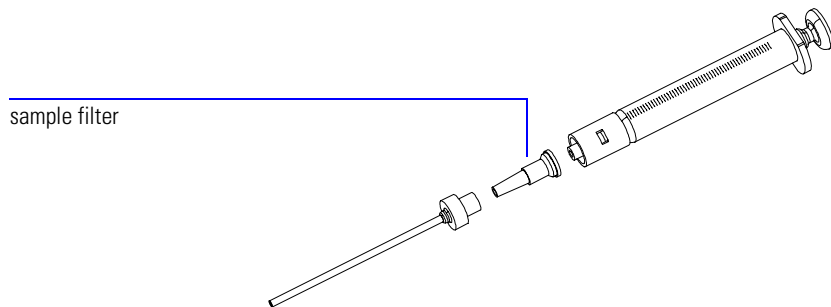

- 7 Lift the needle tip and carefully eject approximately 0.5 ml to remove air out of the syringe and to flush the needle.
- 8 Add the PEEK fitting to the needle tip and fix both at the flow cell inlet.

NOTE

Do not inject the calibration solution without the sample filter.

- 9 Slowly inject about 1.0 ml and wait for about 10 seconds to inject another 0.1 ml. This will assure that the cell is filled properly.

Calibrate Refractive Index

- 1 If the detector response differs from the theoretical response of 512,000 nRIU +/- 5,000 nRIU enter the theoretical value (512,000) in the dialog box. If the detector response is within the theoretical response click *OK* (see Figure 20).

NOTE

Rinse the sample cell with pure water at a minimum of 1.5 ml/min to flush the Sucrose from the cell and the capillaries. When organic solvent is sequentially applied (without rinsing), a blockage of capillaries may occur.

Figure 20

Calibrate Refractive Index

