

GC Firmware Update Utility

Version 2.5.5.0


Installation

1. Download the GC Firmware Update Utility setup to a folder.
2. Open the folder and double click run setup.




GC Firmware Update setup


3. Select “Yes” when asked if Setup can make changes to your device.
4. Select a language and click OK.

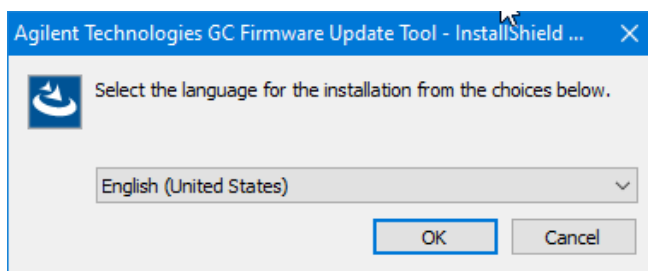


5. Click Next.

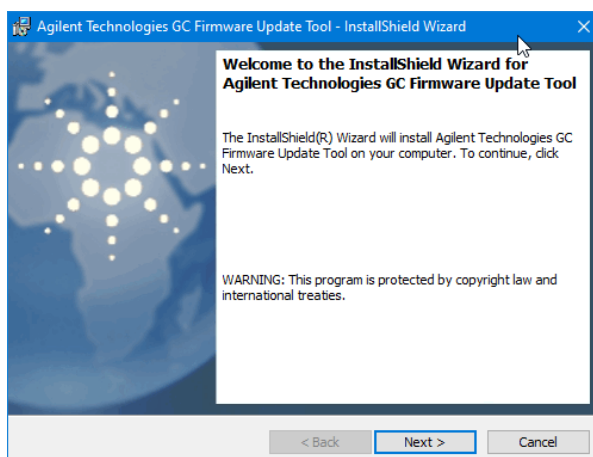


6. Accept the terms and click Next.

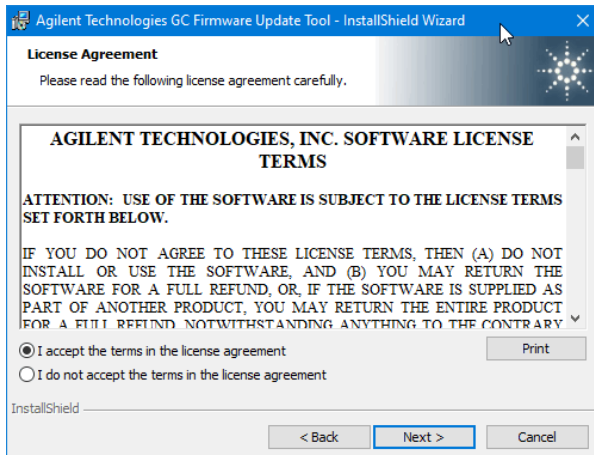

7. Click Install.

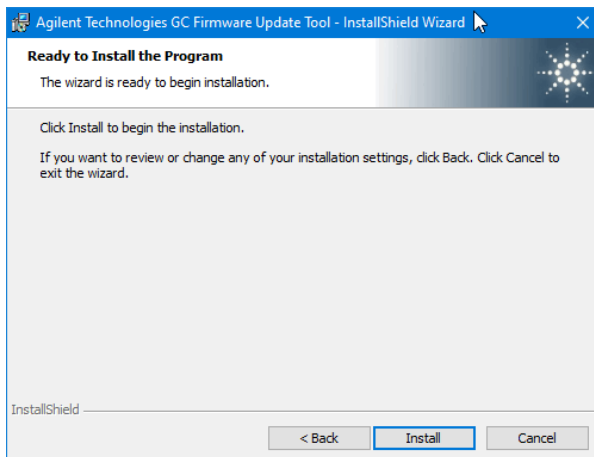

8. Click Finish to complete the installation process.

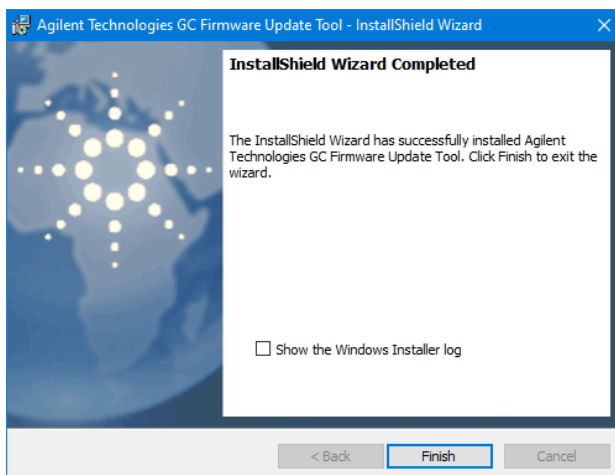

9. Installation is complete. Program icon placed on PC Desktop.


Updating Instrument Firmware

Note: In this example an 8890 Gas Chromatograph will be updated.

1. Double click on the Agilent GC Firmware Update to get started.




2. Click on the GC and ALS box.

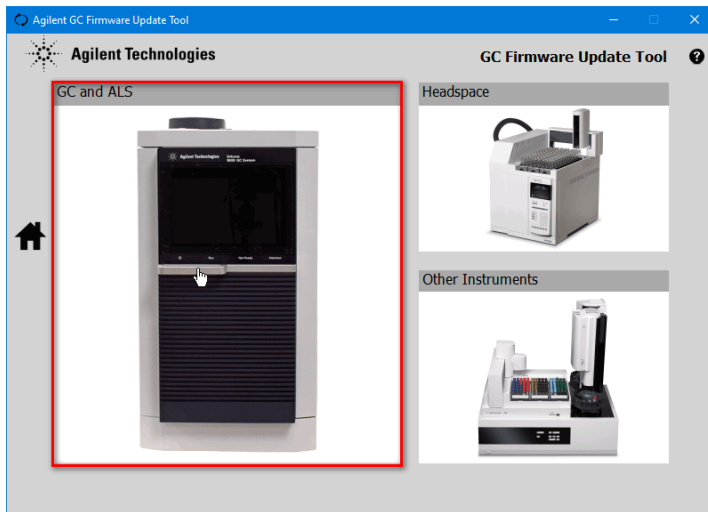


Note: To update the firmware of the Integrated 8697 Headspace Sampler you start by connecting to the GC it is connected to as well.


3. Select 8890 GC.

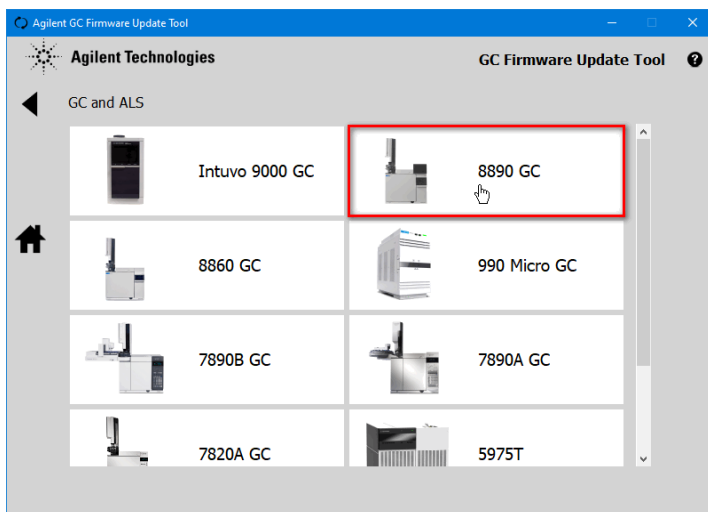


4. Select Firmware update.

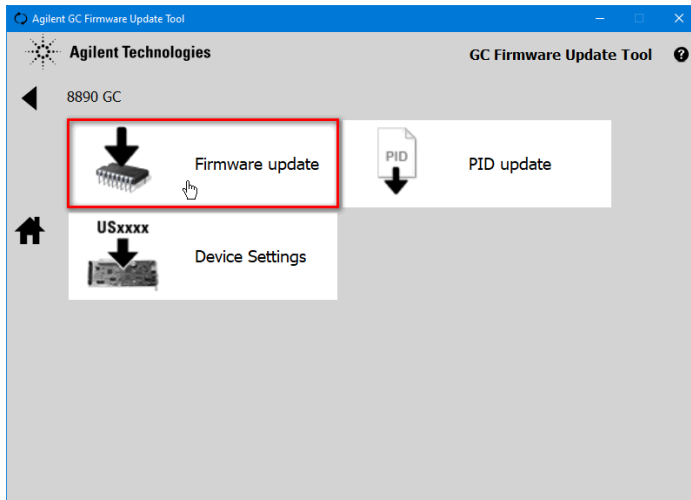


5. Input the IP address or hostname and click Connect.

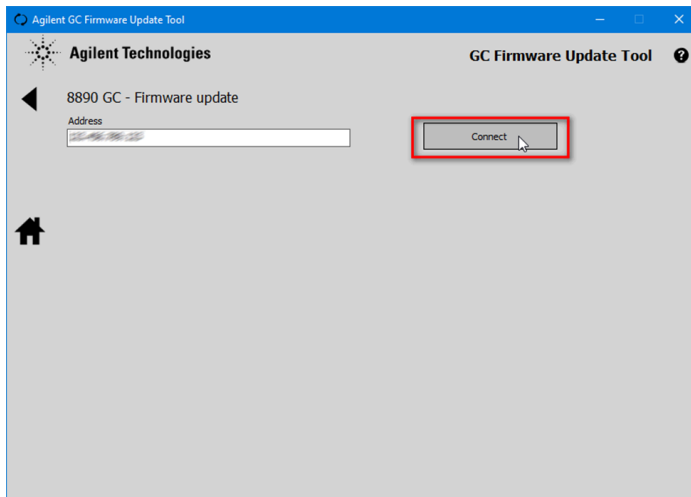

6. Select Update for the 8890 GC.

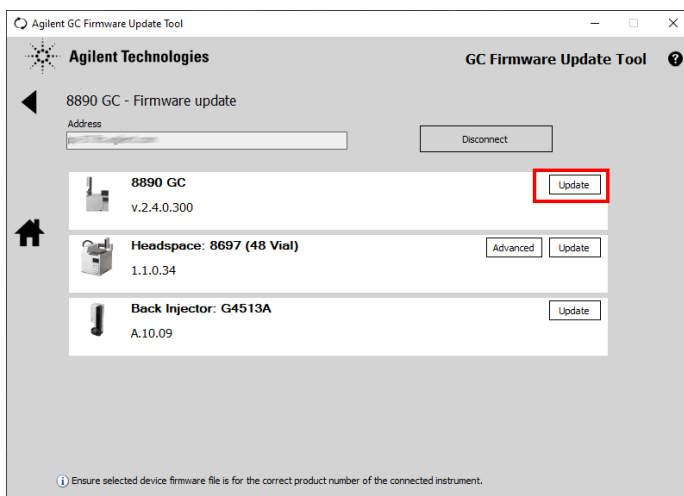


7. Click on Accept and Install.

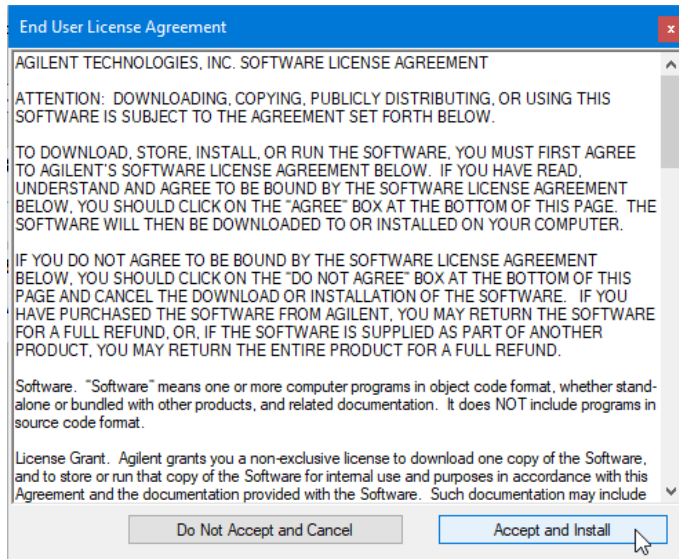

Note: You will be presented with a list of firmware files to choose from. These are files included with the update program. For the 8890 GC it is necessary to download the firmware separately.

Name	Date modified	Type
5975T	11/8/2021 9:43 AM	File folder
6820	11/8/2021 9:43 AM	File folder
6850	11/8/2021 9:43 AM	File folder
6852	11/8/2021 9:43 AM	File folder
6890	11/8/2021 9:43 AM	File folder
7667	11/8/2021 9:43 AM	File folder
7696	11/8/2021 9:43 AM	File folder
7697A	11/8/2021 9:43 AM	File folder
7820	11/8/2021 9:43 AM	File folder
7890A	11/8/2021 9:43 AM	File folder
7890B	11/8/2021 9:43 AM	File folder
8255_8355	11/8/2021 9:43 AM	File folder
8697	11/8/2021 9:43 AM	File folder
35900E	11/8/2021 9:43 AM	File folder
Frit	11/8/2021 9:43 AM	File folder
G1512A	11/8/2021 9:43 AM	File folder
G1513A	11/8/2021 9:43 AM	File folder
G1888	11/8/2021 9:43 AM	File folder
G2612A	11/8/2021 9:43 AM	File folder
G2613A	11/8/2021 9:43 AM	File folder
G2614A	11/8/2021 9:43 AM	File folder
G2880A	11/8/2021 9:43 AM	File folder
G2912A	11/8/2021 9:43 AM	File folder
G2913A	11/8/2021 9:43 AM	File folder
G2916A	11/8/2021 9:43 AM	File folder
G4513A	11/8/2021 9:43 AM	File folder
G4514A	11/8/2021 9:43 AM	File folder
G4515A	11/8/2021 9:43 AM	File folder
G4516A	11/8/2021 9:43 AM	File folder
G4567A	11/8/2021 9:43 AM	File folder
G8130A	11/8/2021 9:43 AM	File folder
Lantronix	11/8/2021 9:43 AM	File folder
pneumatic-EEPROM	11/8/2021 9:43 AM	File folder


8. Navigate to the location of the downloaded 8890 firmware and select open.

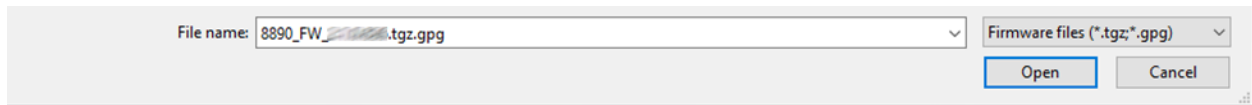

9. Once the file is selected the update will begin.

Note: Do not interrupt the update process.

10. Once the GC update is complete the instrument can be disconnected and the software shutdown unless other connected instruments need to be updated such as the ALS or Integrated Headspace sampler.

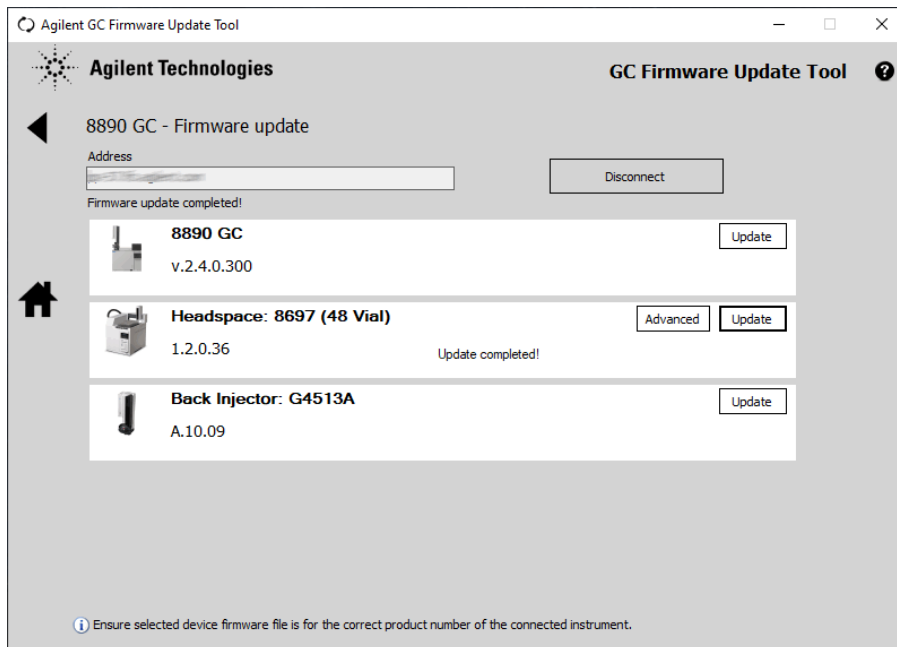

Note: In this picture the 8697 Headspace was updated along with the 8890 GC.